Metrosensor electrodes



Precision is not accidental, but brought about by design!

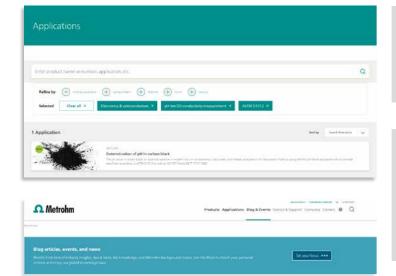


Table of contents

Which electrode for which application?	6-9	Accessories	96-102
Electrodes for pH measurement	6	Accessories for Metrosensors	96-97
Electrodes for titration	7	Ion standards, buffer solutions, electrolytes	98-99
Electrodes for USP monographs	8-9	Electrical connections	100-102
Electrode catalog	10-93	How is a Metrosensor made?	103-104
dTrodes – Digital electrodes for OMNIS	10-11		
iTrodes – Intelligent electrodes for		Theoretical fundamentals	105-126
the Titrando generation	12	1. Fundamentals of potentiometry	105-120
Unitrode and Solvotrode <i>easy</i> Clean –		1.1. Electrode construction	105
Cleaning at the touch of a button	13	1.2. From the measured potential to	
Separate pH glass electrodes	16-17	the ion concentration	105-106
Electrodes for pH measurement	18-21	1.3. Measuring electrodes	107-115
Electrodes for pH titration	22-25	1.3.1. pH glass electrodes	115-111
Special electrodes for pH measurement/pH titration	26-29	1.3.2. Metal electrodes	111-113
Electrodes for pH titration in		1.3.3. Ion-selective electrodes	113-115
HF-containing solutions	30-31	1.4. Reference electrodes	115-120
Separate metal electrodes	34-35	1.4.1. Silver/silver chloride reference electro	de 115
Titrodes – the maintenance-free metal electrodes	s 36-37	1.4.2. Metrosensor «Long Life» reference syst	em 116
Combined metal electrodes	38-39	1.4.3. Diaphragms	116-119
Electrodes for Karl Fischer titration	40-41	1.4.4. Reference electrolytes and	
Ion-selective electrodes and accessories	44-47	bridge electrolytes	120
Electrodes for surfactant titration and accessories	s 48-50	2. Fundamentals of conductometry	121-125
Reference electrodes	54-57	2.1. General	121-124
Working electrodes for trace analysis and CVS	60-61	2.2. Conductivity measurement in accordar	nce
Reference and auxiliary electrodes for		with USP and Pharm. Europe (EP)	124-125
trace analysis and CVS	62-63	3. Temperature measurement	126
Electrodes for Metrohm Autolab RDE and RRDE	64-65		
Microelectrodes	66-67	Appendix	128-141
Electrodes for the electrochemical quartz		Technical specifications	128-141
crystal microbalance	68-69		
Screen-printed electrodes	70-71	Metrohm Electrode Care	142
Interdigitated electrodes/microelectrodes	72		
Custom-made electrochemical sensors	73		
Conductivity measuring cells for			
OMNIS Measuring Module Conductivity	76-77		
Conductivity measuring cells			
for 856 Conductivity Module	78-79		
Conductivity measuring cells for 912/914 meters	80-81		
Conductivity measuring cells	82-83		
Sensors for stability measurement	84-85		
Temperature sensors	86-87		
Sensors for thermometric titration	88-89		
Optical Sensor	90-91		
Sensor for photometry	92-93		

4

Metrohm has long-standing experience in ion analysis – why should you wrack your brains, when we've already done the thinking for you?



Our webpage offers you helpful Metrohm Application Bulletins – instructions that are quaranteed to work.

You can find all applications in the Application Finder at: www.metrohm.com/applications.html

In addition, Metrohm regularly publishes blog posts on current and ubiquitous topics on the webpage.

You can find a lot of useful tips and tricks for chemical analysis, e.g., titration, Karl Fischer titration, pH measurement, and more in the Knowledge section at: www.metrohm.com/discover.html

Are you looking for a suitable electrode for your application? Then the electrode finder on our website will help you. You can refine your search using the filter settings.

Find electrodes for pH/conductivity measurement, titration, electrochemistry, and voltammetry/CVS in the Electrode Finder: www.metrohm.com/discover.html

Electrode catalog

Technical specifications

Detailed information concerning technical specifications can be found in the appendix «Technical specifications». The electrochemical parameters are specified for 25 °C, the outflow rates with a hydrostatic pressure of a 10 cm high water column.

Installation length

The «installation length» is the length from the electrode tip to the upper edge of the standard ground-joint. In the case of flexible SGJ sleeves, this corresponds to the length down to the electrolyte refill opening below. All flexible SGJ sleeves have like their size the standard ground-joint 14/15.

Shaft length

«Shaft length» refers to the length of the electrode tip up to the lower edge of the electrode plug-in head.

LL reference system

The abbreviation «LL reference system» stands for the Metrosensor «Long Life» reference system. More detailed information on this can be found in the theoretical section, Chapter 1.4.2.

Reference electrolyte

Unless otherwise specified, all combined pH electrodes are filled with the referenced electrolytes c(KCI) = 3 mol/L (Order number 6.2308.020).

DJ

«DJ» stands for «double junction». These electrodes are equipped with a bridge electrolyte chamber; the bridge electrolyte is replaceable, which means that it can be adapted to suit the sample.

Materials abbreviations:

LI	громис
EVA	Ethylenvinylacetate
PBT	Polybutylenterephthalate
PCTFE	Polychlortrifluoroethylene
PE	Polyethylene
PEEK	Polyetheretherketone
PMMA	Polymethyl metacrylate
POM	Polyoxymethylene
PP	Polypropylene
PPO	Polyphenylenoxide
PTFE	Polytetrafluoroethylene
PVC	Polyvinyl chloride

Application Specifics Electrode Order number Page Universal Standard laboratory, various samples, Unitrode 6.0258.600 pH 0...14, T = 0...100 °C 6.0221.600 Routine analysis in similar samples, pH 1...11 | Ecotrode Gel 18 Water Demineralized, drinking and sea water, Aquatrode Plus 6.0257.600 20 weakly buffered solutions Waste water 20 General Unitrode 6.0258.600 Sulfide-containing waste water Profitrode 6.0255.100 22 Soil samples Surface measurement or aqueous Flat-membrane 6.0256.100 26 suspensions Agriculture, Culture media, small volume samples Biotrode 6.0224.100 26 plant breeding Fertilizers Unitrode 6.0258.600 20 6.0255.100 22 Liquid manure Profitrode 6.0239.100 Nutrient solutions Viscotrode 28 6.0235.200 Protein-containing solutions Porotrode 26 Food, 6.0258.600 20 General Unitrode stimulants 6.0235.200 26 Protein-containing food, beer Porotrode 6.0226.100 26 Penetration measurement Spearhead (dough, cheese, meat) Drinking water Aquatrode Plus 6.0257.600 20 Juices, wine, spirits Unitrode 6.0258.600 20 Pharmaceutical Creams, liquid formulations, medicinal Viscotrode 6.0239.100 28 syrups mouthwashes, raw materials industry, 6.0258.600 biology 20 Dialysis solutions, urine Unitrode Gastric juice, serum, small-volume samples 6.0224.100 26 Biotrode Infusion solutions 6.0257.600 20 Aquatrode Plus Protein-containing solutions Porotrode 6.0235.200 26 Cosmetics Shampoos, emulsions, shower gels, Viscotrode 6.0239.100 28 lotions, perfumes Microelectrode 6.0234.100 28 Skin (surface measurement) Flat-membrane 6.0256.100 26 Detergents, General Viscotrode 6.0239.100 28 surfactants Profitrode 6.0255.100 22 Samples with pH >10 Leather, Bleaching & dye baths Profitrode 6.0255.100 22 6.0258.600 20 Dampening solutions (offset printing), glue Unitrode textile industry Leather, paper, textiles Flat-membrane 6.0256.100 26 (surface measurement) Washing liquors Viscotrode 6.0239.100 28 Paints, Dye baths, ink, wood stain, lacquers Profitrode 6.0255.100 22 lacquers, Dispersions, emulsions, resins, suspensions Unitrode 6.0258.600 20 solvents Paint (surface) 6.0256.100 Flat-membrane 26 Non-aqueous, polar solvents EtOH-Trode 6.0269.100 20 Electroplating, 6.0255.100 General Profitrode 22 6.0258.600 20 Acidic electroplating baths Unitrode processing Cutting-oil emulsions 6.0239.100 28 Viscotrode Special Concentrated acids Profitrode 6.0255.100 22 applications Photographic baths, fixative solutions Profitrode 6.0255.100 22 6.0258.600 20 Emulsions, suspensions, dispersions Unitrode Polymer dispersions (water-based dispersion Ecotrode Gel 6.0221.600 18 paints and adhesive dispersions on basis of acrylic acid esters and styrene) Samples at pH >12 Unitrode 6.0258.600 20 Temperature 80...100 °C Unitrode (Idrolyte) 6.0258.600 20 20 Ion-deficient solutions, weakly buffered Aquatrode Plus 6.0257.600 Non-aqueous, polar solvents EtOH-Trode 6.0269.100 20 Penetration measurement Spearhead 6.0226.100 26 6.0235.200 Protein-containing solutions Porotrode 26 6.0224.100 Small volume samples Biotrode 26 Flat-membrane 6.0256.100 26 Surface measurements Flat-membrane 6.0256.100 26 Biofuels EtOH-Trode 6.0269.100 20

Electrodes for Titration

Application	Specifics	Electrode	Order number	Page
Aqueous acid/	General	Ecotrode Plus	6.0262.100	22
base titrations		Ecotrode Gel	6.0221.100	22
	Alkaline samples, Bayer liquors. Titrations performed at high temperatures	Unitrode	6.0259.100	22
	Acidity of alcoholic beverages	Unitrode easyClean with Pt1000	6.0260.010	13
	Titrations with small sample volumes	Flat membrane pH electrode	6.0256.100	26
	Titrations in water having a low conductivity	Aquatrode Plus	6.0253.100	24
	Carbonate hardness, acid capacity of water, p- & m-value	Aquatrode Plus with Pt1000	6.0257.600	20
	Electroplating baths, etching baths, phosphate coating	Profitrode	6.0255.100	22
	Electroplating and etching baths containing fluoride or hydrofluoric acid	Solitrode HF	6.0223.100	30
	Protein-containing samples	Porotrode	6.0235.200	26
Non-aqueous acid/base titrations	Determination of the base number (TBN) of mineral oils, titrations with perchloric acid, cyclohexylamine, alcoholic HCl	Solvotrode easyClean with LiClsat in ethanol	6.0229.010	13
	Determination of the acid number (TAN) of mineral oils, titrations with alcoholic KOH, NaOH and TBAOH, potassiummethylate, free acids; hydroxyl number of oils and fats	Solvotrode <i>easy</i> Clean c(TEABr) = 0.4 mol/L in ethylene glycol)	6.0229.010 (+ 6.2320.000)	13
Redox titrations	Titrations at constant pH value	Pt-Titrode	6.0431.100	36
Arsenite, cerium	Titrations with change of the pH value	Combined Pt-ring-electrode	6.0451.100	38
sulfate, iron(III), odine, potassium bromate, sodium	Chemical oxigen demand in water (ferrometry)	Combined Au-ring electrode	6.0452.100	38
nitrite, oxalic acid,	Penicillin, ampicillin	Combined Au-ring electrode	6.0452.100	38
permanganate,	Titrations in I _{pol} mode	Double Pt sheet electrode	6.0309.100	34
thiosulfate, titan(III), Hg(NO ₃) ₂	Bromatometry, iodometry, cerimetry according to Ph. Eur. & USP	Pt-Titrode	6.0431.100	36
Karl Fischer Titrations	Water content according to Karl Fischer	Double Pt-wire electrode (HF-resistant version on request)	6.0338.100	40
Complexometry Titrations with	Back-titration of the Ba2+ excess with EDTA	Combined Ca ²⁺ -ISE polymer membrane	6.0510.100	44
EDTA, Complexon [©] II and IV	Bulletin 125)	Combined Ca ²⁺ -ISE polymer membrane	6.0510.100	44
	Determination of Al, Ba, Bi, Ca, Cd, Co, Fe, Mg, Ni, Pb, Zn (Application Bulletin 101)	Cu ²⁺ ISE – crystal membrane	6.0502.140	44
Precipitation	Chloride in general, sodium chloride in food	Ag-Titrode	6.00430.100	36
titrations Fitrations with	Chloride in dialysis and infusion solutions	Ag-Titrode with Ag ₂ S-coating	6.00430.1005	36
silver nitrate	Titrations according to Ph Eur. & USP	Ag-Titrode with Ag ₂ S-coating	6.00430.1005	36
	Determination of hydrogen sulfide, mercaptans, carbonyl sulfide, sulfides	Ag-Titrode with Ag₂S-coating	6.00430.1005	36
	Chloride, bromide, iodide and cyanide in electroplating baths	Ag-Titrode with Ag₂S-coating	6.00430.1005	36
		FISE crystal membrane	6.0502.150	44
Photometric titrations	Titration in aqueous and non-aqueous media; choice of eight wave lengths (470, 502, 520, 574, 590, 610, 640, 660 mn)	Optrode	6.1115.000	92
Surfactants in non-aqueous media Aromatic/aliphatic,	Titration of anionic and cationic surfactants, titrations in chloroform, surfactant formulations containing oils, pH <10	Surfactrode Resistant	6.0507.130	48
hydrocarbons, ketones, chloroethanes, methylisobutylk- etone	Titration of anionic and cationic surfactants, titration of washing agents, soaps, pH >10	Surfactrode Refill	6.0507.140	48
Surfactants in	Titration of cationic surfactants	Cationic Surfactant	6.0507.150	48
aqueous media	Titration of anionic surfactants	Ionic Surfactant	6.0507.120	48
	Titration of non-ionic surfactants, titration of pharmaceutical ingredients with sodium tetraphenylborate	NIO electrode	6.0507.010	48
Thermometric titrations	Titrations in aqueous and non-aqueous solutions without HF	Thermoprobe	6.9011.020	88
	Titrations in aqueous solutions containing HF	Thermoprobe HF	6.9011.040	88

Electrodes for USP monographs

8

Titrant	Specifics	Electrode Order no.		Page
Acetic acid	All indicators	dEcotrode plus 6.00201.300		10
Ammonium thiocyante	Ferric ammonium sulfate, sample contains silver or silver nitrate VS is added in excess	dAg Titrode	6.00404.300S	10
	Ferric ammonium sulfate, sample contains mercury	Comb. Au ring electrode	6.0452.100	38
	Silver electrode and double-junction Sep. Ag ring electrode reference electrode		6.00350.100S	34
	Other electrodes	dAg Titrode	6.00404.300S	10
Barium perchlorate	All indicators	Cu ²⁺ ISE	6.0502.140	44
Benzethonium chloride	Methylene blue	Surfactrode Resistant	6.0507.130	48
Bismuth nitrate	All indicators	Cu ²⁺ ISE	6.0502.140	44
Bromine	All indicators	dPt Titrode	6.00401.300	10
Cerric ammonium nitrate, cerric	All color indicators, combined platinium electrode, potentiometric indication	dPt Titrode	6.00401.300	10
ammonium sulfate, cerric sulfate	Electrodes with silver-silver chloride or calomel reference system	dPt-ring-electrode	6.00403.300	10
Cupric nitrate	All indicators	Cu ²⁺ ISE	6.0502.140	44
Dichlorophenol- Indophenol	Dichlorophenol - All indicators Double Pt sheet electrode or		6.0309.100 or 6.00353.100	34
Edetate disodium (EDTA)	Hydroxy naphthol blue, calconcarboxylic acid triturate, or calcium ion-selective electrode	dCa ²⁺ ISE	6.00502.300	10
	Other potentiometric indication	Cu ²⁺ ISE	6.0502.140	44
	Other color indicators	Optrode	6.1115.000	92
Ferric and ferrous ammonium sulfate	All indicators	dPt Titrode	6.00401.300	10
Hydrochloric acid	Aqueous titrant, all indicators, solvent is water	dEcotrode plus	6.00201.300	10
	Non-aqueous titrant, all indicators, non-aqueous solvent	dSolvotrode	6.00203.300	10
Iodine	All indicators	dPt Titrode	6.00401.300	10
Lead nitrate and lead perchlorate	All indicators	Pb ²⁺ ISE	6.0502.170	44
Lithium methoxide	All indicators	dSolvotrode	6.00203.300	10
Magnesium	Eriochrome black-T-sodium chloride	Optrode	6.1115.000	24
chloride	Copper ion-selective electrode	Cu ²⁺ ISE	6.0502.140	44
Mercuric nitrate	All indicators	Comb. Au ring electrode	6.0452.100	26
Oxalic acid	All indicators	dPt Titrode	6.00401.300	10
Perchloric acid	All color indicators, all electrodes except below	dSolvotrode	6.00203.300	10
	Electrodes with lithium chloride or lithium per- chlorate in glacial acetic acid as electrolyte	Special double junction Solvotrode		
Potassium arsenite	All indicators	dPt Titrode	6.00401.300	10
Potassium bromate	All indicators	dPt Titrode	6.00401.300	10
Potassium bro- mide - potassium bromate	All indicators	Double Pt-wire electrode for coulometry	6.0341.100	40

Electrodes for USP monographs

Potassium All indicators All indicators All indicators All indicators All indicators All indicators All indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous solvent All indicators	Titrant	Specifics	Electrode	Order no.	Page
Potassium Aqueous titrant, all indicators, solvent is water Aqueous titrant, all indicators, non-aqueous disolvotrode 6.00201.300 10		All indicators	dPt Titrode	6.00401.300	10
Non-aqueous titrant, all indicators, non-aqueous solvent Solvent of the potassium iodate Potassium iodate Potassium All indicators		All indicators	dPt Titrode	6.00401.300	10
Potassium iodate All indicators Silver electrode All indicators Silver electrode Silver electrode All indicators All indicator				6.00201.300	10
Potassium permanganate Potassium permanganate Potassium Potassium permanganate All indicators All indicators Silver electrode and double-junction reference electrode Silver electrode and double-junction reference electrode Sep. Ag ring electrode 6.00350.100S 34	hydroxide		dSolvotrode	6.00203.300	10
Potassium thiocyanate Silver electrode and double-junction reference electrode Silver electrode and double-junction reference electrode Silver electrode and double-junction reference electrode Silver nitrate Potassium chromate is used as indicator All other color indicators, potentiometric indication with all electrodes, except listed below Electrodes with silver-silver chloride or calomel reference system Silver-ion selective electrode Agr/5²- ISE 6.0502.180 44 Silver-ion selectrode and double-junction Sep. Agring electrode 6.00402.3005 10 Silver-ion selectrode and double-junction reference electrode Agr/5²- ISE 6.0502.180 44 Silver-ion selectrode and double-junction reference electrode All indicators described All indicators Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous doolvent Starch, platinum-calomel electrodes dP ring electrode 6.00203.300 10 Solvent Platinum-platinum electrodes Double Pt sheet electrode 6.0309.100 34 Solium nitrite Starch, platinum-calomel electrodes Double Pt sheet electrode 6.0507.010 48 Combined potassium ion selective electrode Rr ISE 6.0510.110 44 Solium thiosulfate Electrodes with calomel reference electrode Agreous titrant, all indicators, non-aqueous dool of the properties of the proper	Potassium iodate	All indicators	dPt Titrode	6.00401.300	10
Silver electrode and double-junction reference electrode February		All indicators	dPt Titrode	6.00401.300	10
Potassium chromate is used as indicator All other color indicators, potentiometric indication with all electrodes, except listed below Electrodes with silver-silver chloride or calomel reference system Silver-ion selective electrode Agr/S²-ISE 6.0502.180 44			dAg Titrode	6.00404.300S	10
All other color indicators, potentiometric indication with all electrodes, except listed below Electrodes with silver-silver chloride or calomel reference system Silver-ion selective electrode Sodium arsenite All indicators Aqueous titrant, all indicators, non-aqueous Sodium nitrite Sodium nitrite Sodium ethoxide Sodium ethoxide Sodium hydroxide Sodium hydroxide Sodium hydroxide Sodium nitrite Sodium ethoxide S	thiocyanate		Sep. Ag ring electrode	6.00350.100S	34
tion with all electrodes, except listed below Electrodes with silver-silver chloride or calomel reference system Silver-ion selective electrode Ag*/5²- ISE 6.0502.180 44 Silver electrode and double-junction reference electrode Ag*/5²- ISE 6.0502.180 Ad Silver electrode and double-junction reference electrode Sodium arsenite All indicators Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous double publication solvent Sodium methoxide All indicators Sodium nitrite Starch, platinum-calomel electrodes Ag* obvolved Sodium electrode Sodium hydroxide All indicators Sodium electrodes Sodium electrodes Sodium electrode Sodium electrodes Sodium electrode Sodium electrodes Sodium electrode Sodium electrodes Solosoficators Sodium electrodes Solosoficators Sodium electrodes Sodium electrodes Sodium electrode	Silver nitrate	Potassium chromate is used as indicator	dAg ring electrode	6.00402.300S	10
calomel reference system Silver-ion selective electrode Silver electrode and double-junction reference electrode Sep. Ag ring electrode Ag*/S²- ISE 6.0502.180 44 Silver electrode and double-junction reference electrode Sep. Ag ring electrode 6.00350.100S 34 Sep. Ag ring electrode 6.00350.100S 34 Sep. Ag ring electrode 6.00401.300 10 Sodium hydroxide Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous solvent Sodium methoxide All indicators All indicators Asilver ring electrode All indicators All indicators Asilver ring electrode All indicators All indicators Asilver ring electrode Asilver ring ring ring ring ring ring ring rin			dAg Titrode	6.00404.300S	10
Silver electrode and double-junction reference electrode Sep. Ag ring electrode 6.00350.100S 34			dAg ring electrode	6.00402.300S	10
Sodium arsenite All indicators All indicators, solvent is water All occupant strant, all indicators, solvent is water All occupant strant, all indicators, solvent is water All occupant strant, all indicators, solvent is water All occupant solvent All indicators				6.0502.180	44
Sodium hydroxide Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous dSolvotrode 6.00203.300 10 10 10 10 10 10 10			Sep. Ag ring electrode	6.00350.100S	34
Non-aqueous titrant, all indicators, non-aqueous dSolvotrode 6.00203.300 10	Sodium arsenite	All indicators	dPt Titrode	6.00401.300	10
Sodium methoxide All indicators All indicators Starch, platinum-calomel electrodes All tring electrode 6.00203.300 10	Sodium hydroxide			6.00201.300	10
Sodium nitrite Starch, platinum-calomel electrodes Double Pt sheet electrode 6.00403.300 10			dSolvotrode	6.00203.300	10
Platinum-platinum electrodes Double Pt sheet electrode 6.0309.100 34 Eterraphenylboron Sodium thiosulfate All color indicators and electrodes except with calomel reference system Electrodes with calomel reference electrode Non-aqueous titrant, all indicators, non-aqueous solvent Tetrabutylammonium hydroxide Tetramethylammonium bromide Tetramethylammonium chloride Titanium trichloride All color indicators All indicator	Sodium methoxide	All indicators	dSolvotrode	6.00203.300	10
Bromophenol blue solution (1 in 2000) NIO surfactant electrode 6.0507.010 48	Sodium nitrite	Starch, platinum-calomel electrodes		6.00403.300	10
tetraphenylboron Combined potassium ion selective electrode All color indicators and electrodes except with calomel reference system Electrodes with calomel reference electrode Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous solvent Tetrabutylammonium hydroxide Tetramethylammonium bromide Tetramethylammonium chloride Titanium trichloride All indicators Combined potassium ion selective electrode K* ISE 6.0510.110 44 APT Titrode 6.00401.300 10 6.00201.300 10 6.00203.300 10 6.00203.300 10 6.00203.300 10 6.00404.300S		Platinum-platinum electrodes	Double Pt sheet electrode	6.0309.100	34
Sodium thiosulfate All color indicators and electrodes except with calomel reference system Electrodes with calomel reference electrode dPt ring electrode 6.00401.300 10		Bromophenol blue solution (1 in 2000)	NIO surfactant electrode	6.0507.010	48
calomel reference system Electrodes with calomel reference electrode dPt ring electrode 6.00403.300 10 Sulfuric acid Aqueous titrant, all indicators, solvent is water dEcotrode plus 6.00201.300 10 Non-aqueous titrant, all indicators, non-aqueous solvent dSolvotrode 6.00203.300 10 Tetrabutylammonium hydroxide Tetramethylammonium bromide Tetramethylammonium chloride Titanium trichloride All indicators dPt Titrode 6.00404.300S 10 Zinc sulfate All color indicators Optrode 6.1115.000 92				6.0510.110	
Sulfuric acid Aqueous titrant, all indicators, solvent is water Non-aqueous titrant, all indicators, non-aqueous Non-aqueous titrant, all indicators, non-aqueous Solvent Solv	Sodium thiosulfate	calomel reference system	dPt Titrode	6.00401.300	10
Non-aqueous titrant, all indicators, non-aqueous dSolvotrode Tetrabutylammo- nium hydroxide Tetramethylammo- nium bromide Tetramethylammo- nium bromide Tetramethylammo- nium chloride Titanium trichloride All indicators dAg Titrode dAg Titrode dAg Titrode 6.00404.300S 10 dAg Titrode 6.00404.300S 10 dAg Titrode 6.00404.300S 10 dAg Titrode 6.00401.300 10 dag Titrode		Electrodes with calomel reference electrode	dPt ring electrode	6.00403.300	10
Solvent Solv	Sulfuric acid		l l	6.00201.300	10
nium hydroxide dag Titrode 6.00404.300S 10 Tetramethylammonium bromide All indicators dag Titrode 6.00404.300S 10 Tetramethylammonium chloride All indicators dag Titrode 6.00404.300S 10 Titanium trichloride All indicators description description 6.00401.300 10 Zinc sulfate All color indicators Optrode 6.1115.000 92			dSolvotrode	6.00203.300	10
nium bromide dAg Titrode 6.00404.300S 10 Tetramethylammonium chloride All indicators dAg Titrode 6.00401.300 10 Titanium trichloride All indicators dPt Titrode 6.00401.300 10 Zinc sulfate All color indicators Optrode 6.1115.000 92		All indicators	dSolvotrode	6.00203.300	10
nium chloride But the control of the cont		All indicators	dAg Titrode	6.00404.3005	10
Zinc sulfateAll color indicatorsOptrode6.1115.00092		All indicators	dAg Titrode	6.00404.3005	10
	Titanium trichloride	All indicators	dPt Titrode	6.00401.300	10
Copper ion-selective electrode Cu ²⁺ ISF 6.0502 140 44	Zinc sulfate	All color indicators	Optrode	6.1115.000	92
CODE TOT SCIENCE CICCIONC CONTROL TO		Copper ion-selective electrode	Cu ²⁺ ISE	6.0502.140	44

dTrodes – Digital electrodes for OMNIS

10 dTrodes – simply easier

The latest generation of intelligent, digital electrodes – the dTrodes – were specifically developed for the OMNIS platform. The dTrodes combine the measuring electrode and the measuring amplifier in one slim and robust sensor.

The electrode head contains an analog-to-digital converter, which converts the measured analog signal into a binary code. The digital data is then optically transmitted to the titrator. This means, the sensor is electrostatically decoupled from the titrator. The signal is thus not affected by electrostatic interferences and is therefore of low noise.

Important sensor data, such as article and serial numbers, calibration data, calibration history, working life, and calibration validity period are automatically saved on the integrated memory chip. Therefore, the sensor can easily be used on another instrument without having to repeat the calibration. This also eliminates the danger of mixing up sensors.

LED for sensor status

When connecting a dTrode to the OMNIS titrator, the sensor is automatically recognized by the software and saved in its sensor list. Immediately, all sensor data is transferred to the software and checked for its validity. A problem with the sensor, e.g., an invalid parameter, can be recognized immediately by the color of the integrated LED. Sensor management at its best!



Ordering information	
dUnitrode with Pt1000	6.00200.300
dEcotrode plus	6.00201.300
dAquatrode plus with Pt1000	6.00202.300
dSolvotrode	6.00203.300
dProfitrode	6.00204.300
dAg ring electrode with Ag₂S coating	6.00402.300S
dAg-Titrode	6.00404.300
dAg-Titrode with Ag₂S coating	6.00404.300S
dPt-Titrode	6.00401.300
dAg ring electrode	6.00402.300
dPt ring electrode	6.00403.300
dFluoride-ISE with Pt1000, combined	6.00500.300
dCalcium-ISE, combined	6.00502.300
dThermoprobe	6.01117.300
dThermoprobe HF	6.01118.300
Digital measuring module	6.02100.010



iTrodes – Intelligent electrodes for the Titrando generation

12 Easy digital identification

The iTrodes contain a memory chip which enables the storage of all relevant sensor data, such as article and serial number, calibration data, calibration history, working live and calibration validity period.

All of the sensor data are uploaded automatically when the iTrode is connected to an instrument. This means, that the possibility of any mix-up or editing error is eliminated and each analysis result is traceable to the used electrode.

Digital data transmission

iPt-ring electrode, combined

iConnect

When connected to the sensor, the analog/digital converter in the 854 iConnect (2.854.0010) converts the analog measuring signal into binary code. Digital data transmission means that the measuring signal is no longer susceptible to electrostatic influences. Interference-free transmission can now be guaranteed, no matter how long the electrode cable is.

iTrodes can be used with the 913 and 914 pH meter, the 916 Ti-Touch or the 888 and 90x Titrandos.

Ordering information iAquatrode Plus with Pt1000 6.0277.300 iUnitrode with Pt1000 6.0278.300 iSolvotrode 6.0279.300 iEcotrode Plus 6.0280.300 6.00470.300 iAg-Titrode iAg-Titrode with Ag₂S coating 6.00470.3005 iPt-Titrode 6.0471.300 iAg-ring electrode, combined 6.00450.300

6.0451.300 2.854.0010

Unitrode and Solvotrode *easy*Clean – Cleaning at the touch of a button

Easy cleaning of the diaphragm: Just press once on the electrode head and the electrolyte flows out. The diaphragm does not need to be touched any more for cleaning.

Highlights

- Easy contact-free cleaning of the diaphragm
- Greater accuracy and reproducibility of the electrolyte flow (glass on glass, the spring in the head returns to the defined starting point)
- Reduced immersion depth of the sensor through optimization of the membrane shape



Ordering information	
Unitrode <i>easy</i> Clean with Pt1000, fixed cable, plug F	6.0260.010
Unitrode <i>easy</i> Clean with Pt1000, fixed cable 2 m, plug F	6.0260.020
Solvotrode <i>easy</i> Clean, fixed cable, plug F	6.0229.010
Solvotrode <i>easy</i> Clean, fixed cable 2 m, plug F	6.0229.020



Electrodes for pH measurement/pH titration



Fine-tune your measurements!

The greatest precision and ease of care – these are the two outstanding properties of Unitrode and Aquatrode Plus. The constant electrolyte outflow of the fixed ground-joint diaphragm (which is largely insensitive to contamination) guarantees a low-noise measuring signal, even in difficult samples and independent of the measuring conditions. Further details can be found in the theoretical section on page 118.

Separate pH glass electrodes

16 Separate pH glass electrode

- Electrically shielded
- Blue T glass for reliable results, e.g. in differential potentiometry in non-aqueous media
- Optimal length for sample changer applications

Technical specifications

pH range 0...14
Temperature range 0...80 °C
Installation length 142 mm
Shaft diameter 12 mm
Minimum immersion depth 15 mm

Electrode plug-in head Metrohm plug-in head G

Differential potentiometry

In addition to the measuring electrode, a reference electrode and an auxiliary electrode are required for differential potentiometry. The shielding of the reference electrode must be identical to that of the measuring electrode.

Reference electrodes for differential potentiometry (see «Reference electrodes» section)

• Ag/AgCl DJ reference electrode, length 100 mm, Metrohm plug-in head G Without electrolyte filling, without cable 6.0729.100

• Ag/AgCl DJ reference electrode, length 138 mm, Metrohm plug-in head G Without electrolyte filling, without cable 6.0729.110

Auxiliary electrodes for differential potentiometry, Metrohm plug-in head B (see «Separate metal electrodes» section)

Separate Pt wire electrode 6.0301.100

Separate Pt rod electrode 6.1241.040 + 6.1248.000



Ordering information

Separate pH glass electrode, without cable

6.0150.100

Electrodes for pH measurement

Solitrode with Pt1000 – robust and reliable, ideal for routine laboratory use

- For solutions that do not contain precipitates, proteins or sulfides
- Long-lasting standard electrode
- Unbrekable plastic shaft

18

- Impact protection for the glass membrane
- LL reference system with long-term stability

Technical specifications

Solitrode

Shaft material PP
pH range 0...14
Temperature range 0...80 °C
Temperature sensor Pt1000
Diaphragm Ceramic pin
Installation length 113 mm
Shaft diameter 12 mm
Minimum immersion depth 15 mm

Ecotrode Gel – the maintenance-free solution

- Ideal for routine measurements in similar samples
- Specialist for measurement in polymer dispersions (water-based dispersion paints and adhesive dispersions on basis of acrylic acid esters and styrene; butyl acrylates)
- Maintenance-free
- Lifetime indicator
- LL reference system with long-term stability
- Variant 6.0228.030 with waterproof plug I for use with the 913/914 pH meter (IP67)

Technical specifications

Ecotrode Gel

Shaft material Glass
pH range 1...11
Temperature range 0...60 °C
Temperature sensor Pt1000
Diaphragm Twin pore
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head U

How to store your electrodes correctly:

Rapid response is not a matter of magic, but rather a question of storage! Metrohm recommends the patented storage solution (6.2323.000) for all combined pH glass electrodes which use c(KCI) = 3 mol/L as the reference electrolyte. This prevents the aging of the glass membrane and, as a result, guarantees response times short as they were on the first day. More information on this can be found in the theoretical section in Chapter 1.3.1. «pH glass electrodes.»







6.0228.030

6.0220.100

6.00221.600

6.0221.100

Ordering information

Solitrode with Pt1000, fixed cable (1.2 m) with plug I (IP67) + 1 x B (2 mm)

Solitrode without temperature sensor, without cable

Ecotrode Gel with Pt1000, without cable, plug-in head U

Ecotrode Gel without temperature sensor, without cable

Unitrode with Pt1000 - high performance in

difficult samples and at high pH values

- For universal use, even in dyes, pigments, inks, suspensions, resins and polymers
- Fixed ground-joint diaphragm insensitive to contamination
- High temperature resistance and very low alkali error
- Rapid response to temperature changes

20

- Outer electrolyte Idrolyte for temperatures of 80...100 °C
- LL reference system with long-term stability

Technical specifications Unitrode

Shaft material pH range

Temperature range

Temperature sensor

Glass 0...14 0...100 °C Pt1000

Diaphragm Fixed ground-joint

Installation length 125 mm Shaft diameter 12 mm Minimum immersion depth 25 mm

Aquatrode Plus with Pt1000 - ideal for weekly buffered aqueous solutions

- Special electrode membrane glass: precise measuring values and very rapid response times, even in weekly buffered solutions such as drinking water, surface water and rain water and other poorly conducting solutions
- Maintenance-free inner reference electrolyte (gel)
- Variable bridge electrolyte for special applications
- Fixed ground-joint diaphragm insensitive to contamination
- Optimized length for sample changer applications
- LL reference system with long-term stability

EtOH-Trode - the specialist for ethanol

- Developed for pHe measurement in ethanol
- Special membrane glass
- Very precise ground-joint diaphragm
- Double-junction system for free choice of electrolytes (e.g. 3 M KCI in ASTM D 6423, 1 M LiCI in EN
- LL reference system with long-term stability

Technical specifications Aquatrode Plus

Shaft material Glass pH range 0...13 Temperature range 0...60 °C Pt1000 Temperature sensor

Diaphragm Fixed ground-joint

Installation length 135 mm Shaft diameter 12 mm Minimum immersion depth 20 mm

Technical specifications

EtOH-Trode

Glass Shaft material pH range 0...13 0...80 °C Temperature range Fixed ground-joint

Installation length 125 mm

Shaft diameter 12 mm Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Diaphragm

15490).

Wellness for the electrode

Reliable measuring results over long periods of time can only be guaranteed if the glass membrane and the diaphragm receive preventative and regular care. Cleaning by means of etching with toxic chemicals or a mechanical treatment of the diaphragm is not only complicated and expensive, it also accelerates the aging of the pH glass electrode as well. The care kit (6.2325.000, pg. 98) was developed for simple, gentle cleaning of pH glass electrodes with a liquid electrolyte. Regular application can considerably prolong its lifetime.



Ordering information

Unitrode easyClean with Pt1000, fixed cable (1.2 m) plug F + 1 x B (2mm) 6.0260.010 Unitrode easyClean with Pt1000, fixed cable (2 m) plug F + 1 x B (2 mm) 6.0260.020 Unitrode without temperature sensor, without cable 6.0259.100 Unitrode with Pt1000, without cable, plug-in head U 6.0258.600 iUnitrode with Pt1000 6.0278.3001 dUnitrode with Pt1000 $6.00200.300^2$ Aquatrode Plus with Pt1000, without cable, plug-in head U 6.0257.600 Aquatrode Plus without temperature sensor, without cable 6.0253.100 iAquatrode Plus with Pt1000 $6.0277.300^{1}$ dAquatrode Plus with Pt1000 6.00202.3002 EtOH-Trode without temperature sensor, without cable 6.0269.100

¹ Further information about iTrodes can be found on pg. 12.

² Further information about dTrodes can be found on pg. 10.

Electrodes for pH titration

Ecotrode Gel - the maintenance-free solution

- Ideal for routine measurements in similar samples
- Maintenance-free

22

- Lifetime indicator
- LL reference system with long-term stability

Technical specifications

Ecotrode Gel

Shaft material Glass
pH range 1...11
Temperature range 0...60 °C
Diaphragm Twin pore
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Ecotrode Plus – high durability in routine use at a fair price

- For acid/base titrations in various kinds of solutions
- Fixed ground-joint diaphragm insensitive to contamination
- Ideal for routine laboratory use
- LL reference system with long-term stability

Technical specifications

Ecotrode Plus

Shaft material Glass
pH range 0...13
Temperature range 0...80 °C
Diaphragm Fixed ground-joint

Installation length 125 mm Shaft diameter 12 mm Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Profitrode – professional working in the most difficult of matrices

- For difficult matrices (galvanic baths, precipitates, samples containing sulfides, etc.)
- Flexible ground-joint diaphragm, particularly easy to clean
- Double-junction construction
- Available in various lengths (113/170/310 mm)
- LL reference system with long-term stability

Technical specifications Profitrode

Shaft material Glass pH range 0...14
Temperature range 0...80 °C

Diaphragm Flexible ground-joint Installation length 113/170/310 mm

Shaft diameter 12 mm Minimum immersion depth 30 mm

Electrode plug-in head Metrohm plug-in head G

Unitrode – high performance in difficult samples and at high pH values

- For universal use, even in dyes, pigments, inks, suspensions, resins and polymers
- Fixed ground-joint diaphragm insensitive to contamination
- High temperature resistance and very low alkali error
- Rapid response to temperature changes
- Outer electrolyte Idrolyte for temperatures of 80...100 °C
- LL reference system with long-term stability

Technical specifications

Unitrode

Shaft material Glass
pH range 0...14
Temperature range 0...100 °C
Diaphragm Fixed ground-joint
Installation length 125 mm

Shaft diameter 12 mm Minimum immersion depth 25 mm



Ordering information	
Ecotrode Gel without temperature sensor, without cable	6.0221.100
Ecotrode Plus, without temperature sensor, without cable	6.0262.100
iEcotrode Plus without temperature sensor	6.0280.300 ¹
dEcotrode Plus, without temperature sensor	$6.00201.300^2$
Profitrode, length 113 mm, without cable	6.0255.100
Profitrode, length 170 mm, without cable	6.0255.110
Profitrode, length 310 mm, without cable	6.0255.120
dProfitrode, without temperature sensor	6.00204.3002
Unitrode easyClean with Pt1000, fixed cable (1.2 m) plug F + 1 x B (2 mm)	6.0260.010
Unitrode easyClean with Pt1000, fixed cable (2 m) plug F + 1 x B (2 mm)	6.0260.020
Unitrode without temperature sensor, without cable	6.0259.100
Unitrode with Pt1000, without cable, plug-in head U	6.0258.600
iUnitrode with Pt1000	6.0278.300 ¹
dUnitrode with Pt1000	$6.00200.300^2$

¹ Further information about iTrodes can be found on pg. 12.

² Further information about dTrodes can be found on pg. 10.

Aquatrode Plus – ideal for aqueous, weekly

buffered solutions

24

- Precise measuring values and very rapid response times in ion-deficient or weekly buffered solutions – such as drinking water, surface water and rain water – thanks to special membrane glass and optimized, fixed ground-joint diaphragm insensitive to contamination
- Maintenance-free inner reference electrolyte (gel)
- Variable bridge electrolyte for special applications
- LL reference system with long-term stability

Solvotrode easyClean – space-saving alternative for titration in non-aqueous media

- For non-aqueous titrations in the pharmaceutical sector
- For determination of TAN/TBN in compliance with ASTM D4739, D2896 and D664 and DIN ISO 3771 and DIN EN 12634
- Reference electrolyte: LiCl(sat) in ethanol
- Rapid response and stable measuring values in organic solvents
- Shielding against electrostatic effects
- EasyClean diaphragm, particularly easy and contactless cleaning
- $\bullet\,$ LL reference system with long-term stability

Technical specifications Aquatrode Plus

Shaft material Glass
pH range 0...13
Temperature range 0...60 °C

Diaphragm Fixed ground-joint

Installation length 135 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Technical specifications Solvotrode easyClean

Shaft material Glass
pH range 0...14
Temperature range 0...70 °C

Diaphragm Flexible ground-joint

Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm

Drinking water analysis – Does it matter at which stirring rate titration is performed?

When stirring in ion-deficient solutions, streaming potentials occur at pH electrodes with ceramic pin diaphragms which falsify measuring values. In the case of a SET titration, e.g. to a defined pH value, a considerable error can be produced if an incorrect value is measured at the start or at the endpoint of the titration. See page 118 to find out why you can forget about this problem when using the Aquatrode Plus.



Ordering information

Aquatrode Plus without temperature sensor, without cable 6.0253.100 6.0257.600 Aquatrode Plus with Pt1000, without cable, plug-in head U iAquatrode Plus with Pt1000 6.0277.3001 dAquatrode Plus with Pt1000 $6.00202.300^2$ Solvotrode easyClean, fixed cable (1.2 m) plug F 6.0229.010 Solvotrode easyClean, fixed cable (2 m) plug F 6.0229.020 iSolvotrode $6.0279.300^{1}$ dSolvotrode $6.00203.300^2$

¹ Further information about iTrodes can be found on pg. 12.

² Further information about dTrodes can be found on pg. 10.

Special electrodes for pH measurement/pH titration

Biotrode - pH measurement in small volumes

- Very low immersion depth and very small diameter of the electrode tip (3 mm), exceptionally suited to small measuring vessels
- For protein-containing samples and solutions with organic components
- Very low electrolyte outflow (Idrolyte)

26

• LL reference system with long-term stability

Technical specifications Biotrode

Shaft material Glass
pH range 1...11
Temperature range 0...60 °C
Diaphragm Platinum wire
Installation length 113 mm
Shaft diameter 12 mm
Shaft diameter bottom 3 mm
Minimum immersion depth 7 mm

Electrode plug-in head Metrohm plug-in head G

Spearhead electrode – pH measurement in semi-solid samples

- Robust electrode tip for measurements in semi-solid samples such as cheese, meat, fruits, etc.
- Maintenance-free reference electrolyte (gel)
- Easy-to-clean diaphragm
- LL reference system with long-term stability
- With or without integrated temperature sensor
- Lifetime indicator

Technical specifications Spearhead electrode

Shaft material Glass 1...11 pH range Temperature range 0...60 °C Temperature sensor none / Pt1000 Diaphragm Twin pore Installation length 98 mm Shaft diameter 12 mm Shaft diameter bottom 6 mm Minimum immersion depth 10 mm

Electrode plug-in head Metrohm plug-in head G

Flat-membrane electrode – pH measurement on surfaces and in small sample volumes

- For pH measurement on surfaces such as paper, textiles, leather or soil samples (aqueous suspensions)
- Measurement/titration in small sample volumes
- Completely made of glass with extremely finegrounded surface
- LL reference system with long-term stability

Technical specifications Flat-membrane electrode

Shaft material Glass
pH range 0...13
Temperature range 0...80 °C
Diaphragm Fixed ground-joint
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 1 mm
Electrode plug-in head G

Porotrode – pH measurement in protein containing samples

- For pH measurement in very contaminated, proteincontaining or viscous samples
- Low-maintenance capillary diaphragm
- Polymer electrolyte Porolyte for uniform electrolyte outflow
- LL reference system with long-term stability

Technical specifications

Porotrode

Shaft material Glass
pH range 0...14
Temperature range 0...80 °C
Diaphragm Ceramic capillaries
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm
Electrode plug-in head Metrohm plug-in head G



Ordering information	
Biotrode, without cable	6.0224.100
Spearhead electrode without temperature sensor, without cable	6.0226.100
Spearhead electrode with Pt1000, without cable	6.00226.600
Flat-membrane electrode, without cable	6.0256.100
Porotrode, without cable	6.0235.200

Special electrodes for pH measurement/pH titration

Microelectrode – routine use with sample changers and small vials

- For simple acid/base titrations in aqueous solutions
- Available in various lengths (113/168 mm)
- LL reference system with long-term stability

Technical specifications Microelectrode

Shaft material

pH range 0...14

Temperature range 0...80 °C
Diaphragm Ceramic pin

Installation length 113/168 mm Shaft diameter 12 mm

Shaft diameter bottom 6.4 mm Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Glass

Viscotrode – universal application in viscous media

- For viscous protein- or sulfides- containing media
- Flexible ground-joint diaphragm, particularly easy to clean
- LL reference system with long-term stability

Technical specifications

Viscotrode

Shaft material Glass pH range 0...14

Temperature range 0...80 °C

Diaphragm Flexible ground-joint

Installation length 113 mm
Shaft diameter 12 mm
Minimum immersion depth 30 mm

Electrode plug-in head Metrohm plug-in head G

Syntrode with Pt1000 – use in synthesis and in bioreactors

- Low-maintenance thanks to storage vessel for reference electrolytes
- Fixed ground-joint diaphragm insensitive to contamination
- High temperature resistance
- Available in various lengths (288/438 mm)
- LL reference system with long-term stability

Technical specifications

Syntrode

Shaft material Glass
pH range 0...14
Temperature range 0...100 °C
Temperature sensor Pt1000

Diaphragm Fixed ground-joint
Installation length 288/438 mm
Shaft diameter 12 mm
Minimum immersion depth 25 mm



Ordering informationMicroelectrode, length 113 mm, without cable6.0234.100Microelectrode, length 168 mm, without cable6.0234.110Viscotrode, without cable6.0239.100Syntrode with Pt1000, length 288 mm, without cable, plug-in head U6.0248.600Syntrode with Pt1000, length 438 mm, without cable, plug-in head U6.00249.600

Electrodes for pH titration in HF-containing solutions

30 Solitrode HF

- pH-glass sensor with good resistance in HF-containing solutions
- Unbreakable plastic shaft
- Fast response time

Technical Specifications

Solitrode HF

Shaft material 1...12 pH range 0...40 °C Temperature range Reference system LL-System c(KCI) = 3 mol/LReference electrolyte Diaphragm Ceramic pin Installation length 113 mm Shaft diameter 12 mm Minimum immersion depth 15 mm

Electrode plug-in head Metrohm plug-in head G


Combined Sb-electrode

- For pH titration in very hygroscopic matrices or in matrices containing hydrofluoric acid
- Unbreakable plastic shaft

Technical Specifications Combined Sb-electrode

Shaft material 2...11 pH range 0...70 °C Temperature range Reference system LL-System Reference electrolyte c(KCI) = 3 mol/LDiaphragm Ceramic pin Installation length 113 mm Shaft diameter 12 mm Minimum immersion depth 10 mm Electrode plug-in head Metrohm plug-in head G

The recommended fields of application of the electrodes are illustrated in the following graph:

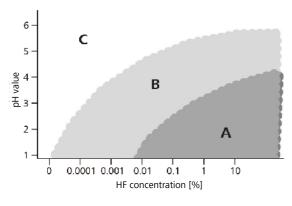

- A Sb-electrode B Solitrode HF
- R 2011ftode HE
- C pH-glass electrode



Ordering information6.0223.100Solitrode HF, without cable6.0223.100Combined Sb-electrode, without cable6.0421.100



Metal electrodes



High-performance metal electrodes for redox and precipitation titration, and water determination according to Karl Fischer.

Separate metal electrodes

34 Separate Pt-wire electrode

- Electrode tip made of Pt-wire (0.8 x 6 mm)
- Auxiliary electrode for differential potentiometry

Technical specifications Separate Pt-wire electrode

Shaft material Glas

Measuring range -2000...2000 mV

Temperature range -20...70 °C

Installation length 125 mm

Shaft diameter 12 mm

Minimum immersion depth 10 mm

Electrode plug-in head Metrohm plug-in head B

Double Pt-sheet electrode

- For bivoltammetric titrations
- Electrode with two polarized Pt sheets (0.15 x 8 x 8 mm)

Technical specifications Double Pt-sheet electrode

Shaft material Gla

Measuring range -2000...2000 mV

Temperature range -20...70 °C

Installation length 101 mm

Shaft diameter 12 mm

Minimum immersion depth 10 mm

Electrode plug-in head Metrohm plug-in head G

Separate Ag-ring electrode

- For precipitation titrations of halides, sulfides, hydrogen sulfide, mercaptans and cyanides
- Available with or without Ag₂S coating (specify when ordering)

Double Au-ring electrode

- For bivoltammetric titrations
- Well suited for Vitamin C determination with DPIP (2,6-Dichlorophenolindophenol)
- Minimum immersion depth 20 mm

Technical specifications

Separate metal-ring electrodesShaft material Glass

Measuring range -2000...2000 mV
Temperature range -20...80 °C
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 7 mm

Electrode plug-in head Metrohm plug-in head G

Separate metal-rod electrodes

 Consisting of separate electrode shaft made of PP and exchangeable metal-rod (76 mm x 2 mm) made of platinum, silver, gold, tungsten or glassy carbon

Degree of purity

Pt 99.90% Ag 99.99% Gold 99.99% W 99.95%

Technical specifications Separate metal-rod electrodes

Total length 162 mm
Installation length 140 mm
Shaft diameter 12 mm
Shaft diameter bottom 8 mm

Electrode plug-in head Metrohm plug-in head B



Ordering information Separate Pt-wire electrode, without cable plug-in head B 6.0301.100 Double Pt-sheet electrode, without cable 6.0309.100 Separate Ag-ring electrode, without cable 6.00350.100 Double Au-ring electrode, without cable 6.00353.100 Shaft for separate metal-rod electrode, without cable, plug-in head B 6.1241.040 Electrode rod Pt 6.1248.000 Electrode rod Ag 6.1248.010 Electrode rod Au 6.1248.030 Electrode rod glassy carbon 6.1248.040 Electrode rod W 6.1248.050

Titrodes – the maintenance-free metal electrodes

36 Pt-Titrode / Pt-Micro Titrode

- For redox titrations without alteration of the pH value
- For bromatometry, iodometry and cerimetry in compliance with Pharm. Europe & USP
- Maintenance-free reference system (pH glass membrane)

Ag-Titrode / Ag-Micro Titrode

- For precipitation titrations without alteration of the pH value
- For precipitation titrations of halides, sulfides, hydrogen sulfide, mercaptans and cyanides
- For titrations in compliance with Pharm. Europe & USP
- Available with or without Ag₂S, AgCl or AgBr coating (specify when ordering)
- Maintenance-free reference system (pH glass membrane)

Au-Micro Titrode

- For ferrometry (determination of the chemical oxygen demand, COD)
- For the determination of penicillin and ampicillin
- For titrations with Hg(NO₃)₂
- For redox titrations in the presence of chromium or iron
- Maintenance-free reference system (pH glass membrane)
- Available in two length (178 mm / 308 mm)

Technical specifications

Titrodes

Shaft material Glass

Measuring range -2000...2000 mV

pH range 0...14
Temperature range 0...80 °C

Reference system pH glass electrode

Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Micro Titrodes

Reference system

Shaft material Glass

Measuring range -2000...2000 mV

pH range 0...14
Temperature range 0...80 °C

Installation length 178 mm

Shaft diameter 12 mm Shaft diameter bottom 6.4 mm

Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

pH glass electrode

Ag-Titrodes: available with or without coating

Depending on the application (see application lists), the use of an Ag-Titrode with or without Ag_2S , AgBr or AgCl coating is recommended.



Ordering information

Ag-Titrode, without cable	6.00430.100
Ag-Titrode, with AgBr coating, without cable	6.00430.100Br
Ag-Titrode, with AgCl coating, without cable	6.00430.100Cl
Ag-Titrode, with Ag₂S coating, without cable	6.00430.100\$
iAg-Titrode, plug-in head K	6.00470.300
iAg-Titrode with Ag₂S coating, plug-in head K	6.00470.300\$
dAg-Titrode, plug-in head Q	6.00404.300
dAg-Titrode with Ag₂S coating, plug-in head Q	6.00404.300\$
Pt-Titrode, without cable	6.0431.100
iPt-Titrode, plug-in head K	6.0471.300 ¹
dPt-Titrode, plug-in head Q	6.00401.3002
Micro Ag-Titrode, without cable	6.0433.110
Micro Pt-Titrode, without cable	6.0434.110
Micro Au-Titrode, length 178 mm, without cable	6.0435.110
Micro Au-Titrode, length 308 mm, without cable	6.00435.120

¹ Further information about iTrodes can be found on pg. 12.

² Further information about dTrodes can be found on pg. 10.

Combined metal electrodes

38 Combined Ag-ring electrode

- For precipitation titrations of halides, sulfides, hydrogen sulfide, mercaptans and cyanides with alteration of the pH value
- Available with or without AgBr, AgCl or Ag₂S coating (specify when ordering)

Combined Pt-ring electrode

• For redox titrations with alteration of the pH value

Combined Au-ring electrode

- For ferrometry (determination of the chemical oxygen demand COD)
- For determination of penicillin and ampicillin

Technical specifications Combined Ag-ring electrode

Shaft material Glass

Measuring range -2000...2000 mVTemperature range 0...70 °CReference system LL systemReference electrolyte $c(KNO_3) = 1 \text{ mol/L}$ Diaphragm Fixed ground-joint

Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 15 mm

Electrode plug-in head Metrohm plug-in head G

Combined Pt-ring electrode / Au-ring electrode

Shaft material Glass -2000...2000 mV Measuring range Temperature range -5...80 °C Reference system LL system c(KCI) = 3 mol/LReference electrolyte Diaphragm Ceramic pin Installation length 113 mm Shaft diameter 12 mm Minimum immersion depth 15 mm

Electrode plug-in head Metrohm plug-in head G



Ordering information

Combined Ag-ring electrode, without cable
iAg-ring electrode, combined
6.00450.300
dAg-ring electrode, combined
6.00402.300²
Combined Pt-ring electrode, without cable
iPt-ring electrode, combined
6.0451.100
dPt-ring electrode, combinded
6.00403.300²
Combined Au-ring electrode, without cable
6.00403.300²
Combined Au-ring electrode, without cable
6.00452.100

¹ Further information about iTrodes can be found on pg. 12.

² Further information about dTrodes can be found on pg. 10.

Electrodes for Karl Fischer titration

Double Pt-wire electrodes

40

Indicator electrode for volumetric KF determination

- For «Ipol» and «Upol» mode titrations
- Thanks to its two fused platinum rings, is more robust and easier to clean than the conventional platinum pin electrode (6.0338.100)

Technical specifications

Measuring range -2000...2000 mV
Temperature range -20...70 °C
Installation length 96 mm
Shaft diameter 8 mm
Shaft diameter bottom 3 mm
Minimum immersion depth 7 mm

Electrode plug-in head Metrohm plug-in head G

Indicator electrode for coulometric KF determination

• With standard ground-joint 14/15

Technical specifications

Measuring range -2000...2000 mV

Temperature range -20...70 °C

Installation length 101 mm

Shaft diameter 12 mm

Shaft diameter bottom 8.75 mm

Minimum immersion depth 10 mm

Electrode plug-in head Metrohm plug-in head G

Indicator electrode for KF sample changers

• Fixed cable (length 2 m with plug F)

Technical specifications

Measuring range -2000...2000 mV
Temperature range -20...70 °C
Installation length 103 mm
Shaft diameter 5.3 mm
Minimum immersion depth 10 mm

Generator electrodes

Generator electrode with diaphragm

- Standard ground-joint 29/22 and Metrohm plug-in head G
- Requires cable 6.2104.120 for connection with KF Coulometers

Generator electrode without diaphragm

- Standard ground-joint 29/22 and Metrohm plug-in head G
- Requires cable 6.2104.120 for connection with KF Coulometers

Technical specifications

Temperature range -20...70 °C Installation length 108 mm Shaft diameter 24 mm Minimum immersion depth 15 mm

Technical specifications

Temperature range -20...70 °C Installation length 108 mm Shaft diameter 24 mm Minimum immersion depth 15 mm



Ordering informationIndicator electrode for KF volumetry (double Pt ring), without cable6.00346.100Indicator electrode for KF volumetry (double Pt wire), without cable6.0338.100Indicator electrode for KF coulometry, without cable6.0341.100Indicator electrode for KF Sample Changers, fixed cable (2 m), plug F6.0340.000Generator electrode with diaphragm, without cable6.0344.100Generator electrode without diaphragm, without cable6.0345.100





Electrodes for ion and surfactant analysis

Ion-selective electrodes

44 Crystal-membrane electrodes

- Robust construction
- Can also be used for brief periods in organic solvents
- Simple cleaning and renewal of electrode surface with polishing set

Polymer-membrane electrodes for K^+ , NO_3^- , Na^+ and Ca^{2+}

- Robust construction
- High selectivity due to ionophores immobilized in the membrane
- Short preparation time after conditioning in a standard solution
- For aqueous solutions

Ammonia-selective gas-membrane electrode

- Robust construction
- Short preparation time after conditioning in a standard solution
- The gas-permeable membrane ensures high selectivity and prevents interferences by the measuring solution

Choose the right sensor for your application:

NH₃-ISE (Low) 6.0506.100 / 6.1255.000

- Clean samples (e.g., drinking water, boiler feed water, etc.)
- Faster response time at the detection limit
- Lower detection limits
- Individually tested and certified
- Complete modules for simple exchange

NH₃-ISE (High) 6.0506.150 / 6.1255.050

- Waste water samples
- For long-term measurements, monitoring
- Faster response times after measuring high ammonium concentrations
- Better signal stability at high ammonium concentrations
- Affordable exchange of contaminated membranes (e.g., oil residues)





Ion	Article no.	Membrane	Min.	Installation-	Shaft-	Tempera-	Measure-	pH range
		material	immersion	length	diameter	ture range	ment range	
			depth (mm)	(mm)	(mm)	(°C)	(mol/L)	
Ag+	6.0502.180	Crystal	1	125	12	080	10 ⁻⁷ 1	28
Br	6.0502.100	Crystal	1	125	12	050	10 ⁻⁶ 1	014
Ca ²⁺	6.0510.100	Polymer ²	10	113	12	040	5*10 ⁻⁷ 1	212
Cl-	6.0502.120	Crystal	1	125	12	050	10 ⁻⁵ 1	014
CN-	6.0502.130	Crystal	1	125	12	080	8*10 ⁻⁶ 10 ⁻²	1014
Cu ²⁺	6.0502.140	Crystal	1	125	12	080	10 ⁻⁸ 10 ⁻¹	212
F-	6.00500.300	Crystal ⁴	20	125	12	050	10 ⁻⁶ sat.	57
F-	6.00500.600	Crystal ⁵	20	125	12	050	10 ^{−6} sat.	57
I -	6.0502.160	Crystal	1	125	12	050	5*10-81	014
K+	6.0510.110	Polymer ³	10	113	12	040	10 ⁻⁷ 1	2.511
Na+	6.0508.100	Polymer	1	125	12	040	5*10 ⁻⁶ 1	212
NH_4^+	6.0506.100	Gas membrane	5	125	12	050	5*10-610-2	11
NH_4^+	6.0506.150	Gas membrane	5	125	12	050	10-41	11
NO_3^-	6.00510.120	Polymer ¹	10	113	12	040	10-61	2.511
Pb^{2+}	6.0502.170	Crystal	1	125	12	080	10-610-1	47
S ²⁻	6.0502.180	Crystal	1	125	12	080	10 ⁻⁷ 1	212

 $^{^{1}}$ Combined sensor. Ceramic pin diaphragm; electrolyte c(KCl) = 3 mol/L.

 $^{^{2}}$ Combined sensor. Ceramic pin diaphragm; electrolyte c(NH₄NO₃) = 1 mol/L.

 $^{^{3}}$ Combined sensor. Ceramic pin diaphragm; electrolyte c(CH $_{3}$ COOLi) = 1 mol/L.

 $^{^4}$ Combined sensor. Ground-joint diaphragm; electrolyte c(KCl) = 3 mol/L.

 $^{^{5}}$ Combined, digital sensor. Ground-joint diaphragm; electrolyte c(KCl) = 3 mol/L.

LL ISE reference 6.0750.100

Double-junction Ag/AgCl reference electrode with fixed ground-joint diaphragm and optimized length for sample changer applications. Standard bridge electrolyte: c(KCl) = 3 mol/L.



LL ISE Reference

A stable, reproducible reference potential is very important at low ion concentrations, low ionic strengths and especially with repeated determinations using sample changer systems. For this reason Metrohm recommends reference electrodes with a fixed ground-joint diaphragm for working with ion-selective electrodes. In addition to a constant electrolyte outflow of approx. 5...30 µL/h, these electrodes are also considerably less influenced by either the ionic strength of the sample solution or the stirring speed than other types of reference electrodes.



47

Spoilt for choice!

Accessories for ion-selective electrodes

To what must I pay particular attention in an ion determination? Precision? Time needed? Costs?

Which method is the most suitable for my application? Titration? Direct measurement? Standard addition?

ISA? TISAB? When is their use advisable? Which solution do I need for my application? You will find the answers to these questions along with many other useful tips for ion determination with ion-selective electrodes from Metrohm in the theoretical part in section 1.3.3. «Ion-selective electrodes.»

49

Surfactant electrodes for two-phase titration

Surfactrode Refill

- Refillable surfactant electrode for the titration of ionic surfactants in non-aqueous solvents
- Renewable electrode surface, therefore practically unlimited working life
- Resistant to virtually all conventional solvents used in surfactant analysis (not to chloroform)
- Particularly suitable for titration of detergents and soap

Technical specifications

Shaft material PEEK
pH range 0...13
Temperature range 0...40 °C
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 1 mm

Electrode plug-in head Metrohm plug-in head G

Surfactrode Resistant

- Durable surfactant electrode for the two-phase titration of anionic and cationic surfactants in non-aqueous solvents
- Easy to clean and low-maintenance, therefore particularly suitable for use in sample changer systems
- Resistant to chloroform and all solvents used in surfactant analysis
- Particularly suitable for samples containing oil such as drilling and cutting oils or cooling lubricants

Polymer-membrane surfactant electrodes for environmentally-friendly surfactant titration

Cationic Surfactant electrode

- For the titration of cationic and anionic surfactants in aqueous matrices
- Optimized for cationic surfactants
- Excellent response due to ionophores immobilized in the membrane
- Long working life with normal use

Ionic Surfactant electrode

- For the titration of anionic and cationic surfactants in aqueous matrices
- Excellent response due to ionophores immobilized in the membrane
- Long working life with normal use

NIO Surfactant electrode

- For the titration of non-ionic surfactants in aqueous matrices
- For the titration of surfactants based on polyoxyethylene adducts
- For the titration of pharmaceutical ingredients
- Long working life with normal use

Technical specifications

Shaft material POM
pH range 0...10
Temperature range 10...50 °C
Installation length 108 mm
Shaft diameter 12 mm
Minimum immersion depth 5 mm

Electrode plug-in head Metrohm plug-in head G

Technical specifications

Cationic Surfactant electrode Ionic Surfactant electrode NIO Surfactant electrode

Shaft material PVC
pH range 0...12
Temperature range 0...40 °C
Installation length 125 mm
Shaft diameter 12 mm
Shaft diameter bottom 2.5 mm
Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G



Ordering informationSurfactrode Refill, without cable6.0507.140Surfactrode Resistant, without cable6.0507.130Cationic Surfactant electrode, without cable6.0507.150Ionic Surfactant electrode, without cable6.0507.120NIO Surfactant electrode, without cable6.0507.010

Accessories for surfactant electrodes

Refill set for Surfactrode Refill

Paste for Surfactrode Refill, 3.5 g 6.2319.000 Filling tool 6.2826.010

Reagents for surfactant titration

TEGO trant A100, titrant for anionic surfactants

18 g 6.2317.030

TEGO add, additive for two-phase titration

250 mL 6.2317.130







Reference electrodes – our best references

Reference electrodes

Double-junction reference electrodes

54

Ag/AgCl reference electrode with Metrohm plug-in head B

- Easy to change reference and bridge electrolytes
- Variable electrolyte outflow at the flexible groundjoint diaphragm
- Available with 125 mm or 162 mm shaft length
- With standard ground-joint 14/15

Ag/AgCl reference electrode with Metrohm plug-in head G

- For differential potentiometry with Metrohm titrators
- Easy to change reference and bridge electrolytes
- Variable electrolyte outflow at the flexible groundjoint diaphragm
- Available with 125 mm or 162 mm shaft length
- With standard ground-joint 14/15

LL ISE Reference

- Double-junction Ag/AgCl reference electrode
- High signal stability thanks to constant, reproducible electrolyte outflow, therefore particularly suitable for sample changer applications
- Fixed ground-joint diaphragm insensitive to contamination
- Minimum immersion depth of 1 mm

Technical specifications

Shaft material Glass
Temperature range 0...80 °C

Diaphragm Flexible ground-joint

Installation length 100/138 mm
Shaft diameter 12 mm
Shaft diameter bottom 12/8 mm
Minimum immersion depth 10 mm

Reference system Ag wire + AgCl

Electrode plug-in head Metrohm plug-in head B

Technical specifications

Shaft material Glass

Temperature range 0...80 °C

Diaphragm Flexible ground-joint Installation length 100/138 mm

Installation length 100/138
Shaft diameter 12 mm
Minimum immersion depth 10 mm

Reference system Ag wire + AgCl

Electrode plug-in head Metrohm plug-in head G

Technical specifications

Shaft material Glass
Temperature range 0...80 °C

Diaphragm Fixed ground-joint

Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 1 mm
Reference system LL system

Electrode plug-in head Metrohm plug-in head B



Reference electrodes

Modular reference system

- Consisting of Ag/AgCl reference system with standard ground-joint 14/15 and exchangeable electrolyte vessel
- Electrolyte vessels without storage vessels, with ceramic diaphragm, various diaphragm diameters

Technical specifications Ag/AgCl reference system

Shaft material Glass
Temperature range 0...80 °C
Shaft length 50 mm

Length to the upper edge

Standard ground-joint 43 mm
Shaft diameter top 12 mm
Shaft diameter bottom 8 mm
Minimum immersion depth 20 mm
Reference system Ag wire

Reference system Ag wire + AgCl

Electrode plug-in head Metrohm plug-in head B

Electrolyte vessels without storage container

Length to the upper edge


Standard ground-joint 101 mm
Shaft material PTFE/glass
Diaphragm Ceramic pin
Shaft diameter 3/5.5 mm

Diaphragm diameter 3 mm (PTFE)/1 mm (glass)



57

Ordering informationAg/AgCl reference system, without cable, plug-in head B6.0724.140Electrolyte vessel made of PTFE without storage vessel, diaphragm diameter 3 mm6.1240.000Electrolyte vessel made of glass without storage vessel, diaphragm diameter 1 mm6.1240.020

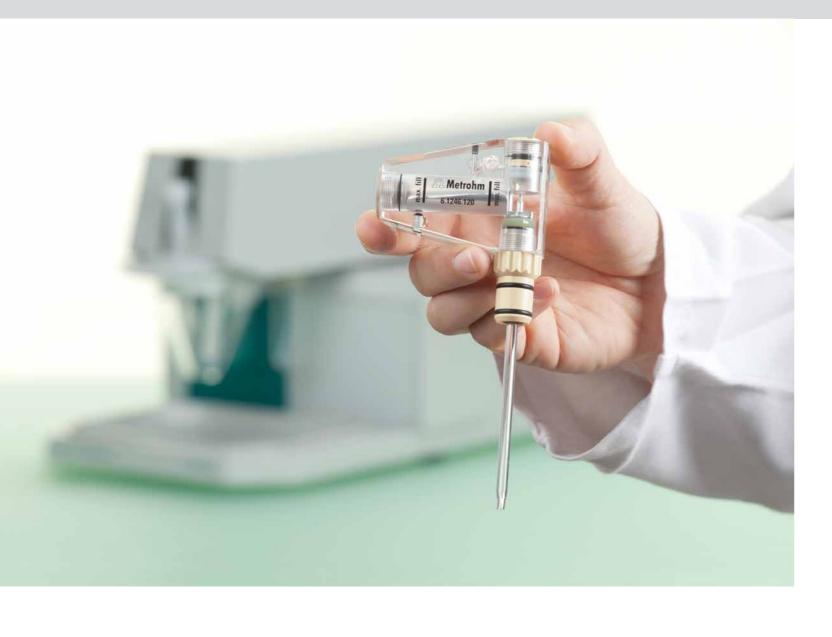

Electrodes for electrochemistry



Electrodes for voltammetric trace analysis, cyclic voltammetric stripping (CVS), cyclic voltammetry and other electrochemical techniques in Metrohm VA instruments or in combination with potentiostats from Metrohm Autolab or Metrohm DropSens.

The electrodes described in the following can be used in various Metrohm voltammetry instruments: 663 VA Stand, 747 VA Stand, 757 VA Computrace, 797 VA Computrace, 884 Professional VA, and 894 Professional CVS.

Multi-Mode Electrode pro

- Universally applicable working electrode for polarography and voltammetry
- Determination of heavy metal ions, organic substances, anions

Glass capillaries for Multi-Mode Electrode pro Non-silanized glass capillaries

- Standard capillary for polarography and stripping voltammetry in alkaline solutions
- For universal use with all pH values in aqueous and non-aqueous solutions

Silanized glass capillaries

- Silanized capillaries for stripping voltammetry in acidic to mildly alkaline solutions
- Long lifetime

Electrodes for mercury-free trace analysis scTRACE Gold

- Combined gold microwire electrode
- For arsenic, mercury, copper, lead and other metals
- For laboratory instruments and the 946 Portable VA Analyzer with scTRACE Gold measuring head

Screen-printed electrodes (SPE)

• Cost effective combined screen-printed electrodes

• Determination of various heavy metals

• For laboratory instruments and the 946 Portable VA Analyzer

Bismuth drop electrode

- Separate working electrode
- Bi drop electrode for Cd, Pb, Ni, Co, Fe
- For laboratory instruments and online analyzers

6.1241.080 +

6.1258.000





Multi-Mode Electrode pro 6.1246.120
Glass capillaries, not silanized, 10 units 6.1226.030
Glass capillaries, silanized, 10 units 6.1226.050



Ordering information

scTRACE Gold (set of 4) 6.1258.000
Holder for scTRACE Gold 6.1241.080
Carbon screen-printed electrode DRP-11L
SPE electrode holder 6.1241.090
Bi drop electrode 6.0346.000

RDE – rotating disk electrodes

An RDE consists of a driving axle and an exchangeable electrode tip.

Ordering information	
Drive for rotating disk electrode (RDE):	
for VA measuring stands	6.1204.210
with mercury contact for VA	
measuring stands	6.1204.220
for Professional VA/CVS systems	6.1204.510
with mercury contact for Professional	
VA/CVS systems	6.1204.520



Electrode tips for the RDE

Order number	Electrode tip	Applications	Determination range
6.1204.600	Glassy carbon Ø 2 mm Shaft made of glass	Determination of heavy metals with anodic stripping voltammetry (mercury film technological kinetic and thermodynamic studies in electrochemistry	
6.1204.130	Silver	Determination of halides and pseudohalid	es ppb to ppt
6.1204.140	Gold	Determination of mercury and other meta with anodic stripping voltammetry	l ions ppb to ppt
6.1204.170	Platinum	Determination of organic additives in elec-	troplating ppm to ppt
	Ø 3 mm	baths with cyclic voltammetric stripping	
	polished	technique (CVS), kinetic and thermodynan	nic
		studies in electrochemistry	
6.1204.190	Platinum Ø 1 mm polished Shaft made of glass	Determination of organic additives in election baths with cyclic voltammetric stripping technique (CVS), kinetic and thermodynam studies in electrochemistry	
6.1204.610	Platinum	Determination of organic additives in elec-	troplating ppm to ppt
	Ø 2 mm	baths with cyclic voltammetric stripping	7 3 11 11
	polished	technique (CVS), kinetic and thermodynan	nic
	Shaft made of glass	studies in electrochemistry	

Reference and auxiliary electrodes for trace analysis and CVS

62

	and CVS			
	Reference electrodes Reference electrode made of plastic with ceramic	Ordering information		
	 diaphragm, filled Universal double-junction reference electrode for voltammetry Inner system filled with c(KCl) = 3 mol/L Aqueous solutions, trace and ultra trace ranges Low carry-over effects, low blank values 	Ag/AgCl reference electrode Electrolyte vessel for reference electrode	6.0728.120 6.1245.010	
	Reference electrode made of plastic with ceramic diaphragm, dry • Double-junction reference electrode. Reference system dry • Studies in organic solvents with any electrolyte solution • Low carry-over effects, low blank values	Ag/AgCl reference electrode (dry) Electrolyte vessel for reference electrode	6.0728.110 6.1245.010	
	 LL reference electrode made of plastic with ceramic diaphragm, filled Double-junction reference electrode for the analysis of electroplating baths with cyclic voltammetric stripping (CVS) Inner system filled with c(KCl) = 3 mol/L Very stable reference potential 	LL-Ag/AgCl reference electrode Electrolyte vessel for reference electrode	6.0728.130 6.1245.010	
	 LL reference electrode made of glass, filled Double-junction reference electrode for the studies of electroplating baths with cyclic voltammetric stripping (CVS) Very stable reference potential Maintenance-free 	LL-Ag/AgCl (gel) reference electrode	6.0730.100	
	Reference electrode made of glass with ground-joint diaphragm Double-junction reference electrode with dry reference system For aqueous as well as non-aqueous solutions Simple electrolyte replacement, low outflow rate	Ag/AgCl reference electrode glass Electrolyte vessel for reference electrode	6.0728.100 6.1245.000	
• Uni • For rota	Auxiliary electrodes Platinum auxiliary electrode Universal auxiliary electrode for voltammetry For all applications with the MME as well as with the rotating platinum disc electrode Robust, easy maintenance	Platinum auxiliary electrode	6.0343.100	
	Glassy carbon auxiliary electrode	Auxiliary electrode holder	6.1241.120	

Glassy carbon rod

• For all applications with rotating disc electrodes as

• Inert surface, no contamination of the measuring

well as with the MME

cell by platinum

6.1247.000



Electrodes for Metrohm Autolab RDE and RRDE

64

Rotating disk electrode and rotating ring disk electrode measurements are powerful experimental methods in which the electrode rotates while immersed in the solution. This creates a well-defined convective drag towards the surface of the electrode as long as a laminar flow is present. Under these conditions, mass transport occurs through convection and diffusion, with both components directly influenced by the angular frequency of the electrode.

Measurements under hydrodynamic conditions provide great benefits with respect to static measurements. Because of the increase in mass transport, the relative contributions from the electron transfer kinetics increase. Electrochemical experiments therefore can provide access to kinetic or mechanistic information.

RDE/RRDE electrode tips

The rotating disk electrode (RDE) consists of an electrode disk inserted in an electrode shaft. The rotating ring disc electrode (RRDE) is based on a similar design with an additional ring electrode located around the disk. The gap between the ring and the disk is typically very small.

Metrohm Autolab provides a range of electrodes for RDE measurements with a wide choice of materials. An empty electrode holder suitable for RDE measurement is also available. RRDE electrodes with a Pt ring are available with a choice of disk material.

Counter electrode

• Pt-sheet counter electrode (8 x 7 mm)



Technical specifications RDE tips

Shaft material PEEK
Length 52.5 mm
Shaft diameter 10 mm
Disk diameter 3 or 5 mm

Disk material Pt, Ag, Au, GC, Cu, Zn,

stainless steel, empty

Connector M4 thread

RRDE tips

Shaft material PEEK
Length 52.5 mm
Shaft diameter 10 mm
Disk diameter 5 mm
Disk material Pt, Au, GC
Ring material Pt
Connector M4 thread

Technical specifications Pt sheet electrode

Shaft material Glass
Temperature range -20...70 °C
Installation length 101 mm
Shaft diameter 12 mm
Minimum immersion depth 10 mm
Electrode plug-in head B (4 mm)



Order number	Electrode tip	Applications
6.09395.014	Glassy carbon disk (3 mm)	Interfacial electrochemistry
6.09395.024	Platinum disk (3 mm)	Interfacial electrochemistry
6.09395.034	Gold disk (3 mm)	Interfacial electrochemistry
6.09395.044	Silver disk (3 mm)	Interfacial electrochemistry
RRDE.PTPT.S	Platinum disk (5 mm) and platinum ring	Electrocatalysis
RRDE.AUPT.S	Gold disk (5 mm) and platinum ring	Electrocatalysis
RRDE.GCPT.S	Glassy carbon disk (5 mm) and platinum ring	Electrocatalysis
3500003900	Glassy carbon disk (5 mm)	Interfacial electrochemistry
RDE.PT50.S	Platinum disk (5 mm)	Interfacial electrochemistry
RDE.AU50.S	Gold disk (5 mm)	Interfacial electrochemistry
RDE.AG50.S	Silver disk (5 mm)	Interfacial electrochemistry
RDE.CU50.S	Copper disk (5 mm)	Interfacial electrochemistry
RDE.ZN50.S	Zinc disk (5 mm)	Interfacial electrochemistry
RDE.STEEL.S	Stainless steel disk (5 mm)	Interfacial electrochemistry
RDE.BLANK.S	Empty disk (5 mm)	_

Microelectrodes

66

Microelectrodes or ultramicroelectrode are electrodes whose dimensions are small enough that their electrochemical properties become a function of size. These working electrodes are much smaller than typical metallic electrodes. When the dimensions are in the range of 25 μm or lower, the term ultramicroelectrode is preferred. The most common geometry of these electrodes is a disk.

Because of their very small size, these electrodes exhibit particularities with respect to their macroscopic counterparts:

- Mass transport of electroactive species changes from a linear to a spherical diffusion (in case of disk electrodes). This results in an enhancement of the mass transport to and from the electrode surface.
- The current measured at such electrodes becomes smaller but is not proportional to the surface area.
- The current density increases.

Moreover, these electrodes are less affected by charging or capacitive currents with the RC constant decreasing with the radius of the disk. This makes them particularly suitable for studying electrochemical reactions with rapid kinetics. These electrodes are also less affected by ohmic drop, which in turn facilitates measurements in low conductivity media.

Metrohm Autolab offers a wide range of glass-sealed disk microelectrodes and ultramicroelectrodes of different sizes and materials. Special electrodes can be manufactured upon request.

Technical specifications Microelectrodes

Shaft material Glass
Temperature range 0...40 °C
Length 51.5 mm
Shaft diameter 5 mm
Disk diameter 8...1000 µm
Disk material Pt, Ag, Au, GC, Ir, Pd
Connector B (2 mm)





Order number	Electrode tip	Applications
PT.10	Platinum disk (10 µm)	Interfacial electrochemistry
PT.20	Platinum disk (20 µm)	Interfacial electrochemistry
PT.25	Platinum disk (25 µm)	Interfacial electrochemistry
PT.50	Platinum disk (50 µm)	Interfacial electrochemistry
PT.100	Platinum disk (100 µm)	Interfacial electrochemistry
PT.200	Platinum disk (200 µm)	Interfacial electrochemistry
PT.300	Platinum disk (300 µm)	Interfacial electrochemistry
PT.500	Platinum disk (500 µm)	Interfacial electrochemistry
AU.10	Gold disk (10 µm)	Interfacial electrochemistry
AU.25	Gold disk (25 µm)	Interfacial electrochemistry
AU.40	Gold disk (40 µm)	Interfacial electrochemistry
AU.50	Gold disk (50 µm)	Interfacial electrochemistry
AU.100	Gold disk (100 µm)	Interfacial electrochemistry
AU.200	Gold disk (200 µm)	Interfacial electrochemistry
AU.300	Gold disk (300 µm)	Interfacial electrochemistry
AU.500	Gold disk (500 µm)	Interfacial electrochemistry
AG.25	Silver disk (25 µm)	Interfacial electrochemistry
AG.30	Silver disk (30 µm)	Interfacial electrochemistry
AG.100	Silver disk (100 µm)	Interfacial electrochemistry
AG.300	Silver disk (300 µm)	Interfacial electrochemistry
IR.75	Iridium disk (75 μm)	Interfacial electrochemistry
PD.25	Palladium disk (25 µm)	Interfacial electrochemistry
PD.100	Palladium disk (100 µm)	Interfacial electrochemistry
PD.300	Palladium disk (300 µm)	Interfacial electrochemistry
PD.500	Palladium disk (500 µm)	Interfacial electrochemistry

Electrodes for the electrochemical quartz crystal microbalance

The electrochemical quartz crystal microbalance (EQCM) is widely used to monitor, simultaneously with the electrochemical signal, the change in frequency which is directly related to the mass changes due to deposition or adsorption of a species to or dissolution of a species from the working electrode.

The principle of operation of the EQCM relies on the piezoelectric properties of the quartz which will oscillate at a given frequency when a sinusoidal external electric field is applied to it. The oscillation frequency of the crystal depends on a number of parameters such as size, the thickness of the crystal, temperature and the oscillating media.

Quartz crystals

68

The crystals provided for the Autolab EQCM are AT-cut crystals (with a small temperature coefficient). They are coated with different metals on top of an adhesion layer (e.g. Ti or Cr oxides). The crystals will also act as working electrodes (WE) being in contact with the electrolyte solution, broadening in this way the application areas of this technique.



Technical specifications

Quartz crystals

Material Quartz
Quartz diameter 13.66 mm
Quartz thickness 250 µm
Electrode diameter 6.80 mm
Electrode material Gold, Platinum
Nominal frequency 6 MHz

Reference electrode

• Single junction reference electrode for EQCM

Technical specifications Reference electrode

Reference system

Shaft material

Shaft length

Shaft diameter

Diaphragm

Temperature range

Connector

LL Ag/AgCl

Glass

55 mm

4 mm

Ceramics

Ceramics

0...40 °C

Counter electrode

• Gold coil counter electrode for EQCM

Technical specifications Counter electrode

Material Gold wire
Wire diameter 1 mm
Length 37.5 mm
Coil diameter 8 mm



Order number	Electrode tip	Applications
EQCM.Au	EQCM 6 MHz Au/TiO2 quartz crystal, polished	Interfacial electrochemistry, electrodeposition
EQCM.CE	EQCM gold counter electrode	Interfacial electrochemistry, electrodeposition
EQCM.REF.EL.S	EQCM Ag/AgCl 3 M KCl reference electrode	Interfacial electrochemistry, electrodeposition

Screen-printed electrodes

Screen-printed electrodes (SPEs) are cost-effective disposable electrodes specially designed to work with microvolumes of sample which is ideal for research purposes and also for teaching electrochemistry.

There is a wide range of SPEs classified below by its working electrode (WE) material and differentiated by the configuration of the electrochemical cell. The reference electrode (RE) material in all of them is silver (unless electrode ref. 11L and C11L which is silver / silver chloride) and, depending on the model, the auxiliary electrode (AUX) material can be carbon, gold or platinum. Also depending on the model, the substrate material can be ceramics, FR4, transparent plastic or white plastic. There are also customized designs available upon request.

Working electrode made of carbon

• Standard type

70

- Work in solution type
- MultiAnalysis configuration dual
- MultiAnalysis configuration 8X electrochemical cells
- MultiAnalysis configuration 96X electrochemical cells
- MultiAnalysis configuration 4WE sharing AUX
- MultiAnalysis configuration 8WE sharing AUX and REF
- Integrated flow-cell

Working electrode made of gold

- Standard type
- Work in solution type
- WE of 1.6 mm
- MultiAnalysis configuration 8X electrochemical cells
- MultiAnalysis configuration 96X electrochemical cells
- Integrated flow-cell

Working electrode made of platinum

- Standard type
- Work in solution type
- MultiAnalysis configuration 96X electrochemical cells
- Integrated flow-cell

Working electrode made of silver

- WE of 1.6 mm
- Standard type

Working electrode made of palladium

Standard type

Transparent working electrode

Materials: ITO, PEDOT, gold, carbon, silver

Standard type

Mediator modified working electrode

Materials: Co-phthalocyanine, meldola's blue, prussian blue, ferrocyanide, manganese (II) phthalocyanine, copper (II) phthalocyanine

- Standard type
- MultiAnalysis configuration dual

Biomodified working electrode

Materials: Streptavidin, extravidin, glucose sensor, lactate sensor, uric acid sensor

- Standard type
- MultiAnalysis configuration dual

Nanomodified working electrode


Materials: Graphene, graphene oxide, reduced graphene oxide, multi-walled carbon nanotubes, single-walled carbon nanotubes, carbon nanofibers, mesoporous carbon, ordered mesoporous carbon, gold nanoparticles, silver nanoparticles, quantum dots, core-shell quantum dots, graphene-gold nanoparticles, carbon nanotubes-gold nanoparticles, carbon nanofibres-gold nanoparticles, gold nanoparticles-streptavidin

- Standard type
- MultiAnalysis configuration dual

Other working electrodes

Materials: Bismuth oxide, nickel oxide, polyaniline, polypyrrole, palladium particles, gold particles, platinum particles, rhodium particles, iridium particles, phenanthroline, poly-L-lysine, potassium ferricyanide, alizarin, anthraquinone, hydrogen ionophore, potassium ionophore, sodium ionophore

- Standard type
- MultiAnalysis configuration dual

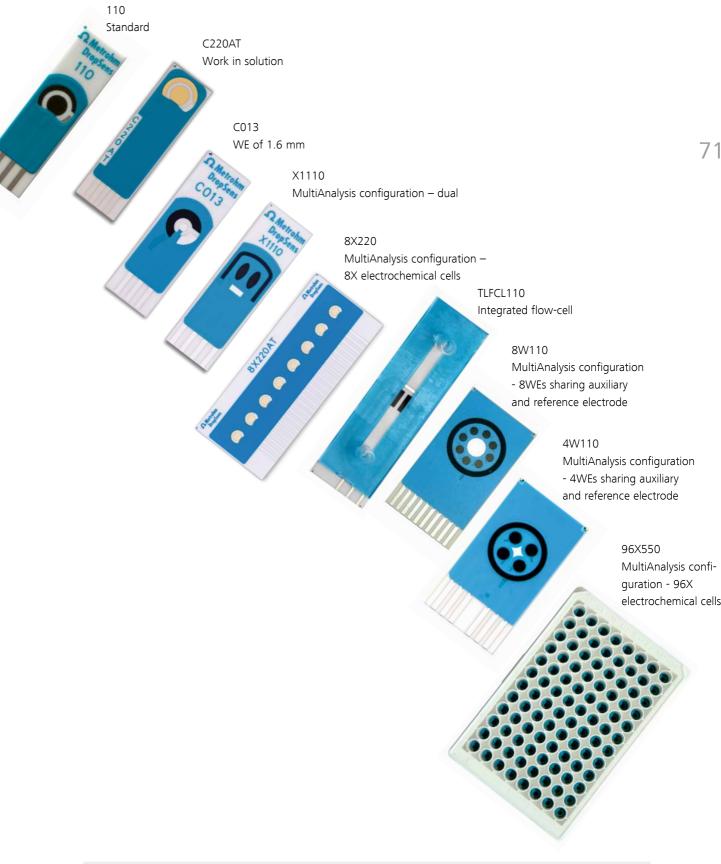

Ordering information

Screen-printed carbon electrode

Screen-printed gold electrode / ink AT / work in solution

Screen-printed silver electrode / d 1.6 mm

Dual screen-printed carbon electrode 8X screen-printed gold electrode

Thin-layer flow-cell integrated screen-printed carbon electrode (Aux.: C; Ref.: Ag)

8 WEs screen-printed carbon electrode (1 AUX.:C; 1 REF.:Ag)

4 WEs screen-printed carbon electrode (1 AUX.:C; 1 REF.:Ag)

96X screen-printed platinum electrode

DRP-110-U75
DRP-C220AT-U75
DRP-C013-U75
DRP-X1110-U75
DRP-8X220AT-U20
DRP-TLFCL110-U10
DRP-8W110-U20
DRP-4W110-U20
DRP-96X550-U4

Interdigitated electrodes/microelectrodes

Interdigitated electrodes and microelectrodes offer several advantages, such as working with low volumes of sample and avoiding tedious polishing of solid electrodes. In addition, the interdigitated configuration typically enhances sensitivity and detection limits.

Depending on the model, substrate materials can be ceramics, transparent plastic, white plastic or glass. AUX and REF electrodes, in models which have them, are made of gold or platinum.

Interdigitated electrodes: gold, platinum

• Squared design

72

- Concentric design
- With auxiliary and reference electrode
- With a platinum heater
- Field effect transistor

Interdigitated electrodes: silver, copper, palladium, ITO

• Squared design

Microelectrodes: gold, platinum

Multi-interdigitated: gold



Ordering information

Interdigitated platinum concentric electrode / 10 microns lines and gaps / glass substrate Interdigitated gold electrode / 100 microns lines and gaps / plastic substrate Interdigitated gold electrode (Aux.:Au; Ref.:Au) / 10 microns lines and gaps / glass substrate Platinum microelectrode array d. 3mm / microholes 10 microns / glass substrate Field-effect transistor gold electrodes / 30 microns lines and gaps Multi-Interdigitated gold electrodes / 40 microns lines and gaps

DRP-G-IDECONPT10-U20 DRP-PW-IDEAU100-U50 DRP-G-IDE222-U20 DRP-G-MEA555-U20 DRP-AUFET30-U50 DRP-PW-4XIDEAU40-U50

Custom-made electrochemical sensors

If you are developing a sensor for a final application, Metrohm DropSens is the experienced partner that will manufacture your customized electrode with a scalable and cost-effective process. Benefit from the multiple possibilities for sensor customization: choose the substrate, place your logo, decide on the geometry and have as many material combinations as needed to suit your electrochemical application.



Key service features:

- Custom production and high-volume manufacturing of electrochemical solutions
- Certified easy-to-use, and best-in-class instruments for your electrochemical research
- Support throughout the entire journey: from the initial idea, to prototyping, and up to the implementation of the final solution.
- Expertise, flexibility, and understanding





Conductivity measuring cells and temperature sensors

Conductivity measuring cells for OMNIS Measuring Module Conductivity

The new conductivity measuring cells are equipped with a fixed cable labelled with a yellow marking and plug N for direct connection to the OMNIS Measuring Module Conductivity.

The measuring cells are based on the already existing sensors for 912 Conductometer or 856 Conductivity Module but have optimized cable length for manual and automated OMNIS systems.

Conductivity measuring cell c = 0.1 cm⁻¹

• With Pt1000 temperature sensor

76

• Available in 0.65 m and 2.0 m cable length

Conductivity measuring cell c = 0.5 cm⁻¹

- With Pt1000 temperature sensor
- Available in 0.65 m and 2.0 m cable length

Conductivity measuring cell $c = 0.7 \text{ cm}^{-1}$

- With Pt1000 temperature sensor
- Available in 0.65 m cable length

Conductivity measuring cell c = 1.6 cm⁻¹

- With Pt1000 temperature sensor
- Available in 0.65 m and 2.0 m cable length

Technical specifications

Shaft material
Ideal measuring range
Temperature range
Installation length
Shaft diameter
Minimum immersion depth
Stainless steel
0...300 µS/cm
1.25 mm
1.25 mm
3.5 mm

Technical specifications

Shaft material PEEK

Ideal measuring range $15...2.5 \times 10^5 \mu \text{S/cm}$

Temperature range 0...70 °C Installation length 125 mm Shaft diameter 12 mm Minimum immersion depth 30 mm

Technical specifications

Shaft material PEEK

Ideal measuring range 5...2 x 10⁴ µS/cm

Temperature range 0...70 °C Installation length 125 mm Shaft diameter 12 mm Minimum immersion depth 35 mm

Technical specifications

Shaft materialGlassIdeal measuring range1...105 μS/cmTemperature range5...70 °CInstallation length125 mmShaft diameter12 mmMinimum immersion depth30 mm



Ordering information

Conductivity measuring cell (stainless steel) $c = 0.1 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 0.65 m) plug N 6.00921.040 Conductivity measuring cell (stainless steel) $c = 0.1 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 2.0 m) plug N 6.00922.040 Four-wire conductivity measuring cell $c = 0.5 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 0.65 m) plug N 6.00923.080 Four-wire conductivity measuring cell $c = 0.5 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 2.0 m) plug N 6.00924.080 Five-ring conductivity measuring cell $c = 0.7 \text{ cm}^{-1}$ with Pt1000 (fixed cable 0.65 m) plug N 6.00925.100 Three-ring conductivity measuring cell $c = 1.6 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 0.65 m) plug N 6.00929.140 Three-ring conductivity measuring cell $c = 1.6 \text{ cm}^{-1}$ with Pt1000 (fixed cable, 2.0 m) plug N 6.00930.140 OMNIS Measuring Module Conductivity 6.02101.020

Five-ring conductivity measuring cells

Modern five-ring conductivity measuring cells have linearity ranges that are wider than those of classic conductivity measuring cells and require no additional platinization. The current applied to the inner electrode generates a current flow to the outer, grounded electrodes, so that external influences and measuring errors are minimized.

Five-ring conductivity measuring cells supply precise measuring values, independent of positioning in the beaker (wall effect). Interferences with the potentiometric measurements are now a thing of the past; conductivity and pH value can now be measured simultaneously in the same beaker.

The measuring cells are equipped with a fixed cable and plug N for direct connection to the 856 Conductivity Module.

Technical specifications 6.0915.100/6.0920.100

Shaft material PEEK

Ideal measuring range $5...2 \times 10^4 \, \mu \text{S/cm}$ Temperature range $0...70 \, ^{\circ}\text{C}$

Temperature range 0...70 °C

Temperature sensor Pt1000

Installation length 125 mm

Shaft diameter 12 mm

Minimum immersion depth 34 mm

Cell constant 0.7 cm⁻¹

Technical specifications 6.0915.130/6.0920.130

Shaft material PEEK

Ideal measuring range 5...105 µS/cm

Temperature range 0...70 °C

Temperature sensor Pt1000

Installation length 142 mm

Shaft diameter 12 mm

Minimum immersion depth 50 mm

Cell constant 1 cm-1





Ordering information

Five-ring conductivity measuring cell c = 0.7 cm⁻¹ with Pt1000, fixed cable (1.2 m) plug N 6.0915.100 Five-ring conductivity measuring cell c = 1.0 cm⁻¹ with Pt1000, fixed cable (1.2 m) plug N 6.0915.130 Five-ring conductivity measuring cell c = 0.7 cm⁻¹ with Pt1000, fixed cable (2 m) plug N 6.0920.100 Five-ring conductivity measuring cell c = 1.0 cm⁻¹ with Pt1000, fixed cable (2 m) plug N 6.0920.130 Conductivity measuring cell (stainless steel) c = 0.1 cm⁻¹ with Pt1000, fixed cable (1.2 m) plug N 6.0916.040

Conductivity measuring cells for 912/914 meters

80 Four-electrode measuring technique

Historically, conductivity measuring cells consists of two platinized platinum electrodes (Kohlrausch cell). If other electrode materials are used instead of platinized platinum, errors in measurement results due to polarization are most likely to occur.

However, platinization has drawbacks, platinized electrodes are susceptible to encrustation, inclusions and also growth of algae, bacteria or mold leading to changes in the cell constant. Additionally, the two-electrode measuring technique is affected by field effects thus falsifying conductivity values measured close to a wall.

Therefore, Metrohm introduced the multi-electrode measurement technique for their instruments. For the 912/914 meters, an electrode based on four-electrode measurement was developed (6.0917.080). This cell consists of two pairs of electrodes, one pair is used for introducing the generator current, and the other pair is recording the measurement voltage.

The current flowing in the solution causes a drop in the measured voltage. By knowing this voltage and the generator current, the conductance and from that the conductivity can be calculated.

In summary, the four-electrode measuring technique allows an almost polarization free recording of the conductance and determinations are more reproducible. Additionally, it enables a very large measuring range. Another benefit of this sensor is its PEEK shaft which makes the sensor more robust.

Technical specifications

6.0918.040

Shaft material Stainless steel Ideal measuring range 0...300 µS/cm
Temperature range 0...70 °C
Temperature sensor Pt1000
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 35 mm
Cell constant 0.1 cm⁻¹

Technical specifications

6.0917.080

Shaft material PEEK

Ideal measuring range 15...2.5*10⁵ µS/cm

Temperature range 0...70 °C
Temperature sensor Pt1000
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 30 mm
Cell constant 0.5 cm⁻¹

Technical specifications

6.0919.140

Shaft material Glass
Ideal measuring range 10²...106 µS/cm
Temperature range 0...70 °C
Temperature sensor Pt1000
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 40 mm
Cell constant 1.6 cm-1



Ordering information

Conductivity measuring cell (stainless steel) with Pt1000, $c = 0.1 \text{ cm}^{-1}$, fixed cable (1.2 m) plug K

Four-wire conductivity measuring cell with Pt1000, $c = 0.5 \text{ cm}^{-1}$, fixed cable (1.2 m) plug K

Three-ring conductivity measuring cell with Pt1000, $c = 1.6 \text{ cm}^{-1}$, fixed cable (1.2 m) plug K

6.0917.080

6.0919.140

Conductivity measuring cells

82 Conductivity measuring cell c = 0.9 cm⁻¹

- Platinized
- With standard ground-joint 14/15
- Metrohm plug-in head G
- Optimum length for sample changer systems

Technical specifications

Shaft materialGlassMeasuring range1...105 μS/cmTemperature range5...70 °CInstallation length120 mmShaft diameter12 mmMinimum immersion depth16 mm



Ordering information

Conductivity measuring cell for sample changer, $c = 0.9 \text{ cm}^{-1}$, without cable

6.0910.120





Accessories for conductivity measuring cells:

Conductivity standard κ = 12.87 mS/cm (25° C), 250 mL 6.2301.060 Conductivity standard κ = 100 μ S/cm (25 °C), 250 mL with DKD certificate 6.2324.010 Conductivity standard κ = 100 μ S/cm (25 °C), 5 x 30 mL with DKD certificate 6.2324.110

Sensors for stability measurement

The oxidation stability characterizes the resistance of oils and fats and fatty foods to oxidation. It is a standard parameter for quality control in the production of oils and fats in the food industry or for incoming goods inspection in processing plants. Biodiesel and PVC are also subject to oxidation, which can be determined by stability measurement with biodiesel Rancimats or PCV

Thermomats.

The determination of the oxidation stability takes place automatically and is measured by means of a conductivity sensor. The exact temperature is of central importance for the measurement. We offer two temperature sensors for determining the temperature compensation in the reaction vessel. A short version for the measurement in the short test tubes and a longer version for the determination of the temperature compensation in longer biodiesel test tubes.

Conductometric measuring cell for 743 Rancimat, 873 Biodiesel Rancimat, 892 Professional Rancimat, 893 Professional Biodiesel Rancimat, 763 PVC Thermomat and 895 Professional PVC Thermomat

Technical specifications

Shaft material

Measuring range 0...400 μS/cm
Cell constant 1.1 cm⁻¹

Temperature sensor Pt100 for 743 Rancimat, 763 PVC Thermomat, 892 Professional Rancimat or 895 Professional PVC Thermomat

- Sensor class B according to EN 60751 (ITS 90), certified
- Fixed cable with mini DIN plug
- Stainless steel shaft

Temperature sensor Pt100 for 873 Biodiesel Rancimat or 893 Professional Biodiesel Rancimat

- Sensor class B according to EN 60751 (ITS 90), certified
- Fixed cable with mini DIN plug
- Stainless steel shaft

Technical specifications

Shaft material Stainless steel (AISI 304)
Measuring range -200...300 °C
Installation length 175 mm
Minimum immersion depth 20 mm
Shaft diameter 2 mm

Technical specifications

Shaft material Stainless steel (AISI 304)
Measuring range -200...300 °C
Installation length 300 mm
Minimum immersion depth 20 mm
Shaft diameter 2 mm



Ordering information	
Conductometric measuring cell for 743, 763, 873, 892, 893 and 895	6.0913.130
Pt100 temperature sensor short for 743, 763, 892, 895	6.1111.010
Pt100 temperature sensor long for biodiesel Rancimats 873, 893	6.1111.020
Accessories	
SET to determine temperature correction for 743, 763, 892, 895	6.5616.100
SET to determine the temperature correction at 873, 893	6.5616.110
Silicone oil for stability measuring instruments (50 mL)	6.2326.000

Temperature sensor

86 Temperature sensor Pt1000

- Rapid, precise temperature setting
- Available in various lengths (90/125/178 mm)

Technical specifications

Shaft material Glass
Temperature range -50...180 °C
Installation length 90/125/178 mm
Shaft diameter 12 mm
Shaft diameter bottom 5/6.4 mm

Minimum immersion depth 20 mm

Electrode plug-in head Metrohm plug-in head G

Temperature sensor Pt1000 steel

- The glass-free alternative
- Shaft made of PEEK
- For use in non-oxidizing media pH 1 13
- For temperature measurement in semi-solid materials such as cheese, not in frozen meat or similar

Technical specifications


Shaft material PEEK
Temperature range -50...100 °C
Installation length 140 mm
Shaft diameter 12 mm
Shaft diameter bottom (75 mm) 3 mm
Minimum immersion depth 10 mm

Temperature sensor Pt1000 for 909 UV Digester

- Shaft made of glass
- Fixed cable with 2 x B (2 mm) plug

Technical specifications

Shaft material Glass
Temperature range -50...180 °C
Installation length 120 mm
Shaft diameter 12 mm
Shaft diameter bottom 5 mm
Minimum immersion depth 20 mm

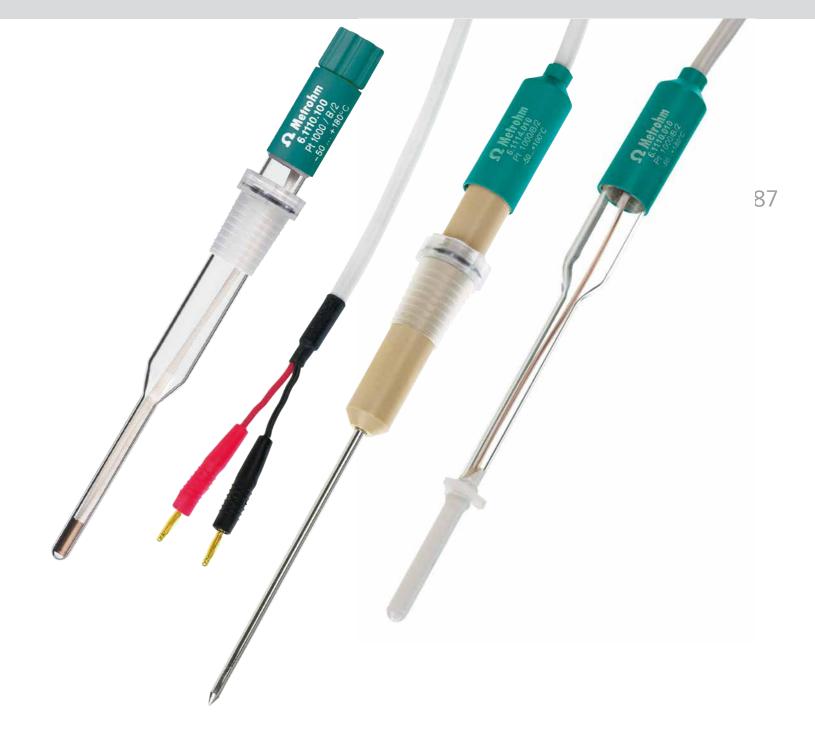

Ordering informationTemperature sensor Pt1000, length 125 mm, without cable6.1110.100Temperature sensor Pt1000, length 178 mm, without cable6.1110.110Temperature sensor Pt1000, length 90 mm, without cable6.1110.120Temperature sensor Pt1000, steel, length 140 mm, fixed cable (1.2 m) plug 2 x B (2 mm)6.1114.010Temperature sensor Pt1000 for 909, length 120 mm, fixed cable (0.5 m) plug 2 x B (2 mm)6.1110.010

Sensor for themormetric titration

Next to potentiometric titration where a suitable electrode is used to measure a potential difference, thermometric titration exists. Thermometric titration monitors the change in temperature caused by the reaction enthalpy of the titrant with the analyte (exothermic or endothermic reaction). For this, a very sensitive temperature sensor is used, a thermistor.

88

The thermistor is manufactured from sintered mixed metal oxide which exhibits a large change of electrical resistance for a small temperature change. The thermistor is then encapsulated in a suitable electrically insulating medium with satisfactory heat transfer characteristics and acceptable chemical resistance. Typically, the encapsulating material is glass, however, when either chemical attack or mechanical stress is expected encapsulation in an epoxy resin might be possible.

Furthermore, a quick electronic circuit is used to maximize the sensitivity for smallest temperature changes. With this, a resolution of 10^{-5} K can be achieved.

Thermometric titration may be the choice where conventional titration sensors are not suitable for the titration environment or when they deliver unsatisfactory titration results.

Technical specifications 6.9011.020

Shaft material Glass
Measuring range 0...60 °C
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 15 mm

6.9011.040

Shaft material PEEK and PVDF coated glass

Measuring range 0...60 °C

Installation length 125 mm

Shaft diameter 12 mm

Minimum immersion depth 15 mm

6.01117.300

Shaft material Glass
Measuring range 0...60 °C
Installation length 125 mm
Shaft diameter 12 mm
Minimum immersion depth 15 mm

Examples for thermometric titration:

- Total acid number (TAN) according to ASTM D8045
- Total base number (TBN)
- Phosphate determination
- Sodium determination
- Titration of plating baths



Ordering informationThermoprobe6.9011.020Thermoprobe HF6.9011.040dThermoprobe6.01117.300

Optical Sensor

This optical sensor for the determination of dissolved oxygen in water or other liquid samples. Compared with conventional, galvanic sensors, the O_2 Lumitrode is much easier to handle and has a longer service life.

Currently, optical sensors for dissolved oxygen measurements are well accepted for the in-situ determination of dissolved oxygen and are already mentioned in the standard ISO 17289 as the method of choice.

O2 Lumitrode

- Low maintenance
- Does not consume oxygen
- Easy to clean
- Can be used for in-situ or continuous dissolved oxygen (DO) measurements
- High accuracy: < 0.15 mg/L at DO concentrations between 0-20 mg/L
- Fast response time, analysis takes < 1 minute

Technical specifications O2 Lumitrode

Shaft material Luran, stainless steel

Ideal measuring range 0...500 % a.s. 0...50 mg/L

Temperature range 0...60 °C Installation length 140 mm Shaft diameter 12 mm

Minimum immersion depth 35 mm

Oxygen measurements with the O2 Lumitrode, examples:

- Dissolved oxygen in tap water / surface water
- Dissolved oxygen in wine
- Dissolved oxygen in liquid dairy products
- Dissolved oxygen in fruit juices
- Dissolved oxygen in acrylic dispersion paint



Ordering information O2 Lumitrode 6.1116.000 Accessories for O2 Lumitrode Calibration vessel O2 Oxygen standard 0% 6.2329.000 O2 Lumitrode replacement set 6.5623.000

Sensor for photometry

Titration with photometric endpoint detection is an integral part of many titration methods. The Optrode is a handy sensor that can be used like any other Metrosensor. Both new and existing titration systems can be equipped with the Optrode. The power supply comes directly via the USB port of a Metrohm instrument (Titrino plus, Ti-Touch, Titrando, USB sample changer). In the case of models without a USB port the power can also be supplied via an optional USB power adapter.

Technical specifications Optrode

Shaft diameter 12 mm
Installation length 135 mm
Shaft material Glass

Measuring range

Dark voltage min. 0 mV
Bright voltage max. 800 mV
Temperature range 0...40 °C
pH range 0...14
Min. immersion depth 30 mm

Optrode

92

- Eight wavelengths: 470, 502, 520, 574, 590, 610, 640 and 660 nm
- 100% solvent resistant (glass shaft)
- Compact and space-saving
- Very easy cleaning
- Easy change of wavelength with a magnet or software controlled (tiamo 2.5 or higher)

Photometric titrations with the Optrode, examples:

- Nonaqueous titrations according to USP and EP
- Determination of the carboxyl end groups (nonaqueous)
- TAN/TBN according to ASTM D974 (nonaqueous)
- Chloride in silicone products (nonaqueous)
- Sulfate determination
- Fe, Al, Ca in cement
- Water hardness (total hardness and Ca/Mg)
- Chondroitin sulfate according to USP



Ordering information

Optrode, fixed cable plug F, USB

6.1115.000

Accessories for Optrode

USB power adapter 5V / 2.1 A (for titrators without USB-connection)
OMNIS adapter cable

6.02117.000 6.02109.000





Accessories for Metrosensors

Accessories for Metrosensors

96 SGJ sleeves for Metrohm electrodes

6.1236.020 SGJ sleeve made of PP, standard

ground-joint 14/15 with O-ring

6.1236.030 SGJ sleeve made of PP, standard

ground-joint 14/15 with O-ring,

for sample changer

6.1236.040 SGJ sleeve made of silicone rubber,

standard ground-joint 14/15

6.1236.050 SGJ sleeve made of EVA, standard

ground-joint 14/15

Other accessories

6.2008.040 Storage vessel made of PP

Length 105 mm Diameter 13 mm

Ground-joint taper or standard

ground-joint 14/15

6.1243.020 Spare ground-joint diaphragm

for Profitrode 6.0255.1XX

(glass sleeve and plastic ring)

6.1243.030 Spare ground-joint for reference

electrodes 6.0726.1XX and

6.0729.1XX

6.2615.050 Electrode holder for 11 electrodes

and 3 \times 50 ml buffer bottles

6.1620.000

6.1620.000 Bottle made of PE

(e.g. for buffer solutions)

Height 95 mm Diameter 31 mm Volume 50 mL













6.1236.040



6.1236.050



6.1236.030



Ion standards, buffer solutions, electrolytes

Ion standards (traceable to NIST)

98

Ion	Concentration	Amount	Ordering number
KCl	0.1000 ± 0.0005 mol/L	250 mL	6.2301.060
Conductivity standard	κ = 12.87 mS/cm (25 °C)	250 mL	6.2301.060
Conductivity standard	for USP <645> and EP 2.2.38	250 mL	6.2324.010
with DKD-certificate	κ = 100 μS/cm (25 °C)	5 x 30 mL (sachets)	6.2324.110

Buffer and calibration solutions (traceable to NIST)

Article		Amount	Ordering number
Ready-to-use buffer solution	pH 4.00	500 mL	6.2307.100
in bottles, colored, with	pH 7.00	500 mL	6.2307.110
paper seal	pH 9.00	500 mL	6.2307.120
Ready-to-use buffer solution	pH 4.00	30 x 30 mL	6.2307.200
in sachets with DKD-certificate	pH 7.00	30 x 30 mL	6.2307.210
	pH 9.00	30 x 30 mL	6.2307.220
	pH 4.00, 7.00, 9.00	10 x 30 mL each	6.2307.230
Redox standard			
yields with reference electrode $Ag/AgCI/c(KCI) = 3 mol/L$		250 mL	6.2306.020
$U = +250 \pm 5 \text{ mV } (20 \text{ °C}); c$	can also be used as buffer pH 7		

Storage solution, cleaning solution, pHit Kit

Electrolyte		Amount	Ordering number
Storage solution	for combined pH glass electrodes with	250 mL	6.2323.000
	reference electrolyte c(KCl) = 3 mol/L		
pHit kit	Care kit for electrodes, containing 50 mL	50 mL each	6.2325.000
	each of cleaning solution, reference		
	electrolyte c(KCl) = 3 mol/L, storage		
	solution, and		
	2 storage vessels		
Cleaning solution	As addition to the pHit Kit to clean	3 x 50 mL	6.2325.100
	combined pH glass electrodes		





Electrolytes

Electrolyte		Amount	Ordering number
KCl 3 mol/L	for Ag/AgCl reference systems	50 mL	6.2308.050
		250 mL	6.2308.020
		1000 mL	6.2313.000
KCl gel 3 mol/L	only for Ag/AgCl reference systems in voltammetry	50 mL	6.2308.060
KCl saturated	for storage of gel electrodes	250 mL	6.2308.000
KNO3 1 mol/L	reference electrolyte for combined Ag-electrodes and bridge electrolyte for Ag/AgCl reference systems	250 mL	6.2310.010
NH ₄ NO₃ 1 mol/L	reference electrolyte for 6.0510.100 combined Ca-ISE	50 mL	6.2327.000
CH₃COOLi 1 mol/L	reference electrolyte for 6.0510.110 combined K-ISE	50 mL	6.2328.000
TEABr (Tetraethylammonium bromide) 0.4 mol/L in ethyl- ene glycol	bridge electrolyte for titrations in nonaqueous solutions and reference electrolyte for Solvotrodes	250 mL	6.2320.000
NH ₄ Cl 0.1mol/L	inner measuring electrolyte for NH ₃ -electrodes	50 mL	6.2316.030
Idrolyte	for 6.0224.100 Biotrode or for pH-measurement > 80 °C with Unitrode/Syntrode	250 mL	6.2308.040
Porolyte	for 6.0235.200 Porotrode	50 mL	6.2318.000



Electrical connections

Connection of pH electrodes, ion-selective electrodes (ISE) and metal electrodes on Metrohm instruments

Electrode plug-in head	Cable Order number		Measuring device
	Plug-in head G – plug F, 1 m	6.2104.020	For pH/ISE and Ind measuring
			inputs, Titrinos and Titrandos,
Plug-in head G	Plug-in head G – plug F, 2 m	6.2104.030	pH-/ion meter from 691
	Plug-in head G – plug F, 3 m	6.2104.040	Ind.
	Plug-in head G – plug E (DIN 19262), 1 m	6.2104.050	For Metrohm pH Meter < 691
	Plug-in head G – plug E (DIN 19262), 2 m	6.2104.060	
	Plug-in head G – plug E (DIN 19262), 3 m	6.2104.070	
	Electrode cables for generator electrodes	6.2104.120	KF Coulometer Gen. EL
	6.0342.110, 6.0344.100 and 6.0345.100		
pH electrodes with fixed cable, plug B (2 mm)	Adapter plug B (2 mm)/4 mm	6.2103.150	Titrinos (Pt1000 only) pH Meter ≤ 744 (Pt1000 only)
pH electrodes with	Adapter plug B (4 mm)/2 mm (red)	6.2103.130	780/781/Titrandos/
fixed cable plug B	Adapter plug B (4 mm)/2 mm (black)	6.2103.140	Ti-Touch/91X meters
(4 mm)			Pt1000 (2 mm)
			Temp.
Plug-in head U	Plug-in head U – plug F + 2 x B (2 mm), 1 m	6.2104.600	Ind.
	Plug-in head U – plug F + 2 x B (2 mm), 2 m	6.2104.610	Temp.

Connection of conductivity measuring cells and temperature sensors to Metrohm instruments

Electrode plug-in head	Cable	Order number	Measuring device
Plug-in head G	Plug-in head G – plug 2 x B (4 mm), 1 m	6.2104.080	712 Conductometer, measuring inputs Pt 100/ Pt1000
	Plug-in head G – plug 2 x B (4 mm), 2 m	6.2104.110	(4 mm) Cond. Cell Cell Cell Cell Cell Cell Cell Cel
	Plug-in head G – plug 2 x B (2 mm), 1 m	6.2104.140	780/781/Titrandos/ Ti-Touch/91X meters
0.0	Plug-in head G – plug 2 x B (2 mm), 2 m	6.2104.150	Pt1000 (2 mm) Temp.

Connection of reference electrodes and separate metal electrodes on Metrohm devices

Electrode plug-in head	Cable	Order number	Measuring device
Plug-in head G	Plug-in head G – plug F, 1 m	6.2104.020	For «Ind.» measuring inputs of Titrinos and for connection
(6.0729.XXX)	Plug-in head G – plug F, 2 m	6.2104.030	to a Metrohm differential amplifier
20.C	Plug-in head G – plug F, 3 m	6.2104.040	Ind.
Plug-in head B	Plug-in head B (4 mm) – plug B (4 mm), 1 m	6.2106.020	For measuring input «Ref.»
20.00 0.00 0.00 0.00	Plug-in head B (4 mm) – plug B (4 mm), 2 m	6.2106.060	
42 Mein 6.0726.0 9.AgCl/0	Plug-in head B (4 mm) – plug B (4 mm), 3 m	6.2106.050	Ref.

Connection of Metrohm electrodes with plug-in head G to devices made by other manufacturers

Electrode plug-in head	Cable	Order number	Measuring device
Plug-in head G	Plug-in head G – BNC plug, 1 m	6.2104.090	Orion, Beckman, Corning, EDT, Fisher, Hanna, Mettler- Toledo, Jenway, Philips, Radiometer, Mitsubishi, SI Analytics, Crison, Kyoto/KEM
	Plug-in head G – LEMO	6.2104.160	Mettler
	Plug-in head G – plug E (DIN 19262), 1 m	6.2104.050	Older Metrohm devices, WTW, Knick, SI Analytics

102 Connection of Metrohm electrodes to OMNIS measuring modules

Electrode plug-in head	Cable	Order number	Measuring device
Plug-in head Q	Plug-in head Q – plug P, 0.55 m	6.02104.300	For a digital measuring module of OMNIS
	Plug-in head Q – plug P, 1.5 m	6.02104.310	
Plug-in head G of pH electrode, ISE, metal electrodes	Plug-in head G – plug P, 0.55 m (green)	6.02104.000	For measuring input "INPUT 1" and "INPUT 2" of an analog measuring module of
	Plug-in head G – plug P, 1.5 m (green)	6.02104.010	OMNIS
	Adapter cable for electrodes with fixed cable	6.02109.000	
Plug-in head G of temperature sensor	Plug-in head G – plug P, 0.55 m (red)	6.02104.020	For measuring input "INPUT 1" and "INPUT 2" of an analog measuring module of
	Plug-in head G – plug P, 1.5 m (red)	6.02104.030	OMNIS
Plug-in head G of electrode used for polarized measure-	Plug-in head G – plug P, 0.55 m (blue)	6.02104.040	For measuring input "INPUT 1" of an analog measuring module of OMNIS
ment	Plug-in head G – plug P, 1.5 m (blue)	6.02104.050	
	Adapter cable for electrodes with fixed cable	6.02109.010	
Plug-in head U	Plug-in head U – plug P, 0.55 m (green/red)	6.02104.600	For measuring input "INPUT 1" and "INPUT 2" of an analog measuring module of
	Plug-in head U – plug P, 1.5 m (green/red)	6.02104.610	OMNIS
Plug-in head B	Plug B (4 mm) – Plug B (2 mm), 0.55 m	6.02105.000	For measuring input "REF" of the analog measuring module of OMNIS
6.0726.100 g./AgCi/U8i'C Ostes Pulland	Plug B (4 mm) – Plug B (2 mm), 1.5 m	6.02105.010	

How is a Metrosensor made?

Our accuracy is not accidental ...

Always keep cool!

Many years of experience and a steady hand with raw materials guarantee that our chemistry is always correct. The right composition of the glass mixture and the greatest possible care during the melting process ensure the perfect quality of the membrane glass.







103

We know how to do it!

Our employees need the right feeling when fusing the membrane glass with the electrode body. That this is not just a matter of luck can be seen from our electrodes at the first glance..









Our employees are only human!

In some manufacturing processes, such as grinding our fixed ground-joint diaphragms, even a practiced eye no longer stands a chance. Such tasks are carried out with unrivalled accuracy by the most modern machines.



Tried and tested electrodes!

Before our electrodes leave our premises they are subjected to a further wet-chemistry computer-supported check. At its conclusion we provide a written confirmation so that you can have complete confidence in our electrodes: each electrode receives its own test certifi-



Certificate of origin:

Precision and guaranteed reliability - Metrohm stands for the highest quality in ion analysis. Just convince yourself!

1. Fundamentals of potentiometry

1.1. Electrode construction

In potentiometry the measuring setup always consists of two electrodes: the measuring electrode, also known as the indicator electrode, and the reference electrode. Both electrodes are half-cells. When placed in a solution together they produce a certain potential. Depending on the construction of the half-cells, the potential produced is the sum of several individual potentials. Potentialdetermining transitions always occur at the phase boundaries, e.g. between the solution and the electrode surface.

pH electrode

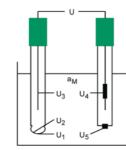


Figure 1: Schematic diagram of a pH electrode

Measuring electrode – glass electrode (left)

- U_1 = Galvani potential between measuring solution and glass membrane
- U₂ = Galvani potential between glass membrane and inner electrolyte
- U_3 = Galvani potential between inner electrolyte and inner reference electrode

Reference electrode – silver/silver chloride (right)

U₄ = Galvani potential of reference electrode

 U_5 = Diaphragm potential (diffusion potential)

 a_M = Activity of measured ion in sample solution

The potentials U_2 , U_3 and U_4 can be kept constant by a suitable electrode construction. Constructive measures and the selection of a suitable reference electrolyte ensure that U₅ is also kept as constant as possible. Ideally the measured potential should depend only on the potential between the glass membrane and the solution. (1) For practical reasons the half-cells of the measuring electrode and the reference electrode are normally contained in a single electrode; this is then known as a combined pH electrode.

Redox electrode

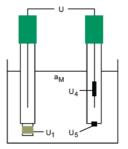


Figure 2: Schematic diagram of a redox electrode

Measuring electrode – metal electrode (left)

 U_1 = redox potential between measuring solution and metal surface

Reference electrode – silver/silver chloride (right)

U₄ = Galvani potential of reference electrode

 U_5 = Diaphragm potential (diffusion potential)

 a_M = Activity of measured ion in sample solution

For metal electrodes the potential forming transitions U₂ and U₃ of the pH electrodes do not exist. Depending on the particular application, it may be possible to use a pH glass electrode as the reference electrode instead of the silver/silver chloride reference electrode. In the combined redox electrodes and Titrodes from Metrohm the halfcells are also contained in a single electrode.

1.2. From the measured potential to the ion concentration

As each ion is surrounded by ions with the opposite charge, it is – to put it simply – no longer as effective as a free ion (see Debye-Hückel law). This affects both the reactivity and the size of the potentials at the measuring electrode. The activity of the measuring ion a_M, which is also used in the Nernst equation, is linked to the normally interesting analytical concentration c_M via the activity

$$a_M = \gamma * c_M$$

The mathematical relationship between the activity a_M of a measuring ion in solution ions and the potential measured between the reference electrode and the measuring electrode is described by the Nernst equation. This applies only for the (ideal) case in which an electrode only responds to a single type of ion. Potentials U_2 to U_5 for pH electrodes and U_4 and U_5 for redox electrodes, which are normally constant, appear as potential U_0 in the Nernst equation.

$$U = U_0 + \frac{2.303 * R * T}{z * F} * log a_M$$

(Nernst equation) (2)

U = measured potential

 U_0 = temperature-dependent standard potential of electrode

R = general gas constant 8.315 J mol-1 K⁻¹

T = temperature in Kelvin

z = ionic charge including sign

F = Faraday constant 96485.3 C mol⁻¹

The term in the Nernst equation in front of the logarithm is known as the Nernst potential U_N (also Nernst slope).

$$U_N = \frac{2.303 * R * T}{z * F}$$

(Nernst potential) (3)

Its value is 0.059 V at T = 298.15 K and z = +1. As a factor in the Nernst equation it represents the theoretical electrode slope. U_N corresponds exactly to the alteration in potential caused by increasing the activity a_M by a factor of ten. From the equation it can be seen that the electrode slope for electrodes that respond to ions with a double charge (z = 2) is only half the size of that for electrodes for ions with a single charge (z = 1). In addition, the sign for cation- and anion-sensitive measuring electrodes is different, as z also takes the charge on the ion into account. The Nernst potential is directly dependent on the temperature (see Equation 3). This is why it is

absolutely necessary to take the temperature into account in all direct potentiometric measurements, as otherwise no correct results will be obtained.

pH value

In practice – particularly when measuring the acid/ base equilibrium – the term pH, introduced by Sörensen in 1909, is frequently used instead of the activity of the measuring ion a_M :

$$pH = -log a_{H^+}$$

(Definition of the pH value) (4)

The pH value is the negative common logarithm of the hydrogen ion activity of a solution. The term p is frequently used for the simplified presentation of very large or small values. In a similar way pNa $^+$ can be used for the activity of sodium ion, or pK $_A$ as acid constant or pK $_B$ as base constant for reaction constants. In each of these cases what is meant is the negative common logarithm of the particular value. If this definition is inserted in the Nernst equation then we obtain for the measured potential U:

$$U = U_0 - \frac{2.303 * R * T}{z * F} * pH$$

(pH value and potential) (5)

Redox potentials (metal electrodes)

In a similar way to the Nernst equation (Equation 2) the equation for the activity-dependent potential is obtained as follows:

$$U = U_0 + \frac{2.303 * R * T}{z * F} * log \frac{a_{ox} * a_{H^+}}{a_{red}}$$

Equation 6 usually allows the potential generated by a redox pair at the measuring electrode to be calculated. As protons are involved in most redox reactions, the measured potential depends on the pH. If proton reactions cannot be excluded then the pH should also be determined or adjusted to a defined value.

1.3. Measuring electrodes

1.3.1. pH glass electrodes

How does a pH glass electrode work?

The glass membrane of a pH glass electrode consists of a silicate framework containing lithium ions. When a glass surface is immersed in an aqueous solution then a thin solvated layer (gel layer) is formed on the glass surface in which the glass structure is softer. This applies to both the outside and inside of the glass membrane. As the proton concentration in the inner buffer of the electrode is constant (pH 7), a stationary condition is established on the inner surface of the glass membrane. In contrast, if the proton concentration in the measuring solution changes then ion exchange will occur in the outer solvated layer and cause an alteration in the potential at the glass membrane. Only when this ion exchange has achieved a stable condition will the potential of the glass electrode also be constant. This means that the response time of a glass electrode always depends on the thickness of the solvated layer. Continuous contact with aqueous solutions causes the thickness of the solvated layer to increase continuously – even if only very slowly – which results in longer response times. This is

why conditioning the electrode in a suitable electrolyte is absolutely necessary to ensure an initial solvated layer condition that is as stationary as possible so that results can be obtained that are as reproducible as possible.

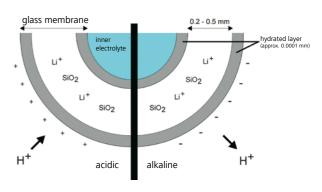


Figure 3: The silicate skeleton of the glass membrane contains lithium ions, among other things. During the formation of the solvated layer at the glass surface these are partly replaced by protons. If the concentration of the protons in the solution changes then a new stationary condition must again be achieved in the solvated layer; this results in a change in potential at the glass membrane.

Table 1: Overview of the different electrode membrane glasses used by Metrohm Ltd $\,$

Application	U glass	T glass	M glass	Aquatrode glass	_
pH range	(green) 014	(blue) 014	(colorless)	(yellow) 013	(yellow) 013
Temperature range	080 °C	080 °C	060 °C	080 °C	080 °C
continuous	0100 °C				
short-term					
Membrane	Electrodes with	Electrodes with	Electrodes with	Large surfaces	Electrodes with
surface	large membrane	medium to large	small membrane		medium to large
	surface	membrane	surface (micro-		membrane
		surface (mini-	electrodes)		surface
		electrodes)			
Special features	For strongly	Measurements in	Measurements in	Responds very	Quick response,
	alkaline solutions,	non-aqueous	small-volume	quickly, so	excellent stability
	long-term meas-	sample solutions	samples		in continuous use
	urements and			for measurements	
	measurements at			in ion-deficient or	
	high temperatures			weakly buffered	
				solutions	
Membrane resist-	< 500	< 150	< 120	< 250	< 250
ance (M Ω)					
With reference to					
sphere membrane 10.5 mm diameter					
10.5 min diameter					

Why are there different types of glass for pH electrodes?

Different demands are placed on a pH glass electrode depending on the particular application. Various properties such as response time, thermal resistance, chemical stability, shape, size and electrical properties must all be taken into account in order to have an optimal electrode available to solve each problem. In order to be able to do justice to the numerous applications, different glasses are available with different properties (see table 1).

Why must a pH glass electrode be calibrated?

The potential of a measuring electrode can always only be given relative to that of a reference electrode. To be able to compare systems, the electrode zero point is defined as being 0 mV for pH 7 and 298.15 K or 25 °C. The electrode slope, i.e. the alteration in the measured value with the pH, is given by the Nernst equation and at 25 °C is 0.059 V per pH 1. These are ideal values from which Metrosensor electrodes only differ slightly. The electrode zero point is \pm 0.015 V. The electrode zero point and the electrode slope may change as a result of the aging of the glass membrane or changes, (e.g contamination) on the diaphragm. For this reason the pH meter must be adapted to the characteristics of the electrode, i.e. calibrated, at regular intervals by using buffer solutions.

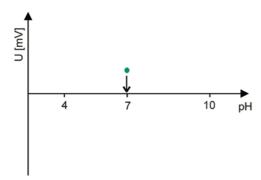


Figure 4: In the first calibration step with buffer pH = 7 the variation from the electrode zero point (= asymmetry potential) is determined and corrected.

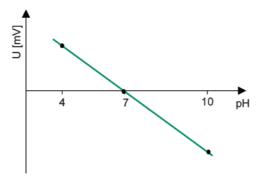


Figure 5: In the second calibration step with another buffer solution the electrode slope is determined and expressed as a percentage of the theoretical value of 0.059 V (at 25 °C).

The electrode zero point is set first (pH 7 corresponding to 0 mV for Metrosensor pH electrodes). The second and further buffer solutions are used to determine the slope of the pH electrode. This slope is expressed as a percentage of the theoretical value (100% = 0.059 V per ΔpH = 1 at 25 °C). In order to minimize subsequent measuring errors, care should be taken that the expected measured value of the sample solution always lies within the pH range covered by the buffer solutions. Modern pH and ion meters such as the 780 pH Meter, the 781 pH/lon Meter, the 913 pH meter and 914 pH/conductivity meter do not require any manual settings to be made. The buffer solutions are recognized automatically and can be presented in any sequence.

Calibration always includes a check of the measuring electrode. The calibration buffers have a medium acid-base concentration and their ionic strength is approximately that of the most common sample solutions. The dependency of the electrode slope on the temperature means that the calibration and measuring temperatures must be known. Information about the electrode condition is provided by the electrode slope, electrode zero point, response time of the signal and its streaming dependency. With the Metrohm 781 pH/lon Meter and 780 pH Meter an automatic electrode test can be carried out; this provides an exact statement of the electrode condition and often allows a source of error to be localized.

pH and temperature - an inseparable couple!

The temperature has a considerable influence on the pH value and the pH measurement. If an electrode is calibrated at 25 °C then it should be capable of linear measurement throughout the whole pH range and provide correct results. However, if the electrode is then used at a different temperature the electrode slope will change - in accordance with the Nernst equation - and possibly the electrode zero point as well. The point at which the two calibration curves (without correction) for different temperatures intersect is known as the isothermal intersection point. Thanks to the optimized inner buffer and «Long Life» reference system precise measurements can be made with Metrosensor pH electrodes at different temperatures. This means that, although calibration is only carried out at a single temperature, measurements can then be made throughout the whole temperature range. The real behavior of Metrosensor pH electrodes varies from the ideal behavior by maximum ± 15 mV. Nevertheless it is still true that the accuracy of the measurement is increased when the electrode is calibrated at the temperature to be used for the subsequent measurements. At T = 298.16 K and z = 1, the Nernst potential U_N is equal to 59.16 mV. For other temperatures it can be corrected in the Nernst equation by using Table 2. Modern pH meters automatically take the temperature dependency of the Nernst potential into account if a temperature sensor is connected. In principle, within the context of GLP/ISO recording and documentation of the temperature is required for all measurements.

However, it must be remembered that a pH meter can only correct the temperature behavior of the electrode and never that of the solution to be measured. For correct pH measurements it is essential that the pH is measured at the temperature at which the sample was taken. For example, sodium hydroxide c(NaOH) = 0.001 mol/L at 0 °C has a pH of 11.94, at 50 °C it is pH = 10.26 and only at 25 °C is it pH = 11.00. This change in pH is caused by the dependency of the ionic product of water on the temperature.

In some conventional electrodes the temperature sensor is not located in the immediate vicinity of the membrane, i.e. in the electrode foot. This means that it cannot measure the temperature of the solution correctly and that the pH compensation will be incorrect as the temperature

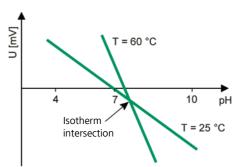


Figure 6: Isothermal intersection point

and pH are not measured at the same location. In modern pH electrodes the temperature sensor should be located within the electrode in the immediate vicinity of the glass membrane. This is the only way in which an accurate pH measurement is possible. If the sensor is located outside the membrane then problems when cleaning the electrode could easily occur.

Table 2: Dependency of the Nernst potential U_N on the temperature

Temperature T (°C)	Slope U _N (mV)	Temperature T (°C)	Slope U _N (mV)
0	54.20	50	64.12
5	55.19	55	65.11
10	56.18	60	66.10
15	57.17	65	67.09
20	58.16	70	68.08
25	59.16	75	69.07
30	60.15	80	70.07
35	61.14	85	71.06
37	61.54	90	72.05
40	62.13	95	73.04
45	63.12	100	74.03

How to store a pH glass electrode?

The swelling of the glass surface is indispensable for the use of glass as membrane for pH glass electrodes; without this solvated layer, no pH measurement would be possible. Glasses for pH glass electrodes are optimized in such a way that only protons can penetrate into the glass membrane. However, because of the very slow but steady swelling of the glass, it is unavoidable that also other ions penetrate into the glass, e.g. sodium and potassium ions. At higher concentrations, these lead to

the so-called alkali error of the glass electrode. This means that the measured value is falsified at comparatively low proton concentrations. If the glass electrode is stored for a very long time in a strong solution of potassium or sodium, this leads to prolonged response times of the glass membrane since the protons must expulse the «added ions» from the solvated layer.

One of the most used electrolytes for pH measurement is c(KCI) = 3 mol/L, since the aequitransferent KCl causes only a very small diffusion potential at the diaphragm and is also economical. Normally a combined pH glass electrode is stored in c(KCI) = 3 mol/L only for this reason, as one wants to have it ready for immediate use without conditioning the diaphragm. However, on a long-term basis the storage in KCl affects the glass, since it leads to ever longer response times. For the membrane glass, storage in distilled water would be optimal, but then the diaphragm would have to be conditioned for several hours. The patented storage solution for combined pH glass electrodes (6.2323.000) solves exactly this problem. If a combined pH glass electrode is kept in this solution, the glass membrane remains unchanged regarding

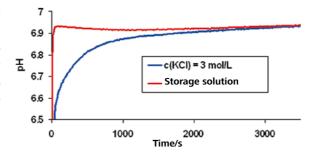


Figure 8: pH measurement in $c(NaHCO_3) = 0.05$ mmol/L. A glass of the Aquatrode stored in the storage solution shows a substantially shorter response time than an electrode glass of the same type stored during the same period in KCl.

response time and alkali error. Moreover, if one uses c(KCI) = 3 mol/L as the reference electrolyte, the optimized composition of the storage solution keeps the pH glass electrode ready for measurement. Conditioning before the measurement is not necessary, no matter for how long the electrode has been stored.

Table 3: The correct storage of pH glass electrodes

Electrode resp. reference electrolyte	Storage
Separate pH glass electrode	Distilled water
Combined pH glass electrode with c(KCl) = 3 mol/L, Porolyte	6.2323.000 Storage solution
Combined pH glass electrode with another reference electrolyte (Idrolyte, non aqueous)	In the respective reference electrolyte
Gel (spearhead electrode), Ecotrode Gel	6.2308.000 Electrolyte solution c(KCl) = sat.

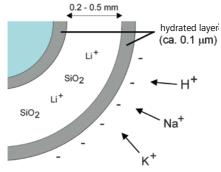


Figure 7: Cross-section of a pH glass membrane. If several kinds of cations are present in the measuring solution, these compete for the free spaces in the solvated layer. Especially potassium and sodium can penetrate into the glass membrane and prolong the response time.

Troubleshooting

The cause of most problems is not to be found in the measuring electrode and its glass membrane, but rather in the reference electrode, as much more critical diaphragm problems can occur there. To avoid incorrect measurements and to increase the working life, attention must still be paid to the following possible sources of error:

Table 4: Possible sources of error and their remedies for pH glass electrodes

Source of error	Effects	Action	Alternatives
HF-containing solutions	Etching and dissolution of the glass membrane → corrosion potential during the measurement/short		Use of the Sb-electrode
High pH value and high alkali content	working life Increased alkali error → pH too low		Use of electrodes with U
High temperatures	Rapid rise in membrane resistance by aging → increased polarizability and drift		Use of electrodes with U glass
Measurements at low temperature	High membrane resistance → polarization effects		Use of electrodes with T glass and Idrolyte as reference electrolyte
Dry storage	Zero point drift	Store in water overnight	Store in storage solution 6.2323.000 or reference electrolyte
Reaction of a solution component with the glass	Slow response, zero point shift, slope reduction		Try other glass types
Non-aqueous media	Reduced sensitivity	Store in water	T glass/non-aqueous electrolyte solution
Deposition of solids on membrane surface	Slow response, zero point shift, slope reduction	Solvent or strong acids	
Electrostatic charging	Slow response	No dab-drying of the electrode	Grounding of measuring instrument
Deposition of proteins on membrane surface	Slow response, zero point shift, slope reduction	5% pepsin in 0.1 mol/L HCl	

Possible sources of error and care information for diaphragm problems are given in Section 1.4. for reference electrodes.

1.3.2. Metal electrodes

How does a metal electrode work?

Metal electrodes have an exposed metal surface. If ions of this metal are contained in the sample solution then an equilibrium is formed at the metal surface that e.g. the Ag/Ag+ equilibrium at a silver surface has a value depends on the concentration of the metal ions in the of E⁰ = 0.7999 V (25°C). If the sample solution does not solution (see «Theory of the electrical double layer» in electrochemistry textbooks). Metal ions are accepted by the metal surface and simultaneously released into the

This concentration-dependent equilibrium is characterized by a corresponding potential E⁰ (Galvani potential), contain any ions of the corresponding metal then metal electrodes can still form a Galvani potential if a redox reaction occurs in the sample solution.

 $S_{ox} + n * e^- \longleftrightarrow S_{red}$

$$Me \longleftrightarrow Me^{n+} + n * e^{-} \qquad E^{0} = \dots$$

metal ions are released from the metal; in this case the metal surface only acts as a catalyst for the electrons. As gold and platinum electrodes are to a large extent $Cl_2(g) + 2e^- \rightarrow 2 Cl^$ chemically inert, they are used for the measurement of redox potentials. Silver electrodes are only used as indicator electrodes for titrations.

Calibrating a metal electrode

Redox-buffer solutions (6.2306.020) are used for quickly checking metal or redox electrodes. As the potential measured in a redox buffer solution is insensitive to the electrode's surface condition, contamination of the metal electrode is often not recognized. For this reason redox-buffer solutions are rather more suitable for checking the reference electrode. If the potential is displaced then the metal electrode is contaminated, the redox buffer partly oxidized or the functioning of the reference electrode is affected. Under no circumstances should the indicated potential be set to the theoretical value.

If measurements are made in weakly redox-buffered solutions then a suitable pretreatment of the metal electrode is recommended to adapt the surface condition as much as possible to the measurement conditions (abrasive pretreatment: carefully clean the electrode with abrasive paste). The reference electrode can either be checked against a second reference electrode that has already been checked in buffer solutions 4 and 7 (response behavior and reference potential) or by using the redox buffer.

The electrode surface is inert to the redox reaction. No In the literature the so-called standard redox potentials E⁰ can usually be found.

$$Cl_2 (g) + 2e^- \rightarrow 2 Cl^ E^0 = + 1.359 \text{ V}$$

 $Fe^{3+} + e^- \rightarrow Fe^{2+}$ $E^0 = + 0.771 \text{ V}$
 $Cd^{2+} + 2e^- \rightarrow Cd^{2-}$ $E^0 = -0.403 \text{ V}$

The zero point of these systems is defined (arbitrarily) with the standard hydrogen electrode (SHE) which is assigned a standard potential of 0 mV. If electrons are released by a redox system to the SHE then this is reduced and the redox pair receives a negative sign; if electrons are accepted then the SHE is oxidized and the result is a redox potential with a positive sign. The standard hydrogen reference electrode is difficult to handle. The specifications of the SHE stipulate that a platinized platinum wire must be used; this is located in a stream of hydrogen gas at a partial hydrogen pressure of 1.0 bar, and that the activity of the hydrogen ions in the solution in which the platinized platinum wire is immersed is to be exactly 1.00 mol/L. The normal alternative is the Ag/ AgCl/KCl reference electrode, which has a potential E⁰ = +207.6 mV at c(KCl) = 3 mol/L and T = 25 °C. The Metrohm redox standard (6.2306.020) can be used for checking separate and combined metal electrodes. Platinum and gold electrodes together with the Ag/AgCl/ KCI reference electrode (c(KCI) = 3 mol/L and T = 20 °C) produce a potential of $\pm 250 \pm 5$ mV.

Table 5: Measuring data for 6.2306.020 redox standard as a function of the temperature

Temp. (°C)	10	20	25	30	40	50	60	70
$mV \pm 5$	+ 265	+ 250	+ 243	+ 236	+ 221	+ 207	+ 183	+ 178
$pH \pm 0.05$	7.06	7.02	7.00	6.99	6.98	6.97	6.97	6.98

If instead of an Ag/AgCl/KCl reference electrode c(KCl) = 3 mol/L an Ag/AgCl/KCl reference electrode c(KCl) = sat. is used for the measurement then at 25 °C a correction of +10 mV must be applied; if the measurement is made using an Hg/Hg₂Cl₂/KCl calomel reference electrode, which for toxicological reasons is no longer available solution.

from Metrohm, the correction to be applied is -37 mV. The Titrodes are checked by a standard titration as no suitable calibration or buffer solutions are available. For example, the certified ion standard c(NaCl) = 0.1 mol/L (6.2301.010) can be titrated with a silver nitrate standard Troubleshooting 113

Table 6: Problems encountered when measuring with metal electrodes

Electrode	Source of error	Effects	Cleaning	Alternatives
Ag	Electrode poisons	Passivation of Ag layer →	Cleaning with abrasives	
	such as S ²⁻ , I ⁻ , Br ⁻	slow response		
Pt/Au	Fats or oils	Isolating layer → slow	Cleaning with solvent	
		response, incorrect potential		
	Weakly redox-	Adsorbed ions on the	Abrasive, oxidative (for	Use of Au or Pt
	buffered solution	surface (e.g. oxides) →	oxidizing solutions) or	
		slow response	reducing (for reducing solu-	
			tions) pretreatment	
	COD determination	Deactivation of Pt		Use of Au

1.3.3. Ion-selective electrodes

How does an ion-selective electrode work?

An ion-selective electrode (ISE) can selectively recognize an ion in a mixture of ions in a solution. There are various types of ion-selective electrodes, the most commonly used ones are:

Glass membrane with	framework of silicate glass interstitial sites for H+ and Na+
Crystal membrane	crystal lattice containing defined gaps for the ion to be measured

Polymer membrane polymer membrane containing a molecule (= ionophore) that only binds the ion to be measured

In contrast to metal electrodes, an ISE does not measure in the sample solution then this ion can penetrate the membrane. This alters the electrochemical properties of the membrane and causes a change in potential. One hundred percent selectivity for exactly one type of ion is only possible on rare occasions. Most ion-selective electrodes have «only» a particular sensitivity for a special type of ion, but also often react with ions with similar chemical properties or a similar structure (see Table 7). This is why the cross-sensitivity to other ions that may be contained in the sample solution must always be taken into consideration when selecting an ISE. One of the

best-known examples of such a cross-sensitivity is the so-called alkali error of pH glass electrodes. With some types of glass the linear range does not extend throughout the whole pH range from 0 to 14 and at high pH values a departure from linear behavior can be observed. The reason for this is that at very low H+-concentrations any alkali ions present in the solution (possibly released from the walls of the vessel) will falsify the measured value. Unfortunately there are only a very few ion-selective electrodes that have a linear range similar to that of pH glass electrodes. The use of an ISE is normally restricted to a concentration range of 6 to 8 powers of ten. If an ISE is used for a measurement right at the limit of the linear range then the Nernst equation (Eq. (5), Section 1.2.) must be extended by the contribution made by the particular interfering ion for the evaluation of the measured potential:

(Nikolsky equation) (9)

K_S is the so-called selectivity coefficient of the ion-selective electrode for interfering ion S. This is a factor that describes the influence of the interfering ion in relationship to the ion to be measured. These selectivity coefficients are known for the most important interfering ions for an ISE and therefore a simple estimation can be made as to whether an interfering ion contained in the sample solution will influence the measured value or not.

Direct measurement or standard addition?

The question often arises as to which determination method is most suitable for a particular sample. In principle there are three different ways of carrying out an ion measurement with ion-selective electrodes:

Direct measurement

Direct measurement is chiefly of benefit with high sample throughputs or with a known sample solution of a simple composition. The ion-selective electrode is calibrated with special standard solutions of the ion to be equation. measured before the measurement itself in a similar way to the calibration of a pH glass electrode and can then be ISA and TISAB – when and why? used for several determinations in series.

Standard addition

Standard addition is recommended whenever a determination only needs to be carried out occasionally or when the composition of the sample is unknown. Defined volumes of a standard solution of the ion to be measured are added to the sample solution in several steps. The concentration in the original solution can then be calculated from the initial potential and the individual potential steps after the addition of the standard. The advantheir addition. tage of standard addition is that the ISE is calibrated directly in the sample solution, which eliminates all matrix effects.

Sample addition

Similar to standard addition, with the difference that defined volumes of the sample solution are added to a defined amount of an ion standard.

Modern ion meters such as the 781 pH/lon Meter from Metrohm can carry out these addition methods automatically. The addition of the standard or sample solution is automatically controlled from the ion meter – by pressing a single key – and evaluated by using the Nikolsky

The activity coefficient of an ion (Section 1.2.) is a function of the ionic strength. For this reason care must be taken that ion-selective measurements are always carried out in solutions with approximately the same ionic strength. In order to achieve this, the so-called ISA solutions (Ionic Strength Adjustor) or TISAB solutions (Total Ionic Strength Adjustment Buffer) should be added to the sample solution (see Table 7). These are chemically inert and have such a high ionic strength that the ionic strength of the sample solution can be neglected after

Table 7: Interfering ions and recommended ISA and TISAB solutions for ion-selective electrodes

lon	Membrane material	pH range ¹	ISA or TISAB ²	Most important interfering ions ³	Remarks
Ag+	Crystal	28	$c(KNO_3) = 1 \text{ mol/L}$	Hg ²⁺ , Proteins	1) The given pH range only
Br-	Crystal	014	$c(KNO_3) = 1 \text{ mol/L}$	Hg ²⁺ , Cl ⁻ , I ⁻ , S ²⁻ , CN ⁻	applies to ion-selective
Ca ²⁺	Polymer	212	c(KCI) = 1 mol/L	Pb ²⁺ , Fe ²⁺ , Zn ²⁺ , Cu ²⁺ , Mg ²⁺	electrodes from Metrohm
Cl-	Crystal	014	$c(KNO_3) = 1 \text{ mol/L}$	Hg ²⁺ , Br ⁻ , I ⁻ , S ²⁻ , S ₂ O ₃ ²⁻ , CN ⁻	AG 2) Alternatives or more de-
CN-	Crystal	1014	c(NaOH) = 0.1 mol/L	Cl-, Br-, I-,	tailed compositions can be
Cu^{2+}	Crystal	212	$c(KNO_3) = 1 \text{ mol/L}$	Ag+, Hg ²⁺ , S ²⁻	found in the manual «Ion
F-	Crystal	57	NaCl/glacial acetic acid/ CDTA	OH-	Selective Electrodes (ISE)», order number 8.109.8042
I-	Crystal	014	$c(KNO_3) = 1 \text{ mol/L}$	Hg^{2+} , S^{2-} , $S_2O_3^{2-}$,	2) 14
K+	Polymer	2.511	c(NaCl) = 0.11 mol/L	TRIS+, NH ₄ +, Cs+, H+	3) More detailed information about interfering ions
Na+	Polymer	312	$c(CaCl_2) = 1 mol/L$	SCN ⁻ , K ⁺ , lipophilic ions	and other interferences
NH_4^+	Gas membrane	11	_	_	can be found in the
NO ₃ -	Polymer	2.511	$c((NH_4)_2SO_4) = 1 \text{ mol/L}$	Cl-, Br-, NO ₂ -, OAC-	manual «Ion Selective
Pb^{2+}	Crystal	47	$c(NaClO_4 \cdot H_2O) = 1 \text{ mol/L}$	Ag+, Hg ²⁺ , Cu ²⁺	Electrodes (ISE)», order
S ²⁻	Crystal	212	c(NaOH) = 2 mol/L	Hg ²⁺ , Proteins	number 8.109.8042

Troubleshooting

Table 8: Possible sources of interference and remedies for ion-selective electrodes

Electrode Source of inter-		Effects	Action
	ference		
Ion-selective crystal	Dissolution processes,	Rough surface → slow response,	Polish with polishing cloth
membrane	oxidation processes	poor detection limits	
	Electrode poisons	Formation of more sparingly soluble	Polish with polishing cloth,
		salts on the electrode surface than with	mask interfering ion
		the ion to be measured \rightarrow zero point	
		shift, reduced linearity range	
Ion-selective	Dissolution processes	Diffusion into the membrane or	Elimination of interfering
polymer membrane		dissolution of membrane component	components
NH₃ sensor	Volatile bases	Electrolyte becomes contaminated →	Change electrolyte
	(amines)	displacement of calibration line, limited	
		linearity	
	Surfactants	Membrane becomes wetted →	Replace membrane
		slow response	

1.4. Reference electrodes

Reference electrodes are usually electrodes of the second kind. In this type of electrode a metal electrode is in contact with a sparingly soluble salt of the same metal. The potential depends only on the solubility of the salt. As a first approximation, electrodes of the second kind do not themselves react with the solution and therefore supply a constant potential.

The most frequently used reference electrode is the silver/silver chloride electrode (Ag/AgCl/ KCl). The calomel electrode (Hg/Hg₂Cl₂/KCl), which was formerly widely used, is hardly used at all today as mercury and its salts are extremely toxic and all the applications can also be carried out with the silver/ silver chloride reference electrode. The standard hydrogen electrode SHE is also an electrode of the second kind. It is only used for calibra-

tion purposes. Some titrations offer the possibility of using pH glass electrodes as reference electrodes. Even if protons are transferred during the titration it is usually still possible to make an accurate determination of the 115

1.4.1. Silver/silver chloride reference electrode

The reference element of the silver/silver chloride reference electrode is the silver/silver chloride/potassium chloride solution system: Ag/AgCl/KCl. The reference electrode is usually filled with c(KCl) = 3 mol/L or saturated KCl solution. Tables 9 and 10 show the potentials of the reference electrode as a function of the reference electrolyte and temperature. Each of these values has been measured against the standard hydrogen electrode under isothermal conditions.

Table 9: Standard redox potentials of the silver/silver chloride reference electrode as a function of the temperature and concentration

Temp. (°C)	0	+10	+20	+25	+30	+40	+50	+60	+70	+80	+90	+95
E ⁰ (mV) with	+224.2	+217.4	+210.5	+207.0	+203.4	+196.1	+188.4	+180.3	+172.1	+163.1	+153.3	+148.1
c(KCI) = 3 mol/L												
E ⁰ (mV) with	+220.5	+211.5	+201.9	+197.0	+191.9	+181.4	+170.7	+159.8	+148.8	+137.8	+126.9	+121.5
c(KCI) = sat.												

Table 10: Standard redox potentials of the silver/silver chloride reference electrode as a function of the concentration

c(KCl) / mol/L (25 °C)	0.1	1.0	3.0	3.5	sat.
E ⁰ (mV)	+291.6	+236.3	+207.0	+203.7	+197.0

1.4.2. The Metrosensor «Long Life» reference system

Most electrodes are equipped with the silver/silver chloride reference system. The solubility product of silver chloride in water is very small (10⁻¹⁰ mol²/L²). In the concentrated, chloride-containing solution of the reference The advantages of the «Long Life» reference systems at a electrolyte soluble complexes of the series (AgCl₂)-, (AgCl₃)²⁻, (AgCl₄)³⁻ are formed. This means that the reference system poses several problems. Outside the electrode the chloride concentration is frequently lower and the complexed silver chloride precipitates in the region surrounding the diaphragm («liquid junction»). The result: precipitated silver chloride blocks the diaphragm, and the response time of the pH electrode increases. A further problem is presented by the dependency of the solubility product of AgCl on the temperature. If the electrode is used at a different temperature then the equilibrium that determines the potential of the reference electrode must be reestablished. The larger the surface with solid AgCl in As in the «Long Life» reference system the silver chloride relationship to the electrolyte volume, the shorter the time required. The «Long Life» reference system prevents high concentrations of complexed AqCl from occurring in the outer electrolyte, as the silver chloride reservoir is connected with the outer electrolyte by a highly effective diffusion barrier. The concentration of the silver complex in the reference electrolyte remains low. Even after one 1.4.3. Diaphragms

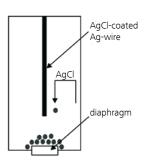


Figure 9: Conventional Ag/AgCl/ KCl system. The chloride concentration outside the electrode is usually lower than in the electrolyte chamber. The soluble silver chloride complexes precipitate out in the region surrounding the diaphragm and may block it.

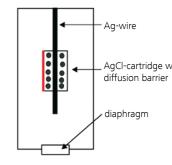


Figure 10: The Metrohm «Long Life» reference system AgCl-cartridge with The dissolved AgCl is retained in the AqCl cartridge and can no longer block the diaphragm.

year the concentration of complexated silver ions in the outer electrolyte has only reached 5% of the saturation

- Long working life of the electrode
- Rapid response to changes in pH
- Rapid response to temperature changes
- Less sensitive to electrode poisons, e.g. S²⁻

Blocking the diaphragm by crystallized AgCl also affects the electrolyte flow. If the «Long Life» reference system is used then the flow of the KCl solution through the diaphragm into deionized water only decreases slightly.

is present in a smaller volume of potassium chloride solution, the thermodynamic equilibrium between silver, silver chloride (solid) and silver chloride (dissolved) is established very quickly and the potential of the reference electrode becomes stable after a very short time.

Faulty measurements, unstable measured values and very long response times usually have their source in the «liquid junction» between the sample solution and the reference electrode. The diffusion, streaming and Donnan potentials that occur there – which are normally known together as the diaphragm potential - have various causes and can result in a very incorrect measured value.

The measuring error may assume vast proportions if measurements are made under the following conditions:

- with a blocked, virtually impermeable diaphragm,
- in ion-deficient solutions with an unsuitable
- in strong acids and bases with an unsuitable diaphragm,
- in colloidal solutions.

In all such cases errors may occur that cannot be tolerated. This is why the following questions must be in the foreground whenever an electrode and therefore the optimal type of diaphragm are to be selected:

- Does the reference electrolyte react with the sample solution to form a precipitate in the diaphragm?
- Does the electrolyte flow alter the composition of the sample solution in an unacceptable way?
- Is there a risk of depositing sample solution components on the diaphragm?
- Is the chemical resistance assured?
- Can physical parameters such as flow, pressure or temperature cause measuring errors?
- Does the process allow cleaning/maintenance of the electrode at certain intervals?
- Is a short response time and/or high reproducibility necessary?

The time required for cleaning and maintenance can usually be considerably reduced if the correct choice of electrode is made. The most frequent cause of measuring problems is contamination of the diaphragm. This is why with pH electrodes the chief attention is paid to the diaphragm during maintenance with the pH membrane being of secondary importance. If existing means cannot be used to determine whether the indicator electrode or the reference electrode requires cleaning/regeneration, then it is usually best to treat the reference electrode. Various types of diaphragm are available to satisfy the diverse requirements. These requirements have already been taken into consideration for the electrode recommendations in the application lists on pages 6 and 7.

Ceramic pin diaphragms

Ceramic pin diaphragms are frequently used diaphragms. They are primarily suitable for clear, aqueous sample solutions. They normally have pore diameters of up to 1 µm with a length and diameter each of about 1 mm. This results in an electrolyte flow rate of up to 25 µL/h, depending on the condition of the diaphragm. This means that the reference electrolyte only requires refilling at long intervals; this is why electrodes with ceramic pin diaphragms are particularly suitable for long-term measurements. On the other hand, because of their small pores and large polar surface (>>500 mm²), ceramic diaphragms tend to become blocked and therefore should not be used in solutions containing precipitates. An important advance with regard to the prevention of diaphragm blockages by silver chloride and silver sulfide has been achieved by the introduction of the «Long Life» reference system (see Section «The Metrosensor «Long Life» reference system»).

Ground-joint diaphragms with fixed or separable ground-joint

Ground-joint diaphragms with fixed or separable groundjoint are used in ion-deficient media, among others, as they produce a steady signal that is almost independent of sample flow conditions. The risk of blockage by silver chloride or by precipitates formed in the sample solution is relatively low because of the large surface area. Streaming potentials, which may occur in measurements in flowing or stirred solutions, remain negligibly small. These properties are particularly important for a SET titration to a defined pH or potential value. For example: the determination of the carbonate alkalinity by a SET titration to pH = 5.4 according to ISO 9963-2 is a widely used method in the routine analysis of drinking water. During a titration it is not possible to dispense with stirring, i.e. with an incorrectly measured pH or potential at the start of the titration an incorrect endpoint is the inevitable result. Figures 11 and 12 clearly show the difference between the Aguatrode Plus (6.0253.100), which was specially developed for this application, and a conventional pH glass electrode with ceramic pin diaphragm.

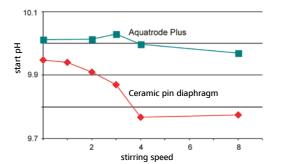


Figure 11: Measured pH of a solution with $c(Na_2CO_3) = 0.14$ mmol/L. Even under vigorous stirring the Aquatrode Plus deviates by only approx. 0.05 pH units (corresponding to approx. 3 mV) from the unstirred value, in contrast the pH glass electrode with ceramic pin diaphragm deviates by approx. 0.2 pH units.

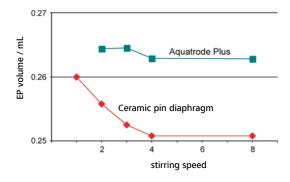


Figure 12: Endpoint volumes of a SET titration of a solution with $c(Na_2CO_3) = 0.14$ mmol/L with the titrant $c(H_2SO_4) = 0.035$ mol/L to pH 5.4. The endpoints of the Aquatrode Plus are virtually independent of the stirring speed. At higher stirring speeds the deviation from the theoretical value of the pH electrode with ceramic diaphragm amounts to approx. 5%.

Fixed ground-joint diaphragms have a uniform and reproducible electrolyte flow and are therefore particularly suitable for use with sample changers.

Separable ground-joint diaphragms are easy to clean and therefore particularly suitable for applications where contamination of the diaphragm cannot be prevented. The electrolyte flow may reach up to 100 μ L/h and is normally considerably higher than the amount of electrolyte flowing from a ceramic or fixed ground-joint diaphragm.

The ring-shaped geometry and the small polar surface of the ground-joint diaphragm have a favorable effect on the measurement. The increased electrolyte flow influences the sample solution more than if a ceramic pin diaphragm was to be used, the reference electrolyte normally needs refilling on a daily basis during long-term measurements.

An alternative is the easyClean diaphragm. It allows easy, contact-free cleaning just by pressing once on the electrode head. The spring in the electrode head returns to the defined starting position thereby ensuring greater accuracy and reproducibility of the electrolyte outflow.

Capillary diaphragms

In pH measurements in critical samples the very small pores of conventional ceramic diaphragms are easily blocked. The concept that has been realized in the Porotrode (6.0235.100), with two capillaries and a flow rate of 15...25 µL/h ensures unhindered contact between the reference electrolyte and the sample solution (liquid/ liquid phase boundary), while the two capillaries of the Porotrode are practically insensitive to contamination. The reference electrode is filled with Porolyte, which has been specially developed for this electrode. The constant flow of Porolyte ensures that the potential is established quickly and reproducibly. The flow rate and therefore the refilling intervals are comparable to those of conventional electrodes. Extra maintenance work is not necessary. Measurements in problematic samples can be carried out easily and reproducibly thanks to the concept that has been realized in the Porotrode. The pH of samples containing protein, such as milk and beer, can now be determined without any diaphragm problems. In contrast to traditional pH electrodes the Porotrode measures correctly even at high surfactant concentrations.

Twin pore

Measuring the pH in semi-solid samples such as cheese, meat and fruit places special demands on an electrode. Proteins, fats and carbohydrates and other semi-solid substances in foodstuffs tend to block the fine pores of the ceramic diaphragms used in most pH electrodes, as such substances adhere extremely well to the fine-pore ceramic surface. With the development of the spearhead electrode (6.0226.100) and the polymer electrolytes this problem has been elegantly eliminated: two pinhole diaphragms take over the function of the «liquid junction» between the sample and the reference electrode. The polymer electrolyte adjacent to the openings, which is spiked with potassium chloride and thickened, is to a large extent insensitive to contamination by media containing proteins and fats. This insensitivity to contamination, the efficient protection of the reference electrode against the penetration of electrode poisons and the optimized inner buffer of the measuring electrode ensure that the new spearhead electrode has an outstanding

long-term behavior: even when used in difficult media the electrode zero point retains its long-term stability. The use of polymer electrolytes means that refilling a liquid reference electrolyte is no longer necessary.

The new Ecotrode Gel electrodes (6.0221.x00) are equipped with this diaphragm which keeps maintenance effort low

Plied platinum wire

In combination with the reference electrolyte Idrolyte, which contains glycerol, the plied-platinum-wire diaphragm is outstandingly suitable for applications in biological media. The precipitation of proteins is suppressed by using an electrolyte with a low KCl content. The multi-capillary system (channels between the platinum wires) reduces contamination effects and the electrically conductive platinum reduces the response time and the diaphragm resistance. However, cross-sensitivity may occur in strongly redox-buffered solutions.

Cleaning and care of diaphragms

Table 11: Recommended ways of cleaning diaphragms

Type of diaphragm	Type of contamination	Cleaning
General	Preventiv and regular care	pHit kit (6.2325.000) according to instructions
	Precipitates of silver halides and silver sulfides	Immerse diaphragm for several hours in a solution of 7% thiourea in 0.1 mol/L HCl.
	Proteins, polypeptides	Immerse diaphragm for several hours in a solution of 5% pepsin in 0.1 mol/L HCl.
	Suspensions, solids, resins, glues, oils, fats	Clean electrode with suitable solvent
Fixed ground-joint	All types of contamination	Aspirate off reference electrolyte and immerse electrode in the corresponding cleaning solution.
Separable ground-joint	All types of contamination	Loosen the ground-joint sleeve (using hot water if necessary) and clean according to the type of contamination.
Capillary	Electrolyte flow interrupted	Apply slight counterpressure to electrolyte refilling opening

1.4.4. Reference electrolytes and bridge electrolytes

The reference or bridge electrolyte is in electrical contact with the sample solution via the diaphragm. The sample solution and electrolyte form a phase boundary with different ion concentrations on each side. This difference in concentration causes diffusion of the ions to the other side and, because of the different ion mobilities, a so-called diffusion potential occurs. In order to achieve a high degree of measuring accuracy the electrolyte composition must be selected so that any diffusion potentials formed are as negligible as possible; this is to a large

extent achieved by the use of c(KCI) = 3 mol/L. On the one hand the ionic mobilities of K^+ and CI^- are practically the same, on the other hand the ionic concentration in the sample solution is normally negligibly low in comparison to c(KCI) = 3 mol/L. This is why the equal-transference KCI electrolyte is used as standard in all combined Metrohm electrodes and reference electrodes. However, certain media require the use of other electrolyte compositions in order to suppress effects that occur in addition to the diffusion potential.

Table 12: Alternatives to the standard reference electrolyte c(KCl) = 3 mol/L

Medium	Problems with standard electrolytes c(KCl) = 3 mol/L	Alternative electrolyte
Silver ions	Reaction with Cl⁻ with precipitation of AgCl → slow response	$c(KNO_3) = 1 \text{ mol/L (or Titrode for more or less constant pH value)}$
Non-aqueous	Precipitation of KCl, solutions and electrolyte immiscible → unsteady signal	c(LiCl) = 2 mol/L in ethanol or LiCl saturated in ethanol
Ion-deficient water	Contamination of the medium by salt → drift	KCl solution of lower concentration
Proteins/polypeptides	Precipitation of the proteins with KCl and AgCl → zero point shift/reduced slope	Idrolyte ¹
Semi-solid substances	Contamination of diaphragm → zero point shift/slow response	Solid electrolyte in combination with pinhole diaphragm
Surfactants (proteins)	Adsorption on diaphragm → zero point shift/ reduced slope	Porolyte ²

¹ Idrolyte is a glycerol-based electrolyte whose chloride ion activity corresponds to that of a KCl solution with c(KCl) = 3 mol/L. This means that the latter can also be readily replaced by Idrolyte. Idrolyte is excellent for use with solutions containing proteins and aqueous solutions with an organic fraction.

Table 13: Electrolyte flow rates and viscosities

Electrolyte	Viscosity	Flow rate µL/h (10 c	low rate μL/h (10 cm water column)						
	(25 °C) (mPas)	Ceramic pin	Flexible ground- joint	Fixed ground- joint	Ceramic capillary	Plied Pt wire			
c(KCl) =	~1	Standard electrode	Ø 10 mm: 20100	530	_	_			
3 mol/L		525							
		Microelectrode 515	Ø 5 mm: 530						
$c(KNO_3) =$	~1	1025	Ø 10 mm: 20100	-	-	_			
1 mol/L			Ø 5 mm: 530						
Idrolyte	810	_	_	_	_	325			
Porolyte	12001500	_	_	_	530	_			

2. Fundamentals of conductometry

2.1. General

Conductometry means measuring the conductivity – a conductometer measures the electrical conductivity of ionic solutions. This is done by applying an electric field between two electrodes. The ions wander in this field. The anions migrate to the anode and the cations to the cathode. In order to avoid substance conversions and the formation of diffusion layers at the electrodes (polarization), work is carried out with alternating voltage. The rule of thumb is that the frequency of the alternating voltage must be increased as the ion concentration increases. Modern conductometers automatically adapt the measuring frequency to the particular measuring conditions.

Ion migration in an electric field depends on many factors. The temperature has a decisive influence on the viscosity of the solution and therefore on the mobility of the ions. As the temperature increases the viscosity decreases and the conductivity increases. Dissociation constants are also temperature-dependent quantities. This is why it is important to make measurements at a constant temperature or to compensate for changes of temperature by using the so-called temperature coefficient. The temperature coefficient of most salt solutions is approx. 2%/°C, but depends on the temperature in very dilute solutions.

The measuring unit used in conductivity measurements is the electrical resistance of the solution. This means that the conductivity is a sum parameter which includes all dissolved ions. Conductivity cannot be used for the determination of a single type of ion, unless the sample is a solution of a single salt or the concentrations of the other ions are known. The reciprocal value of the measured resistance of the solution, the so-called conductance L with the unit Siemens (S = $\Omega^{\text{-1}}$) is by itself less meaningful, as the shape of the measuring cell must be taken into account. The cell constant c of a conductometric measuring cell

$c = \frac{distance\ between\ Pt\ sheets}{electrode\ surface\ area}[cm^{-l}]$

Cell constant (10)

must be known. The result of the measurement is therefore always given as the specific conductivity K with the unit Siemens per cm ($S \cdot cm^{-1}$).

$$\kappa = L * c /S cm^{-1}$$

Specific conductivity (11)

This means that the conductometer must be calibrated before each measurement by determining the cell constant in a solution of known specific conductivity. The specific conductivity for various concentrations of many salts is given in tables. The specific conductivity K is linked with the concentration ci of the individual ion i via the concentration-dependent equivalent conductivity Λi . The equivalent conductivity Λi is similar to the activity coefficient γ (see Section 1.2.) and is also a quantity that depends on the concentration.

$$\kappa = \sum_{i} (\Lambda_i * z_i * c_i)$$

Specific conductivity and concentration (12)

At great dilutions, i.e. $ci \le 0.001$ mol/L, the equivalent conductivity Λi can be equated with the equivalent conductivity shown in the tables for an infinite dilution.

² Porolyte is a KCl solution that has been gelled by polymerization and is used in electrodes with a capillary diaphragm (Porotrode).

Table 14: Conductivity κ of various substances and solutions

Conductor	T (K)	Conductivity due to	Conductivity κ (μS cm ⁻¹)
Metallic copper	273	Electron conduction	645,000,000,000
Potassium hydroxide	291	Ionic conduction resulting from the complete	184,000
solution (c = 1 mol/L)		dissociation of KOH	
KCl solution	293	Ionic conduction resulting from the complete	11,660
(c = 0.1 mol/L)		dissociation of KCl	
Brackish water	273	lonic conduction resulting from the dissociation of salts	20,000 to 1,000,000
		and carbonic acid	
Acetic acid	291	lonic conduction resulting from the partial dissociation	1300
(c = 1 mol/L)		of CH₃COOH	
Drinking water	298	lonic conduction resulting from the dissociation of salts	10 to 2000
		and carbonic acid	
Graphite	273	Electronic conduction	1200
Distilled water	273	Ionic conduction resulting from contamination	0.0610
		by salts, dissociation of water and carbonic acid	
Ultrapure water	273	Ionic conduction resulting from low self-dissociation	0.056
Pure benzene	273	Ionic conduction resulting from the dissociation of	0.00000005
		traces of water	

Conductometry is used for direct measurements and in titration. The theory is identical for both methods. Whereas in direct measurements it is the absolute value that is of interest, in titrations it is the change in the measured value. Direct measurement is often used for monitoring surface waters, waterworks, water desalination plants and in the preparation of ultrapure water, where particular limits must not be exceeded. Conductivity detection is mostly used for precipitation titrations, where the equivalent point is recognized by the conductivity reaching a minimum value. The absolute value is of secondary importance.

Conductivity measurements

Whereas the instruments used for potentiometry have been standardized (input impedance >1012 Ω , zero point at pH 7), this is not the case with conductometers. The influence of the cable capacity, the measuring frequency level, the conductivity range and the adjustable cell constant, the method used for conductivity measurements (phase-sensitive, frequency-dependent, bipolar pulse, etc.) vary and depend on the type of instrument. This means that the instrument must be taken into account for solving application problems. Important parameters are:

- Platinizing quality (platinum black) → high series capacity C_S
- Electrode area A \rightarrow high series capacity C_S
- Cell constant c
- Measuring frequency f
- Cable capacity C_P
- Cable resistance R_C
- Instrument measuring range (resistance range)

Selecting the right cell constant

The cell constant c is defined for conductometric measuring cells. A measuring cell with two parallel electrodes at a distance of 1 cm and each with an area of 1 cm² theoretically has a cell constant $c = I \cdot A^{-1} = 1$ cm³. The cell constant is never exactly $I \cdot A^{-1}$, as the electric field is not strictly homogeneous. The rules of thumb given in Table 15 are used for selecting the correct measuring cell:

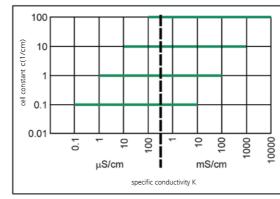


Figure 13: Cell constants and recommended conductivity intervals.

Interferences, care

Conductivity measuring cells with Pt sheets

Conductivity cells have a very porous black platinum coating in order to avoid polarization effects in media with a high conductivity. However, the properties of this coating may change in time (contamination, abrasion of the platinum coating, etc.); this could alter the cell constant. This is why it is absolutely necessary to calibrate the conductometric measuring cell before making a measurement in order to avoid measuring errors. For the exact determination of cells with a cell constant of <1 cm⁻¹ a solution with a conductivity of about 100 mS/ cm is recommended. If measurements are made in well conducting media then a check of the activity of the platinum layer is additionally recommended, e.g. in a 0.1 mol/L KCl solution. If a lower specific conductivity is shown then cleaning with a suitable oxidizing agent or solvent is indicated.

If measurements are made in ion-deficient water then frequent calibration is unnecessary, as in this case the activity (series capacity) of the platinum layer is not very important and the deposition of highly isolating substances is not to be expected. The measuring cell must be thoroughly cleaned after calibration in order to avoid incorrect measuring results caused by adherent KCl solution.

Table 15: Recommended cell constants

Cell constant	Sample
$c = 0.1 \text{ cm}^{-1}$	For very poorly conducting solutions such as distilled water, deionized or partly deionized water, etc. For applications according to USP 645 and EP 2.2.38
$c = 1 \text{ cm}^{-1}$	For moderately conducting solutions such as drinking water, surface water, wastewater, etc.
$c = 10 \text{ cm}^{-1}$	For solutions with good conductivity such as seawater, rinsing water, physiological solutions, etc.
$c = 100 \text{ cm}^{-1}$	For solutions with very good conductivity such as electroplating baths, salt solutions, etc.

Conductivity measuring cells (stainless steel) 124

tamination or corrosion. However, also these measuring cells should be calibrated before the measurements in

These are usually substantially more insensitive to con-order to achieve the highest possible measuring accuracy. For cleaning, water and/or ethanol alone should be used.

Table 16: Conductivity measurement – interferences

Source of	Effects	Measures
interference		
Low conductivity	Values too high, drift	Drive off atmospheric CO ₂ with inert gas
with open vessel		(Ar, N_2) or use flow-through cell, avoid carryover of salt solutions (e.g. too frequent calibration,
		inadequate rinsing)
Oils, precipitation	Isolating layer on electrode →	Clean with solvent or oxidizing agent
products	cell constant increases, measuring range	
	limited to higher values	
Unstable	Unstable values	Temperature compensation, if temperature
temperature		coefficient is known, or thermostatting
		(temperature coefficient generally approx. 2%/°C)
Conductance	Stray fields outside electrode shaft	Watch distance from vessel during
depends on	(particularly for cells with constants >1 cm ⁻¹)	calibration and measurement or select
electrode position	→ measured value displaced	cell constant 1 cm ⁻¹
		Use of 5-ring conductivity measuring cells
Foreign salts	Carryover of residual salts when changing	Thorough previous rinsing of electrode
	to solutions with low conductivity → drift	
	to higher values	
Air bubbles	Air bubble located between	Remove air bubble by tapping
	electrode plates → unsteady signal	

2.2. Conductivity measurement in accordance with USP and Pharm. Europe (EP)

There are special requirements for the conductivity measlatest EP -4.8-07/2004:0169. Besides a precision conduc- following three conditions is met: tometer whose temperature compensation can be disabled, a conductivity measuring cell and a conductivity

standard are to be used that allow to determine the cell constant with a maximum measuring error of 2%. To prevent the uptake of carbon dioxide, the measurement urement in water for pharmaceutical use («water for should be carried under exclusion of air and/or in a flow injections») according to USP 645, EP 2.2.38 resp. the cell. The sample fulfills the specification if one of the

Stage 1:

The sample is measured directly without further pretreatment and without temperature compensation. If the test is considered as passed.

Stage 2:

If the conditions are not fulfilled by stage 1, continue as cannot be used for pharmaceutical purposes. follows: The conductivity of at least 100 mL sample is measured under strong agitation at 25 °C ±1 °C as soon as the drift caused by the uptake of carbon dioxide is smaller than 0.1 μ S/cm per five minutes. If the measured value is smaller than 2.1 μ S/cm, the test is considered as passed.

Stage 3:

However, if the sample does not fulfill the specification of stage 2, a sample of exactly 100 mL is mixed with water fulfills the specification indicated in table 17, the 0.3 mL saturated KCI solution. Then the pH value of this solution is measured exactly to 0.1 pH units. Only if the conductivity fulfills the conditions specified in table 18 the test is considered as passed. Otherwise the water

Table 17: First step of the conductivity measurement according to USP 645 and EP -4.8-07/2004:0169

Temperature	Conductivity	Temperature	Conductivity
	not larger than (µS/cm)		not larger than (µS/cm)
0	0.6	55	2.1
5	0.8	60	2.2
10	0.9	65	2.4
15	1.0	70	2.5
20	1.1	75	2.7
25	1.3	80	2.7
30	1.4	85	2.7
35	1.5	90	2.7
40	1.7	95	2.9
45	1.8	100	3.1
50	1.9		

Table 18: pH and conductivity criteria for stage 3

pH	Conductivity	рН	Conductivity
	not larger than (µS/cm)		not larger than (µS/cm)
5.0	4.7	6.1	2.4
5.1	4.1	6.2	2.5
5.2	3.6	6.3	2.4
5.3	3.3	6.4	2.3
5.4	3.0	6.5	2.2
5.5	2.8	6.6	2.1
5.6	2.6	6.7	2.6
5.7	2.5	6.8	3.1
5.8	2.4	6.9	3.8
5.9	2.4	7.0	4.6
6.0	2.4		

3. Temperature measurement

126

Only a few of the electrodes offered are fitted with a pensation. Under such circumstances it is possible to built-in temperature sensor. The decision whether temperature measurement/compensation is necessary depends on the required accuracy. Differing diffusion However, if a high degree of reproducibility of the measglass can result in measuring errors that are far in excess absolutely necessary. of the errors caused by the absence of temperature com-

potentials, e.g. in highly concentrated or very dilute solu- ured values is demanded or if GLP requirements have to tions, or changes to the diaphragm or the membrane be met then temperature measurement/compensation is

Table 19: Temperature measurement or temperature compensation: yes or no?

Measurement requirements	Temperature compensation or measurement
Are GLP requirements to be met?	Yes: temperature measurement
Is high measuring accuracy required?	Yes: temperature compensation (see Nernst potential)
Direct measurement?	Yes: temperature compensation (see Nernst potential)
Titration?	No: relative measurement
Is the pH of the sample about 7?	Yes: temperature compensation not absolutely necessary
	(low influence, as electrode zero point is at $pH = 7$),
	possibly temperature measurement
Does the pH value differ greatly from pH 7?	Yes: temperature compensation (see Nernst potential)
Are measurements made at different temperatures?	Yes: temperature compensation/measurement
	(see Nernst potential)
Is the pH of the sample solution very temperature-	Yes: temperature measurement (measurement
dependent?	temperature must be mentioned)
Does the application require a different type of	Yes: separate temperature sensor (electrodes with
diaphragm than a ceramic pin?	built-in temperature sensors are only available with
	ceramic pin and fixed ground-joint)
Is the working life of the electrode very short?	Yes: for cost reduction use separate temperature sensor

	1				1			1							1			I		1	
6.01 - 6.02 pH glass electrodes	Max. installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Temperature sensor	Temperature range short-term (°C)	Temperature range long-term (°C)	Shape	pH range	Membrane glass	Membrane resistance (M요)	Electrode zero point (mV)	Electrode slope	Asymmetry potential (mV)	Diaphragm	Reference electrolyte	Electrolyte outflow (µL/h)	Reference system	Max. Reference resistance (kΩ)
6.A2:W360	142	12	12	15	Glass	G		080	080	Sphere	014	Т	40150	±15	>0.97	±15		 <u> </u>	_	_	_
6.0233.100	101	12				G			080		014	T	150400	±15	>0.97	±15	Coromic	c(KCl)=3 mol/L	1025	II system	5
6.0253.100	135	12	12 12	20	Glass	G		080	060	Hemisphere Sphere	014	A	80200	±15	>0.97	±15	Ceramic Fixed ground-joint	c(KCI)=3 mol/L c(KCI)=3 mol/L (gel)	530	LL system	50
6.0257.600	135	12	12	20	Glass	U U	Pt1000	060	060	Sphere	013		80200	±15	>0.97	±15		_		LL system	50
6.00257.000	135	12	12	20	Glass	Fixed cable with plug F		060	060	Sphere	013	A	80200	±15	>0.97	±15	Fixed ground-joint Fixed ground-joint	c(KCL)=3 mol/L (gel) c(KCl)=3 mol/L (gel)	530	LL system	50
6.0257.020	260	12	12	20	Glass	Fixed cable with plug F		060	060	Sphere	013	A	80200	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L (gel)	530	LL system	50
6.00202.300	135					, ,		060	060	Sphere	013	A	80200	±15	>0.97		Fixed ground-joint	_	530	LL system	
6.00202.300		12	12	20	Glass Glass	Q			060	Sphere	013		80200	±15	>0.97	±15	,	c(KCI)=3 mol/L (gel)		LL system	50
6.0277.300	135 113		3	7	Glass	G	Pt1000	060			111	A	300600	±15	>0.97	±15	Fixed ground-joint Platinum wire	c(KCI) = 3 mol/L	530	LL system	50
		12							060	Hemisphere		Spec.				±15		Idrolyt		LL system	30
6.0262.100	125	12	12	20	Glass	G G		080	080	Hemisphere	013	E	150400	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	20
6.0221.100	125	12	12	20	Glass	-	D+1000	060	040	Hemisphere	111	E	<400	±20	>0.97	±15	Twin pore	Gel	-	LL system	20
6.00221.600	125	12	12	20	Glass	U	Pt1000	060	040	Hemisphere	111		<400	±20	>0.97	±15	Twin pore	Gel	- - -	LL system	20
6.00201.300	125	12	12	20	Glass	Q		080	080	Hemisphere	013	E	150400 <400	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0280.300	125	12	12	20	Glass	K C		080	080	Hemisphere	013	E		±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	10
6.0269.100	125	12	12	20	Glass	G		080	080	Sphere	013	A	80200	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	10
6.0256.100	125	12	12	20	Glass	G		060	060	Flat membrane	013	Spec.	<2000	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0234.100	113	12	6.4	20	Glass	G		080	080	Hemisphere	014	T	200500	±15	>0.97	±15	Ceramic	c(KCl)=3 mol/L	525	LL system	5
6.0234.110	168	12	6.4	20	Glass	G		080	080	Hemisphere	014	T	200500	±15	>0.97	±15	Ceramic	c(KCl)=3 mol/L	525	LL system	5
6.0235.200	125	12	12	20	Glass	G		080	080	Hemisphere	014	T	200400	±15	>0.97	±15	Two capillaries (ceramic)	Porolyt	530	LL system	15
6.0255.100	113	12	12	30	Glass	G		0100	080	Cylinder	014	U	150500	±15	>0.97	±15	Ground-joint	c(KCl)=3 mol/L	5100	LL system	5
6.0255.110	170	12	12	30	Glass	G		0100	080	Cylinder	014	U	150500	±15	>0.97	±15	Ground-joint	c(KCl)=3 mol/L	5100	LL system	5
6.0255.120	310	12	12	30	Glass	G		0100	080	Cylinder	014	U	150500	±15	>0.97	±15	Ground-joint	c(KCl)=3 mol/L	5100	LL system	5
6.00204.300	125	12	12	30	Glass	Q		0100	070	Cylinder	014	U	150500	±15	>0.97	±15	Ground-joint	c(KCl)=3 mol/L	5100	LL system	5
6.0220.100	113	12	12	15	PP	G		080	080	Hemisphere	014	T	200500	±15	>0.97	±15	Ceramic	c(KCl)=3 mol/L	310	LL system	5
6.0223.100	113	12	12	15	PP	G	D:4000 (2	040	040	Hemisphere	112	Spec.	100650	±15	>0.97	±15	Ceramic	c(KCI)=3 mol/L	310	LL system	5
6.0228.000	113	12	12	15	PP	Fixed cable with plug F	Pt1000 (2 mm)	080	080	Hemisphere	014	T	200400	±15	>0.97	±15	Ceramic	c(KCl)=3 mol/L	310	LL system	5
6.0228.030	113	12	12	15	PP	Fixed cable with plug I (IP67)	Pt1000 (2 mm)	080	080	Hemisphere	014	T	200400	±15	>0.97	±15	Ceramic	c(KCl)=3 mol/L		LL system	5
6.0229.010	125	12	12	20	Glass	Fixed cable with plug F		070		Sphere	014	T	40150	0100	>0.90	±15	EasyClean	LiCl/EtOH		LL system	250
6.0229.020	125	12	12	20	Glass	Fixed cable with plug F		070	070		014	T	40150	0100	>0.90	±15	EasyClean	LiCI/EtOH		LL system	250
6.0229.100	113	12	12	30	Glass	G		070	0 70		014	T	40150	0100	>0.90	±15	Ground-joint	LiCI/EtOH	350	LL system	250
6.00203.300	125	12	12	30	Glass	Q		070		Sphere	014	Т	40150	0100	>0.90	±15	Ground-joint	LiCI/EtOH	350	LL system	250
6.0279.300	113	12	12	30	Glass	K		070		Sphere	014	T	40150	0100	>0.90	±15	Ground-joint	LiCl/EtOH	350	LL system	250
6.0226.100	98	12	6	10	Glass	G		060		Needle	111	Spec.	200500	±20	>0.97	±15	Twin pore	Gel	-	LL system	20
6.00226.600	98	12	6.4	10	Glass			060	040	Needle	111	Spec.	200500	±20	>0.97	±15	Twin pore	Gel	F 30	LL system	20
6.0248.600	288	12	12	25	Glass		Pt1000	0100	080	Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.00249.600	438	12	12	25	Glass	U	Pt1000	0100	080	Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.00249.540	112	12	12	25	Glass	S		0100		Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0259.100	113	12	12	25	Glass	G		0100		Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0258.600	125	12	12	25	Glass		Pt1000	0100	1	Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L		LL system	5
6.0258.010	125	12	12	25	Glass	Fixed cable with plug F		0100		Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L		LL system	5
6.0260.010	125	12	12	20	Glass	Fixed cable with plug F			0100		014	U	150600	±15	>0.97	±15	EasyClean	c(KCl)=3 mol/L		LL system	5
6.0260.020	125	12	12	20	Glass	Fixed cable with plug F	Pt1000			·	014	U	150600	±15	>0.97	±15	EasyClean	c(KCl)=3 mol/L	350	LL system	5
6.00200.300	125	12	12	20	Glass	Q		0100		Hemisphere	014	U	150600	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0278.300	125	12	12	25	Glass	K		0100		Cylinder	014	U	150500	±15	>0.97	±15	Fixed ground-joint	c(KCl)=3 mol/L	530	LL system	5
6.0239.100	113	12	12	30	Glass	G		080	080	Hemisphere	014	T	150400	±15	>0.97	±15	Ground-joint	c(KCl)=3 mol/L	5100	LL system	5

6.03 Separate metal electrodes	Max. installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Туре	Temperature range short-term (°C)	Temperature range long-term (°C)	Shape	Measuring range	pH range
6.0301.100	125	12	12	10	Glass	В	Pt	-2070	-2070	Wire	-20002000 mV	014
6.0309.100	101	12	12	10	Glass	G	Pt	-2070	-2070	Sheet	-20002000 mV	014
6.0338.100	96	8	8	5	Glass	G	Pt	-2070	-2070	Wire	-20002000 mV	014
6.0340.000	103	5.3	5.3	10	Glass	Fixed cable / Plug F	Pt	-2070	-2070	Wire	-20002000 mV	014
6.0341.100	101	12	8.8	10	Glass	G	Pt	-2070	-2070	Wire	-20002000 mV	014
6.0343.100	86	8	8	10	Plastic	Plug-in connector 5.5 mm	Pt	080	080	Pin	-20002000 mV	014
6.0331.000	100	NS14/15	3	5	PP	Socket B negative with cable (I = 1 m)	Pt	080	080	Pin	-20002000 mV	014
6.0331.010	100	NS14/15	3	5	PP	Socket B negative with cable (I = 1 m)	Ag	080	080	Pin	-20002000 mV	014
6.0331.010S	100	NS14/15	3	5	PP	Socket B negative with cable (I = 1 m)	Ag	080	080	Pin	-20002000 mV	014
6.0344.100	108	24	24	15	Glass	G	Pt	-2070	-2070	Grid/Sheet		014
6.0345.100	108	24	24	15	Glass	G	Pt	-2070	-2070	Grid/Sheet		014
6.00346.100	96	8	3	7	Glass	G	Pt	-2070	-2070	Ring	-20002000 mV	014
6.00353.100	125	12	6	20	Glass	G	Au	-2080	-2080	Ring	-20002000mV	014
6.0351.100	125	12	12	7	Glass	G	Pt	-2080	-2080	Ring	-20002000 mV	0 14
6.0352.100	125	12	12	7	Glass	G	Au	-2080	-2080	Ring	-20002000 mV	0 14
6.00350.100	125	12	6	7	Glass	G	Ag	-2080	-2080	Ring	-20002000 mV	014
6.00350.100S	125	12	6	7	Glass	G	Ag	-2080	-2080	Ring	-20002000 mV	014
6.0343.110	86	8	8	10	Stainless steel	Plug connector 5.5 mm	Steel	080	080			
6.0346.000	86	12	1.5	5	Glass	Plug connector 5.5 mm		060	060			

6.04 Combined metal electrodes	Max. installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Туре	Temperature range short-term (°C)	lemperature range long-term (°C)	Shape	Measuring range	pH range	Diaphragm	Reference electrolyte	Electrolyte outflow (µL/h)	Reference system	Reference resistance (kΩ)	Reference resistance (kΩ)
					S Noryl/					N	-20002000			c(KCl)=				
6.0421.100	113	12	12	10	PP	G	Sb rod	070	070		mV -20002000	211	Ceramic	3 mol/L	310	LL system	5	5
6.00430.100	125	12	6	20	Glass	G	Ag	080	080	Cylinder	mV -20002000	014				pH	150400	150400
6.00430.100BR	125	12	6	20	Glass	G	Ag	080	080	Cylinder	mV	014				pH	150400	150400
6.00430.100CL	125	12	6	20	Glass	G	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	200500
6.00430.100S	125	12	6	20	Glass	G	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	200500
6.00404.300	125	12	6	20	Glass	Q	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	200500
6.00404.300S	125	12	6	20	Glass	Q	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	5
6.00470.300	125	12	6	20	Glass	K	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	5
6.00470.300S	125	12	6	20	Glass	K	Ag	080	080	Cylinder	-20002000 mV	014				рН	150400	5
6.0431.100	125	12	12	20	Glass	G	Pt	080	080	Ring/ hemisphere	-20002000 mV	014				рН	150400	5
6.00401.300	125	12	12	20	Glass	Q	Pt	080	080	Ring/ hemisphere	-20002000 mV	014				рН	150400	5
6.0471.300	125	12	12	20	Glass	K	Pt	080	080	Ring/ hemisphere	-20002000 mV	014				рН	150400	150400
6.0433.110	178	12	6.4	20	Glass	G	Ag	080	080	Ring/ hemisphere	-20002000 mV	014				рН	200500	200500
6.0434.110	178	12	6.4	20	Glass	G	Pt	080	080	Ring/ hemisphere	-20002000 mV	014				рН	200500	150400
6.0435.110	178	12	6.4	20	Glass	G	Au	080	080	Ring/ hemisphere	-20002000 mV	014				рН	200500	< 5 kΩ
6.00435.120	308	12	6.4	20	Glass	G	Au	080	080	Ring/	-20002000 mV	014				рН	200500	< 5 kΩ
6.00450.100	125	12	6	15	Glass	G	Ag	080	080	hemisphere Ring	-20002000 mV	014	Fixed	c(KNO ₃)= 1 mol/L	1025	LL system	5	150400
6.00450.100BR	125	12	6	15	Glass	G	Ag	070	070	Ring	-20002000	014	ground-joint Fixed	c(KNO ₃)=	1025	LL system	5	200500
6.00450.100CL	125	12	6	15	Glass	G	Ag	070	070	Ring	mV -20002000	014	ground-joint Fixed	1 mol/L c(KNO ₃)=	1025	LL system	5	
6.00450.100S	125	12	6	15	Glass	G	Ag	070	070	Ring	mV -20002000	014	ground-joint Fixed	1 mol/L c(KNO ₃)=	1025	-	5	
6.00402.300	125	12	12	20	Glass	Q	Ag	080	080	Ring	mV -20002000	014	ground-joint Fixed	1 mol/L c(KNO3)=	330	,	5	
6.00402.3005	125	12	6	20	Glass	Q	Ag	070	070	Ring	mV -20002000	014	ground-joint Fixed	1 mol/L c(KNO ₃)=	330	LL system		
										_	mV -20002000		ground-joint Fixed	1 mol/L c(KNO ₃)=				
6.00450.300	125	12	6	15	Glass	K	Ag	080	080	Ring	mV -20002000	014	ground-joint	1 mol/L c(KCl)=	1025	LL system	5	
6.0451.100	113	12	12	15	Glass	G	Pt	-580	-580	Ring	mV -20002000	014	Ceramic	3 mol/L c(KCl)=	1025	LL system		
6.00403.300	125	12	12	20	Glass	Q	Pt	-580	-580	Ring	mV -20002000	014	ground-joint	3 mol/L c(KCl)=	330		5	
6.0451.300	113	12	12	15	Glass	K	Pt	-580	-580	Ring	mV	014	Ceramic	3 mol/L	1025	,	5	
6.0452.100	113	12	12	15	Glass	G	Au	-580	-580	Ring	-20002000 mV	014	Ceramic	c(KCl)= 3 mol/L	1025	LL system	5	

6.05 Ion-selective electrodes	Max. installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Туре	Temperature range short-term (°C)	Temperature range long-term (°⊂)	Shape	Measuring range	pH range
6.0502.100	125	12	12	1	EP	G	Crystal (Br)	050	050	Flat	1x10 ⁻⁶ 1 mol/L	014
6.0502.120	125	12	12	1	EP	G	Crystal (CI)	050	050	Flat	1x10 ⁻⁵ 1 mol/L	014
6.0502.130	125	12	12	1	EP	G	Crystal (CN)	080	080	Flat	8x10 ⁻⁶ 10 ⁻² mol/L	1014
6.0502.140	125	12	12	1	EP	G	Crystal (Cu)	080	080	Flat	1x10 ⁻⁸ 0.1 mol/L	212
6.0502.150	125	12	12	1	EP	G	Crystal (F)	080	080	Flat	1x10 ⁻⁶ sat. mol/L	57
6.00500.600	125	12	6	20	Glass	U	Crystal (F)	050	040	Flat	1x10 ⁻⁶ sat. mol/L	57
6.00500.300	125	12	6	20	Glass	Q	Crystal (F)	050	040	Flat	1x10 ⁻⁶ sat. mol/L	57
6.0502.160	125	12	12	1	EP	G	Crystal (I)	050	050	Flat	5x10 ⁻⁸ 1 mol/L	014
6.0502.170	125	12	12	1	EP	G	Crystal (Pb)	080	080	Flat	1x10-60.1 mol/L	47
6.0502.180	125	12	12	1	EP	G	Crystal (Ag/S)	080	080	Flat	1x10 ⁻⁷ 1 mol/L	212
6.0506.100	125	12	12	5	PEEK/POM	G	NH ₃ -permeable membrane	050	050	Flat	5x10 ⁻⁶ 10 ⁻² mol/L	11
6.0506.150	125	12	12	5	PEEK/POM	G	NH₃-permeable membrane	050	050	Flat	10 ⁻⁴ 1 mol/L	11
6.0507.010	125	12	2.5	20	PVC	G	Polymer (Non- ionic surfactants)	040	040	Pin	surfactant- dependent	012
6.0507.120	125	12	2.5	20	PVC	G	Polymer (Ionic surfactants)	040	040	Pin	surfactant- dependent	012
6.0507.130	125	12	12	5	POM	G	Graphit (Ionic surfactants)	1050	1050	Flat	surfactant- dependent	010
6.0507.140	125	12	12	1	PEEK	G	Graphit (Ionic surfactants)	040	040	Flat	surfactant- dependent	013
6.0507.150	125	12	2.5	20	PVC	G	Polymer (Ionic surfactants)	040	040	Pin	surfactant- dependent	012
6.0501.100	100	12	12	20	Glass	G	Glass (Na)	080	080	sphere	5x10 ⁻⁵ 1 mol/L	59
6.0508.100	125	12	12	1	PVC	G	Polymer (Na)	040	040	Flat	5x10 ⁻⁷ 1 mol/L	312
6.0508.110	125	12	12	1	PVC	G	Polymer (Ca)	040	040	Flat	5x10 ⁻⁶ 1 mol/L	212
6.0510.100	113	12	12	10	PMMA/PP	G	Polymer (Ca)	040	040	Flat	5x10 ⁻⁷ 1 mol/L	212
6.00502.300	113	12	12	10	PMMA/PP	Q	Polymer (Ca)	040	040	Flat	5x10 ⁻⁷ 1 mol/L	212
6.0510.110	113	12	12	10	PMMA/PP	G	Polymer (K)	040	040	Flat	1x10 ⁻⁷ 1 mol/L	2.511
6.00510.120	113	12	12	10	PMMA/PP	G	Polymer (NO ₃)	040	040	Flat	1x10 ⁻⁶ 1 mol/L	2.511

6.07 Reference electrodes	Max. installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Temperature range short-term (°C)	Temperature range long-term (°C)	Diaphragm	Reference electrolyte	Bridge electrolyte	Outflow (microL/h) in relation to 3 mol/L KCl	Reference system	Max Reference resistance (kOhm) in relation to 3 mol/L KCl
6.0724.140	43	12	8	20	Glass	В	080	080		variable		0		1
6.0726.100	100	12	12	10	Glass	В	080	080	Ground-joint	variable	variable	550	Ag wire/AgCl	3
6.0726.107	100	12	12	10	Glass	В	080	080	Ground-joint	c(KCl)=3 mol/L	c(KCl)=3 mol/L	550	Ag wire/AgCl	3
6.0726.110	138	12	8	10	Glass	В	080	080	Ground-joint	variable	variable	550	Ag wire/AgCl	3
6.0729.100	100	12	12	10	Glass	G	080	080	Ground-joint	variable	variable	550	Ag wire/AgCl	3
6.0729.110	138	12	12	10	Glass	G	080	080	Ground-joint	variable	variable	550	Ag wire/AgCl	3
6.0736.110	178	12	6.4	10	Glass	В	080	080	Ground-joint	variable		550	Ag wire/AgCl	3
6.0733.100	100	12	6	5	Glass	В	080	080	Ceramic	c(KCl)=3 mol/L		525	LL system	3
6.0750.100	125	12	12	1	Glass	В	080	080	Fixed ground-joint	c(KCl)=3 mol/L (Gel)	c(KCl)=3 mol/L	530	Ag wire/AgCl	40
6.0728.100	53	15	7	3	PCTFE	5.5 mm	060	060	Ceramic	variable		12.5	Ag wire/AgCl	3
6.0728.120	53	15	7	3	PCTFE	5.5 mm	060	060	Ceramic	c(KCl)=3 mol/L		12.5	Ag wire/AgCl	3
6.0728.130	53	15	7	3	PCTFE	5.5 mm	060	060	Ceramic	c(KCl)=3 mol/L		12.5	LL system	5
6.0730.100	65	12	6	3	Glass	5.5 mm	040	040	Ceramic	c(KCl)=3 mol/L (thickened)			LL system	20

6.09 Conductivity cells	cell constant (cm-1)	Installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Plug-in head	Туре	Temperature range short-term (°C)	Temperature range long-term (°C)	Measuring range	Temperature sensor
6.0914.040	0.1	125	12	12	35	Steel, stainless	Fixed cable 4xB (4 mm)	Steel, stainless	070	070	0300 μS/cm	Pt1000
6.0916.040	0.1	125	12	12	35	Steel, stainless	Fixed cable, plug N	Steel, stainless	070	070	0300 μS/cm	Pt1000
6.0918.040	0.1	125	12	12	35	Steel, stainless	Fixed cable, plug O	Steel, stainless	070	070	0300 μS/cm	Pt1000
6.00921.040	0.1	125	12	12	35	Steel, stainless	Fixed cable, plug N	Steel, stainless	070	070	0300 μS/cm	Pt1000
6.00922.040	0.1	125	12	12	35	Steel, stainless	Fixed cable, plug N	Steel, stainless	070	070	0300 μS/cm	Pt1000
6.0917.080	0.5	125	12	12	30	PEEK	Fixed cable, plug O	4-wire, Pt	070	070	0.015250 mS/cm	Pt1000
6.00923.080	0.5	125	12	12	30	PEEK	Fixed cable, plug N	4-wire, Pt	070	070	0.015250 mS/cm	Pt1000
6.00924.080	0.5	125	12	12	30	PEEK	Fixed cable, plug N	4-wire, Pt	070	070	0.015250 mS/cm	Pt1000
6.0915.100	0.7	125	12	12	34	Glass, PCTFE / PEEK	Fixed cable, plug N	5-ring, Pt	070	070	0.00520 mS/cm	Pt1000
6.00925.100	0.7	125	12	12	35	Glass, PCTFE / PEEK	Fixed cable, plug N	5-ring, Pt	070	070	0.00520 mS/cm	Pt1000
6.0920.100	0.7	125	12	12	34	Glass, PCTFE / PEEK	Fixed cable, plug N	5-ring, Pt	070	070	0.00520 mS/cm	Pt1000
60912110	0.8	125	12	12	35	PP	Fixed cable 4xB (4 mm)	Pt platinized	570	570	0.001100 mS/cm	Pt1000
6.0910.120	0.9	120	12	12	16	Glass	G	Pt platinized	570	570	0.001100 mS/cm	
6.0915.130	1	142	12	12	50	Glass, PCTFE / PEEK	Fixed cable, plug N	5-ring, Pt	070	070	0.005100 mS/cm	Pt1000
6.0920.130	1	142	12	12	50	Glass, PCTFE / PEEK	Fixed cable, plug N	5-ring, Pt	070	070	0.005100 mS/cm	Pt1000
6.0913.130	1.1					PP					0400 μS/cm	
6.0919.140	1.6	125	12	12	40	Glass	Fixed cable, plug O	3-ring, Pt	070	070	0.11000 mS/cm	Pt1000
6.00929.140	1.6	125	12	12	40	Glass	Fixed cable, plug N	3-ring, Pt	570	570	0.11000 mS/cm	Pt1000
6.00930.140	1.6	125	12	12	40	Glass	Fixed cable, plug N	3-ring, Pt	570	570	0.11000 mS/cm	Pt1000

6.11 Temperature sensors / Other sensors»	Installation length (mm)	Shaft diameter top (mm)	Shaft diameter bottom (mm)	Min. immersion depth (mm)	Shaft material	Shaft material (completion)	Plug-in head	Temperature range short-term (°C)	Temperature range long-term (°C)	Measuring range	Temperature sensor
6.1110.010	120	12	5	20	Glass		Fixed cable 2xB (2 mm)	-50180	-50180	-50180 °C	Pt1000
6.1110.100	125	12	5	20	Glass		G	-50180	-50180	-50180 °C	Pt1000
6.1110.110	178	12	6.4	20	Glass		G	-50180	-50180	-50180 °C	Pt1000
6.1110.120	90	12	5	20	Glass		G	-50180	-50180	-50180 °C	Pt1000
6.1111.010	175	2	2	20	Stainless steel (AISI 304)		mini-DIN			-200 300 °C	Pt100
6.1111.020	300	2	2	20	Stainless steel (AISI 304)		mini-DIN			-200 300 °C	Pt100
6.1114.010	140	12	3	10	PEEK	Tip (75 mm) stainless steel	Fixed cable 2xB (2 mm)	-50100	-50100	-50100 °C	Pt1000
6.1115.000	140	12	12	30	Glass		Fixed cable (I = 1 m) with plug F	040	040		
6.1116.000	140	12	12	35	Luran, stainless steel		Fixed cable (I = 1 m) with mini-USB	060	060	0500 % a.s.	Pt100
6.9011.020	125	12	12	15	Glass		Fixed cable, plug O			060 °C	
6.01117.300	125	12	12	15	Glass		Q			060 °C	
6.9011.040	125	12	12	15	Glass	PVDF coated glass	Fixed cable, plug O			060 °C	

RDE/RRDE tips/ microelectrodes (6.12, PT., RDE., RRDE.)	Length (mm)	Shaft diameter (mm)	Shaft material	Disk diameter (mm)	Disk material	Ring width (mm)	Ring material	Gap between ring and disk (mm)	Connector	Temperature range short-term (°C)	Temperature range long-term (°C)
6.1204.130	52.5	7	PEEK	2	Ag				M3	040°C	040°C
6.1204.140	52.5	7	PEEK	2	Au				M3	040°C	040°C
6.1204.170	52.5	7	PEEK	3	Pt				M3	040°C	040°C
6.1204.190	52.5	7	PEEK	2	Pt				M3	040°C	040°C
6.09395.014	52.5	10	PEEK	3	GC				M4	040°C	040°C
6.09395.024	52.5	10	PEEK	3	Pt				M4	040°C	040°C
6.09395.034	52.5	10	PEEK	3	Au				M4	040°C	040°C
6.09395.044	52.5	10	PEEK	3	Ag				M4	040°C	040°C
6.1204.600	52.5	8	Glass	2	GC				M3	050°C	050°C
6.1204.610	52.5	8	Glass	1	Pt .				M3	050°C	050°C
AG.100	52	2	Glass	0.1	Ag				B (2 mm, male)	040°C	040°C
AG.25	52	2	Glass	0.025	Au				B (2 mm, male)	040°C	040°C
AG.300 AG.300	52 52	2	Glass	0.03	Ag				B (2 mm, male)	040°C 040°C	040°C
AU.10	52	2	Glass	0.3	Ag Au				B (2 mm, male) B (2 mm, male)	040°C	040°C
AU.100	52	2	Glass	0.01	Au				B (2 mm, male)	040°C	040°C
AU.200	52	2	Glass	0.2	Au				B (2 mm, male)	040°C	040°C
AU.25	52	2	Glass	0.025	Au				B (2 mm, male)	040°C	040°C
AU.300	52	2	Glass	0.3	Au				B (2 mm, male)	040°C	040°C
AU.40	52	2	Glass	0.04	Au				B (2 mm, male)	040°C	040°C
AU.50	52	2	Glass	0.05	Au				B (2 mm, male)	040°C	040°C
AU.500	52	2	Glass	0.5	Au				B (2 mm, male)	040°C	040°C
IR.75	52	2	Glass	0.075	Ir				B (2 mm, male)	040°C	040°C
PD.100	52	2	Glass	0.1	Pd				B (2 mm, male)	040°C	040°C
PD.25	52	2	Glass	0.025	Pd				B (2 mm, male)	040°C	040°C
PD.300	52	2	Glass	0.3	Pd				B (2 mm, male)	040°C	040°C
PD.500	52	2	Glass	0.5	Pd				B (2 mm, male)	040°C	040°C
PT.10	52	2	Glass	0.01	Pt				B (2 mm, male)	040°C	
PT.100	52	2	Glass	0.1	Pt				B (2 mm, male)	040°C	
PT.20	52	2	Glass	0.02	Pt				B (2 mm, male)	040°C	
PT.200	52	2	Glass	0.2	Pt				B (2 mm, male)	040°C	
PT.25	52	2	Glass	0.025					B (2 mm, male)	040°C	
PT.300	52	2	Glass	0.023	Pt				B (2 mm, male)	040°C	
PT.50	52	2	Glass	0.05	Pt				B (2 mm, male)	040°C	
PT.500	52	2	Glass	0.03	Pt				B (2 mm, male)	040°C	
RRDE.PTPT.S	51.5	11.5	PEEK	5	Pt	0.375	Pt	0.375	M4	040°C	
RRDE.AUPT.S	51.5	11.5	PEEK	5	Au	0.375	Pt	0.375	M4	040°C	
RRDE.GCPT.S	51.5	11.5	PEEK	5	GC	0.375	Pt	0.375	M4	040°C	
3500003900	52.5	10	PEEK	5	GC				M4	040°C	
RDE.PT50.S	52.5	10	PEEK	5	Pt				M4	040°C	
RDE.AU50.S	52.5	10	PEEK	5	Au				M4	040°C	
RDE.AG50.S	52.5	10	PEEK	5	Ag				M4	040°C	
RDE.CU50.S	52.5	10	PEEK	5	Cu				M4	040°C	
RDE.ZN50.S	52.5	10	PEEK	5	Zn				M4	040°C	040°C
RDE.STEEL.S	52.5	10	PEEK	5	Stainless steel				M4	040°C	040°C
RDE.BLANK.S	52.5	10	PEEK	5	Empty				M4	040°C	040°C

Screen-printed electrodes and interdigitated electrodes/micro-electrodes/micro-electrodes	WE material	WE dimension (mm)	AUX material	Cell configuration	Substrate size (mm)	Substrate material	Units/box
110	Carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 units
11L	Carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 units
C110	Carbon	d. 4	Carbon	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
C110-NTC	Carbon	d. 4	Carbon	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
C11L	Carbon	d. 4	Carbon	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
150	Carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 unit
220AT	Gold AT	d. 4	Platinum	Standard	34 x 10 x 0.5	Ceramic	75 unit
220BT	Gold BT	d. 4	Gold BT	Standard	34 x 10 x 0.5	Ceramic	75 unit
C220AT	Gold AT	d. 4	Gold AT	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
C220BT	Gold BT	d. 1.6	Gold BT	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
C223AT	Gold AT	d. 1.6	Gold AT	1.6 d	34 x 10 x 0.5	Ceramic	75 unit
C223BT	Gold BT	d. 4	Gold BT	1.6 d	34 x 10 x 0.5	Ceramic	75 units
250AT	Gold AT	d. 4	Platinum	Standard	34 x 10 x 0.5	Ceramic	75 unit
250BT	Gold BT	d. 4	Platinum	Standard	34 x 10 x 0.5	Ceramic	75 unit
410	Co-phthalocyanine/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 unit
550	Platinum	d. 4	Platinum	Standard	34 x 10 x 0.5	Ceramic	75 unit
550BT	Platinum BT	d. 4	Platinum BT		34 x 10 x 0.5	Ceramic	75 unit
C550	Platinum	d. 4	Carbon	Work in solution	34 x 10 x 0.5	Ceramic	75 unit
610	Meldola's blue/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 unit
710	Prussian blue/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 units
F10	Ferrocyanide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 units
810	Ruthenium oxide	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
010	Silver	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	75 unit
C013	Silver	d. 1.6	Carbon	1.6 d	34 x 10 x 0.5	Ceramic	75 unit
110AGNP	Silver nanoparticles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110ALI	Alizarin/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110AQU	Anthraquinone/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110AUP	Gold particles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110BI	Bismuth oxide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CNF	Carbon nanofibres/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CNF-GNP	Carbon nanofibres-gold Nanoparticles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CNT	Multi-walled carbon nanotubes/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CNT-GNP	Multi-walled carbon Nanotubes-gold nanoparticles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CSQD	Core-shell quantum dots ZnS/CdSe/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110CUPH	Copper (II) phthalocyanine/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110FEPH	Iron (II) phthalocyanine/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110FERRI	Potassium Ferricyanide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110GNP 110GNP-STR	Gold nanoparticles/carbon Streptavidin modified gold	d. 4 d. 4	Carbon	Standard Standard	34 x 10 x 0.5	Ceramic	50 unit
	Nanostructured/carbon						
110GPH	Graphene/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110GPH-GNP	Graphene-gold nanoparticles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110GPHOX	Graphene oxide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
110IRP	Iridium Particles /carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 unit
140146	1						
110MC 110NI	Mesoporous carbon/carbon Nickel oxide/carbon	d. 4 d. 4	Carbon Carbon	Standard Standard	34 x 10 x 0.5 34 x 10 x 0.5	Ceramic Ceramic	50 unit

Screen-printed electrodes and interdigitated electrodes/micro- electrodes	WE material	WE dimension (mm)	AUX material	Cell configuration	Substrate size (mm)	Substrate material	Units/box
110PANI	Polyaniline/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110PDP	Palladium particles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110PHEN	Phenanthroline/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110PPYR	Polypyrrole/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110PTP	Platinum particles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110PLYS	Poly-L-Lysine/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110QD	Core quantum dots CdSe/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110RGPHOX	Reduced graphene oxide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110RHP	Rhodium Particles/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110SFT	Surface treated/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110STR	Streptavidin/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110SWCNT	Single-walled carbon nanotubes/ Carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
110XTR	Extravidin/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
4W110	Carbon	d. 2.95	Carbon	4WE sharing AUX and REF	38 x 20 x 1.0	Ceramic	20 units
8W110	Carbon	d. 2.95	Carbon	8WE sharing AUX and REF	38 x 20 x 1.0	Ceramic	20 units
8X110	Carbon	d. 2.56	Carbon	Array of 8X ectrochemical cells	34 x 79 x 1.0	Ceramic	20 units
8X110STR	Streptavidin/Carbon	d. 2.95	Carbon	Array of 8X ectrochemical cells	34 x 79 x 1.0	Ceramic	10 units
8X220AT	Gold AT	d. 2.56	Gold	Array of 8X ectrochemical cells	34 x 79 x 1.0	Ceramic	20 units
96X110	Carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	4 plates
96X110CNT	Multi-walled carbon nanotubes/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110CNT-GNP	Multi-walled carbon Nanotubes-gold nanoparticles/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110GNP	Gold nanoparticles/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110GNP-STR	Streptavidin modified gold Nanostructured/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110STR	Streptavidin/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110SWCNT	Single-walled carbon nanotubes/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X110XTR	Extravidin/carbon	d. 3	Carbon	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	2 plates
96X220	Gold	d. 3	Gold	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	4 plates
96X550	Platinum	d. 3	Platinum	ELISA 96X electrochemical cells	74 x 110 x 5.0	PCB	4 plates
X1110	Carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	75 units
X4410	Co-phthalocyanine/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	75 units
XFF10	Ferrocyanide/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	75 units
X1110AGNP	Silver nanoparticles/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110ALI	Alizarin/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110AQU	Anthraquinone/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110AUP	Gold particles/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110BI	Bismuth oxide/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CNF	Carbon nanofibres/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CNF-GNP	Carbon nanofibres-gold Nanoparticles/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CNT	Multi-walled carbon nanotubes/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CNT-GNP	Multi-walled carbon nanotubes- Gold nanoparticles/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CUPH	Copper (II) phthalocyanine/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110CSQD	Core-shell quantum dots ZnS/CdSe/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110FEPH	Iron (II) phthalocyanine/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110FERRI	Potassium Ferricyanide/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units

Screen-printed electrodes and interdigitated electrodes/micro- electrodes	WE material	WE dimension (mm)	AUX material	Cell configuration	Substrate size (mm)	Substrate material	Units/box
X1110GNP	Gold nanoparticles/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110GNP-STR	Streptavidin modified Gold nanostructured/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110GPH	Graphene/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110GPH-GNP	Graphene-gold nanoparticles/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110GPHOX	Graphene oxide/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110IRP	Iridium Particles /carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110MC	Mesoporous carbon/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110NI	Nickel oxide/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X11100MC	Ordered mesoporous carbon/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PANI	Polyaniline/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PDP	Palladium particles/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PHEN	Phenanthroline/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PPYR	Polypyrrole/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PTP	Platinum particles/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110PLYS	Poly-L-Lysine/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110QD	Core quantum dots CdSe/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110RGPHOX	Reduced graphene Oxide/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110RHP	Rhodium Particles/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110SFT	Surface treated/carbon	d. 4	Carbon	WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110STR	Streptavidin/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110SWCNT	Single-walled carbon nanotubes/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
X1110XTR	Extravidin/carbon	d. 4	Carbon	2WE sharing AUX and REF	34 x 10 x 0.5	Ceramic	50 units
GLU10	Glucose	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
LACT10	Lactate	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
JA10	Uric Acid	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	50 units
ITO10	ITO	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent Plastic	20 units
P10	PEDOT	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent Plastic	75 units
AGTR10	Transparent silver	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent plastic	50 units
AUTR10	Transparent Gold	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent Plastic	20 units
COTE10	Transparent Carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent Plastic	20 units
AL10	Aluminium	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
BIOFV1	PEDOT	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent Plastic	75 units
CR10	Chromium	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
MO10	Molybdenum	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
PB10	Lead	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
SB10	Antimony	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
5N10	Tin	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
TA10	Tantalum	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
W10	Tungsten	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	20 units
PW-AU10	Gold	d. 4	Carbon	Standard	34 x 10 x 0.5	White Plastic	20 units
** 7010	Silver	d. 4	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units

Screen-printed electrodes and interdigitated electrodes/micro-electrodes	WE material	WE dimension (mm)	AUX material	Cell configuration	Substrate size (mm)	Substrate material	Units/box
TLFCL110	Carbon	12.6 mm ²	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
TLFCL110-CIR	Carbon	d. 4	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
TLFCL110S	Carbon	2 mm ²	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
TLFCL1110	Carbon Gold AT	2 mm ² x 2	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
TLFCL210AT-CIR TLFCL210BT-CIR	Gold BT	d. 4 d. 4	Carbon Carbon	Thin layer flow-cell incorporated Thin layer flow-cell incorporated	80 x 25 x 0.1 80 x 25 x 0.1	Ceramic Ceramic	10 units
		0.8 x 5 (external)					
TLFCL2222AT	Gold AT	0.4 x 5 (internal)		Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
TLFCL510-CIR	Platinum	d. 4	Carbon	Thin layer flow-cell incorporated	80 x 25 x 0.1	Ceramic	10 units
SPESMIX	Mix: Carbon, gold, platinum	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	100 units
AUMIX	Mix: Gold	d.4 and d.1.6	Gold	Standard, 1.6 d and work in solution	34 x 10 x 0.5	Ceramic	100 units
MEDIATORSPES	Mix: Co-phthalocyanine, meldola's blue,prussian blue, ferrocyanide/carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Ceramic	100 units
OTEMIX	Mix: ITO, PEDOT, transparent gold and transparent carbon	d. 4	Carbon	Standard	34 x 10 x 0.5	Transparent plastic	40 units
AUFET30	Gold	30 microns lines and gaps	Gold	Field effect transistor	34 x 10 x 0.7	White plastic	50 units
G-IDE222	Gold	10 microns lines and gaps	Gold	Interdigitated with AUX and REF	22 x 7.0 x 0.7	Glass	20 units
G-IDE555	Platinum	10 microns lines and gaps	Platinum	Interdigitated with AUX and REF	22 x 7.0 x 0.7	Glass	20 units
G-IDEAG5	Silver	5 microns lines and gaps	Silver	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDEAU10	Gold	10 microns lines and gaps	Gold	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDEAU5	Gold	5 microns lines and gaps	Gold	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDECU5	Copper	5 microns lines and gaps	Copper	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDEPT10	Platinum	10 microns lines and gaps	Platinum	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDEPT5	Platinum	5 microns lines and gaps	Platinum	Interdigitated squared format	22 x 7.0 x 0.7	Glass	20 units
G-IDEMIX	Mix: Gold, platinum	5 and 10 microns lines and gaps	Mix: gold, platinum	Interdigitated:concentric; squared	22 x 7.0 x 0.7	Glass	20 units
G-IDECONAU10	Gold	10 microns lines and gaps	Gold	Interdigitated concentric format	22 x 7.0 x 0.7	Glass	20 units
G-IDECONPT10	Platinum	10 microns lines and gaps	Platinum	Interdigitated concentric format	22 x 7.0 x 0.7	Glass	20 units
G-MEA222	Gold	d. 3 / Microholes 10 microns	Gold	Interdigitated microelectrode array d. 3mm	22 x 7.0 x 0.7	Glass	20 units
G-MEA555	Platinum	d. 3 / Microholes 10 microns	Platinum	Interdigitated microelectrode array d. 3mm	22 x 7.0 x 0.7	Glass	20 units
G-MEAB222	Gold	Microband 10 microns	Gold	Interdigitated microelectrode array	22 x 7.0 x 0.7	Glass	20 units
G-MEAB555	Platinum	Microband 10 microns	Platinum	Interdigitated microelectrode array	22 x 7.0 x 0.7	Glass	20 units
IDEAU200	Gold	200 microns lines and gaps	Gold	Interdigitated squared format	22 x 7.0 x 0.7	Ceramic	20 units
IDEAU200-HPT-WB	Gold	200 microns lines and gaps	Gold	Squared design / Pt heater	22 x 7.0 x 0.7	Ceramic	50 units

Screen-printed electrodes and interdigitated electrodes/micro-electrodes	WE material	WE dimension (mm)	AUX material	Cell configuration	Substrate size (mm)	Substrate material	Units/box
P-IDEAG50	Silver	50 microns lines and gaps	Silver	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
P-IDEAG100	Silver	100 microns lines and gaps	Silver	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
P-IDEAU50	Gold	50 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
P-IDEAU100	Gold	100 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
P-IDEITO50	ITO	50 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
P-IDEITO100	ITO	100 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	Transparent plastic	50 units
PW-IDEAU50	Gold	50 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	White plastic	50 units
PW-IDEPD50	Palladium	50 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	White plastic	50 units
PW-IDEAU100	Gold	100 microns lines and gaps	Gold	Interdigitated squared format	23 x 7.0 x 0.2	White plastic	50 units
PW-IDEPD100	Palladium	100 microns lines and gaps	Palladim	Interdigitated squared format	23 x 7.0 x 0.2	White plastic	50 units
PW-4XIDEAU20	Gold	20 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units
PW-4XIDEAU30	Gold	30 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units
PW-4XIDEAU40	Gold	40 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units
PW-4XIDEAU50	Gold	50 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units
PW-4XIDEAU60	Gold	60 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units
PW-4XIDEAU70	Gold	70 microns lines and gaps	Gold	Multi-interdigitated	16 x 16 x 0.2	White plastic	50 units

Metrohm Electrode Care

Be sure that your electrode is still up to its job

Enjoy peace of mind with 142 **Metrohm Electrode Care**

whether your pH electrodes perform according to specifications. To this end, we check crucial parameters such and precision of the equivalence point, the response as response time, drift behavior, and slope of your electime, and the potential jump are compared against tartrode against target values. Based on the results, the get values. If one or several parameters fail this test, a performance of your electrode is rated «excellent», replacement of the electrode is highly recommended. «good», «usable», or «poor». In case of a «usable» or

«poor» result, a replacement is highly recommended in order to reduce the risk of OOS results. Other test meth-«Electrode Care» is a service we provide to confirm ods, which are based on standardized titrations, are available for metal and surfactant electrodes. Accuracy



Highlights

- Recognize poor electrode quality before your analysis fails
- Omit tedious error tracking
- Get a replacement recommendation based on real data
- Increase accuracy and the precision of your results
- Reduce sample waste and tedious analysis repetition due to incorrect results
- Reduce downtime of your system

Nobody can afford Out-of-Specification (OOS) investigations!

Ask your Metrohm Sales Rep for Metrohm Electrode Care and be sure your electrodes are always up to their job.

For further information, please consult our dedicated Application Bulletins.

- Metrohm Application Bulletin AB-048 check of metal electrodes
- Metrohm Application Bulletin AB-188 pH measurement technique
- Metrohm Application Bulletin AB-305 check of surfactant electrode

www.metrohm.com

