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CO₂-Adsorberkartusche CW	1
Übersicht	1
Vorbereitung	1
Installation	2
Pflege und Wartung	3
CO₂ adsorption cartridge CW	4
Overview	4
Preparation	4
Installation	5
Care and maintenance	5
Cartouche d'adsorption de CO₂ CW	6
Aperçu	6
Préparation	6
Installation	7
Entretien et maintenance	8
Cartucho de adsorción de CO₂ CW	8
Visión conjunta	8
Preparación	8
Instalación	9
Conservación y mantenimiento	10

DE

CO₂-Adsorberkartusche CW

Übersicht

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CO₂ adsorption cartridge CW

Overview

In order to efficiently remove disturbing CO₂ from the eluent, the Metrohm CO₂ Suppressor (MCS) aspirates air through the CO₂ adsorption cartridge CW (6.2837.100). The CO₂ adsorption cartridge CW is filled with cotton and a CO₂ adsorber material.

You can find the CO₂ adsorption cartridge CW and all necessary parts in the accessory kit: Vario/Flex SeS (6.5000.020).

Preparation

CAUTION

Insufficient CO₂ adsorption

In order to achieve an optimal CO₂ adsorption, the CO₂ adsorption cartridge CW must be put into operation correctly.

Proceed in accordance with the following instructions.


Required accessories

- CO₂ adsorption cartridge CW with dust filter (6.2837.100)


Preparing the CO₂ adsorption cartridge CW

Prepare the CO₂ adsorption cartridge CW for use as follows:


1. Remove the protective cap from the tip of the CO₂ adsorption cartridge CW.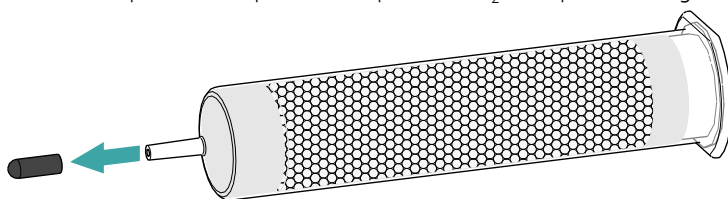

2. Attach the dust filter to the tip of the CO₂ adsorption cartridge CW.

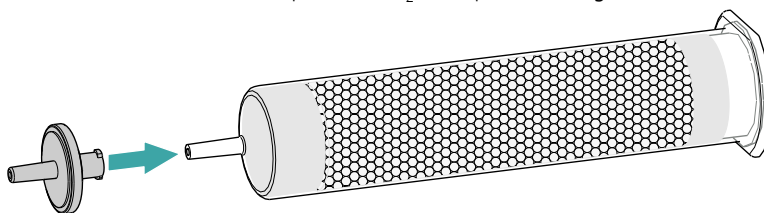3. Remove the label from the lid of the CO₂ adsorption cartridge CW.

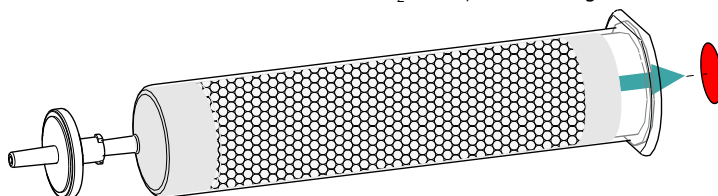This uncovers the small opening in the lid of the CO₂ adsorption cartridge CW, through which air is aspirated. The CO₂ adsorption cartridge CW is now ready for installation.

NOTE

The new CO₂ adsorption cartridge works **without** attached H₂O adsorption cartridge.

Installation


Required accessories

- Adsorption cartridge holder (6.2057.080)
- Prepared CO₂ adsorption cartridge CW (6.2837.100)


Installing the CO₂ adsorption cartridge CW

Install the CO₂ adsorption cartridge CW as follows:


1. Push two clips into the slots on the adsorption cartridge holder.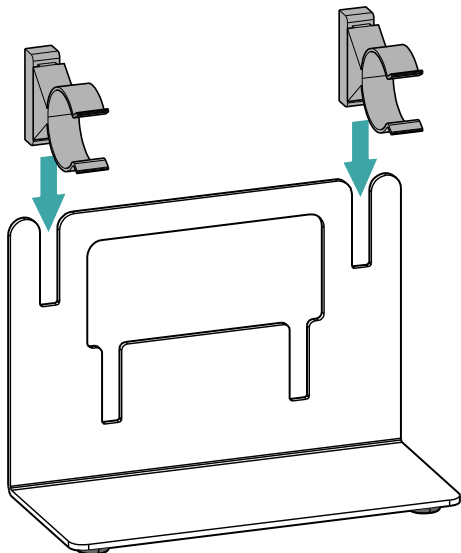

2. Attach the capillary connected to the Metrohm CO₂ Suppressor's (MCS) **Air in** connection to the tip of the CO₂ adsorption cartridge CW.

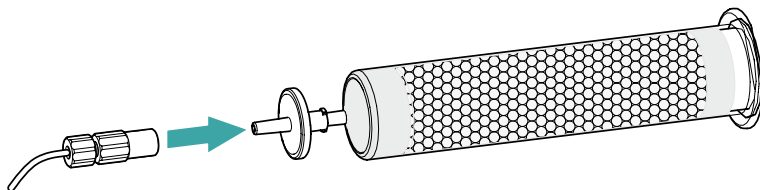3. Clamp the CO₂ adsorption cartridge CW in place using the clips on the adsorption cartridge holder.

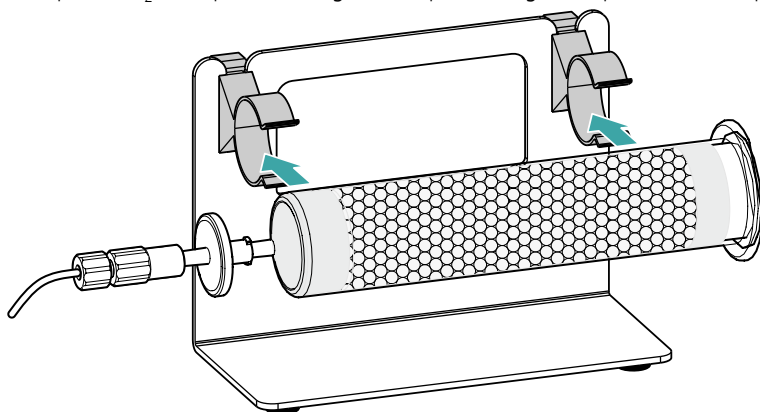4. Place the adsorption cartridge holder with the CO₂ adsorption cartridge CW into the instrument's detector chamber.

Care and maintenance

The CO₂ adsorption cartridge CW needs to be optically checked at regular intervals. Depleted cartridges change their color to violet. New cartridges are white.

Maintenance interval

We recommend replacing the CO₂ adsorption cartridge CW at least once a year, irrespective of the degree of color change.

The order number for a new cartridge is 6.2837.100.

Removing the depleted CO₂ adsorption cartridge CW

1. Remove the depleted CO₂ adsorption cartridge CW from the adsorption cartridge holder.
2. Remove the **Air in** capillary of the Metrohm CO₂ Suppressor (MCS).
3. Professionally dispose of the depleted CO₂ adsorption cartridge CW.

Installing the new CO₂ adsorption cartridge CW

1. Prepare the new CO₂ adsorption cartridge CW (see title "Preparation").
2. Install the new CO₂ adsorption cartridge CW (see title "Installation").

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