

1 Purpose

This document describes the innovations and fixed bugs in the following new firmware version:

- 5.885.0030

in comparison with the firmware version:

- 5.885.0013

2 What's new

Topic	Description
New processor	The device runs with a new processor. Devices with firmware version 5.885.0013 or older can therefore not be updated to firmware version 5.885.0030. A new device with a new processor is required for using firmware version 5.885.0030.
Methods and backups	Methods and backups which were created on program version 5.885.0030 are not reverse-compatible.
USB Plug & Play	USB flash drives and printers can be connected or removed at any time. The device does not need to be switched off before a USB flash drive or a printer is connected or removed.
RS-232/USB Box Version 2.0	There is a new version of the RS-232/USB Box 6.2148.030. Starting with program version 5.885.0030, Version 2.0 of the RS-232/USB Box must be used. Version 2.0 of the RS-232/USB Box is identified with a sticker accordingly.
Languages installed	All available languages are installed on the device. No additional language files need to be loaded. The following languages are available: <ul style="list-style-type: none">▪ German▪ English▪ French▪ Spanish▪ Portuguese▪ Italian

Topic	Description
Portuguese dialog texts	The dialog texts are now available in Portuguese.
Program files directory	New program files must be stored in the 885 ► SwUpdates directory.

3 Fixed bugs

Topic	Description
Changing the contrast	If the contrast was changed, it would be reset back to the default value when the device was restarted.

4 Known issues

Topic	Description
Restoration after backup	In firmware version 5.885.0030, if a backup from a lower firmware version is restored, then the settings under System ► External devices and System ► Lift will not be applied.