



Application Note AN-COR-019

# Determining the corrosion rate with INTELLO

## Tafel analysis and more

Tafel analysis is an important electrochemical technique used to understand reaction kinetics. Studying the Tafel slope reveals the rate-determining steps in electrode reactions, aiding fields like corrosion and fuel cell research. This method helps industries optimize processes and improve device performance by tailoring materials and conditions for greater efficiency. Regarding corrosion, Tafel analysis provides insights into the corrosion rate and mechanisms of different metals in various environments. By examining the Tafel slope, researchers can

determine the polarization resistance and corrosion rate as well as the corrosion current and potential, shedding light on how materials interact with their surroundings. This analysis aids industries in selecting the appropriate coatings, inhibitors, and materials to mitigate corrosion, extend the lifespan of structures, and ensure the integrity of metal components in challenging environments.

This Application Note illustrates Tafel analysis in INTELLO with aluminum in artificial seawater.

## EXPERIMENTAL


A 250 mL three-electrode corrosion cell was constructed for this study. An Ag/AgCl reference electrode was used along with two stainless steel rod-shaped counter electrodes, and an aluminum disc mounted on a sample holder served as the working electrode. The electrolyte was artificial seawater (3.5% NaCl).

The default procedure in INTELLO (Linear Polarization – Tafel Analysis) was chosen. First, an open circuit potential (OCP) measurement was made, and then -30 mV vs OCP was applied. A linear sweep to +20 mV vs OCP was programmed. A scan rate of 50 mV/s was used.

## RESULTS


The result of the linear polarization experiment ( $j$  vs  $E$  plot) is shown in **Figure 1**. Note that the

normalized current (current density) is plotted by default.

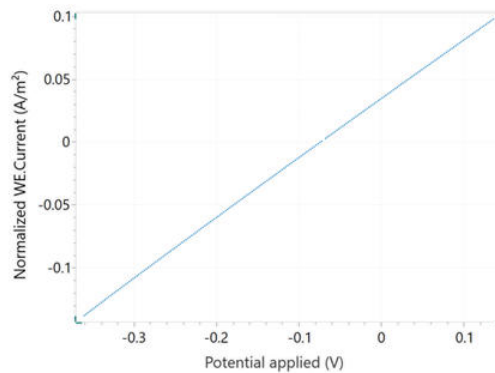

**Figure 1.** Current vs potential plot of aluminum in seawater.


A Tafel plot,  $\text{Log}(I)$  vs  $E$ , is also shown in **Figure 2**.

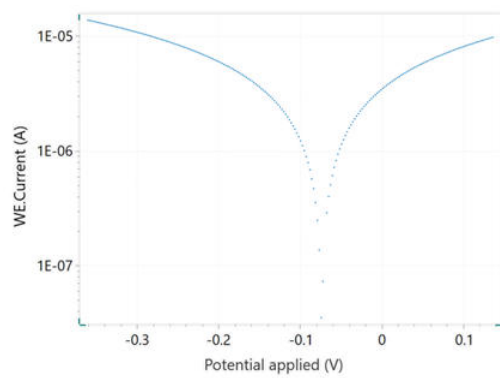

**Figure 2.** Tafel plot which corresponds to the  $j$  vs  $E$  curve shown in Figure 1.

When selecting the Tafel analysis command, it is possible to specify the density of the sample, its equivalent weight, and the surface area (**Figure**


3). If these variables are already known, they can also be added in the main parameters window prior to beginning the experiment.

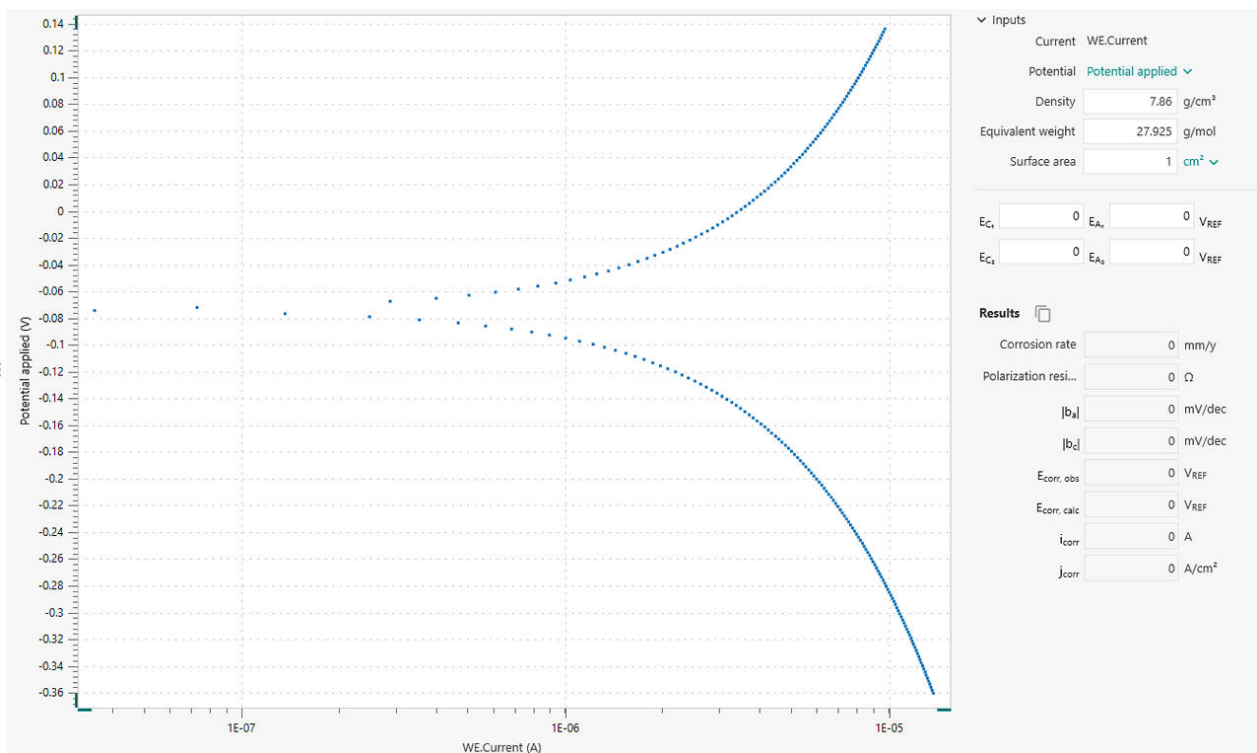

**Figure 3.** Tafel analysis window in INTELLO.

In order to perform a corrosion rate analysis, the linear regions of the anodic and cathodic side of the Tafel plot must be specified. This is accomplished with the help of four markers: two each for the anodic side and the cathodic side. Such markers define the lines for the linear regressions, according to the Tafel equations below.

Once the regions are selected, the linear regressions appear in the plot. The positions can be readjusted by dragging the markers to new positions on the plot. The corrosion potential and the corrosion current correspond to the X and Y coordinates, respectively, of the intersection of the two regression lines (**Figure 4**).

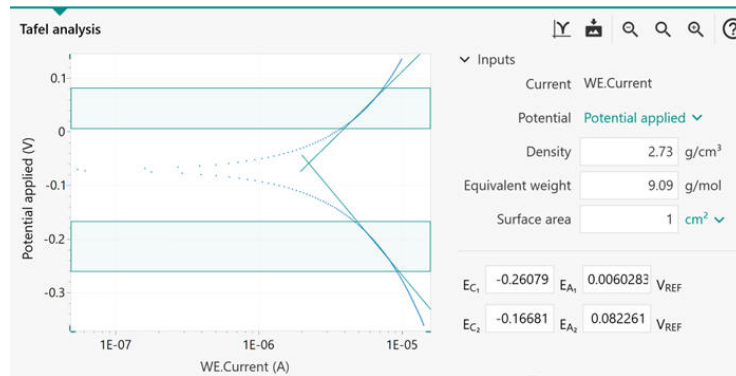

**Figure 4.** Tafel analysis window with the linear regions identified in the shaded area between the horizontal markers. The green angled lines indicate linear regressions.

The results of the Corrosion Rate Analysis command are listed in **Figure 5**. The Tafel

analysis provides a quick estimation of the corrosion rate and the polarization resistance.

Results	
Corrosion r...	0.024289 mm/y
Polarization...	28037 $\Omega$
$ b_a $	0.26199 mV/dec
$ b_c $	0.31993 mV/dec
$E_{corr, obs}$	-0.070577 $V_{REF}$
$E_{corr, calc}$	-0.058915 $V_{REF}$
$i_{corr}$	2.2308E-06 A
$j_{corr}$	2.2308E-06 A/cm <sup>2</sup>

**Figure 5.** Results of the corrosion rate analysis.

The corrosion rate ( $R_M$ , mm/year) is calculated in the following equation using the corrosion

current,  $i_{corr}$ .

$$R_M = 3.17E - 9 \frac{M}{nF\rho A} i_{corr}$$

3.17E - 9: conversion factor from  $\text{cm s}^{-1}$  to  $\text{mm year}^{-1}$  **M** ( $\text{g mol}^{-1}$ ): atomic weight of the sample  
**n**: number of electrons exchanged in the reaction **p** ( $\text{g cm}^{-3}$ ): density of the sample **F** ( $96485 \text{ C mol}^{-1}$ ): Faraday constant **A** ( $\text{cm}^2$ ): area

of the sample

The ratio  $M/n$  is also named equivalent weight.

To calculate  $i_{corr}$  the Butler-Volmer equation is employed, as shown here.

$$i_{corr} \left( e^{2.303 \frac{\eta}{b_a}} - e^{-2.303 \frac{\eta}{b_c}} \right)$$

$i_{corr}$ : corrosion current (i.e., the rate of charge transfer at the corrosion potential  $E_{corr}$ ) **2.303**: conversion factor between  $\log$  and  $\log_e$  and  $\log_{10}$   **$\eta$**  (V): overpotential, defined as the difference between applied potential  $E$  and the

corrosion potential  $E_{corr}$   **$b_a$**  (V): Tafel slope of the anodic branch  **$b_c$**  (V): Tafel slope of the cathodic branch

For large anodic overpotentials,

$$\left(\frac{\eta}{b_a} \gg 1\right)$$

---

the Butler-Volmer equation simplifies to the Tafel equation for the anodic reaction:

$$\eta = b_a \cdot \log\left(\frac{i}{i_{corr}}\right)$$

---

Analogously, for large cathodic overpotentials,

$$\left(\frac{\eta}{b_c} \ll -1\right)$$

---

the Tafel equation for the cathodic reaction is given by:

$$\eta = -b_c \cdot \log \left( \frac{i}{i_{corr}} \right)$$

The Tafel equations predict a straight line for the variation of the logarithm of the current with the potential. Therefore, currents are often shown in semi-logarithmic plots, known as Tafel plots. Care must be taken when choosing the regions for Tafel slope analysis. A correct estimate of the Tafel slopes is possible only if the linear Tafel region (the shaded regions between each pair of markers in **Figure 3**) covers at least one decade in current.

In certain cases, the Tafel slope analysis is not possible. For example, under the diffusion control, the reactants are hampered in reaching the sample surface (e.g., in stagnant solutions). As a consequence, the cathodic reaction does not occur, resulting in  $b_c = \infty$ . Similarly, under passivation conditions, the surface of the sample is coated with a protective layer. This prevents the anodic reaction from occurring, resulting in  $b_a = \infty$ .

## CONTACT

瑞士万通中国  
北京市海淀区上地路1号院  
1号楼7702  
100085 北京

marketing@metrohm.com.cn

## CONFIGURATION



### VIONIC

VIONIC 是我的新一代恒位/恒流,由 Autolab 的新 INTELLO 件提供力。

相比目前市上任何一器,VIONIC 提供了**最全面**的合格

。

- 从:  $\pm 50$  V
- 准流  $\pm 6$  A
- EIS 率: 10 MHz
- 采隔: 低至  $1 \mu$ s

VIONIC 的价格中通常包括大多数其他器生的外用,例如:

- 化学阻抗 (EIS)
- 可浮配件
- 第二量 (S2)
- 模描